



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Data Sheet

BRD3 (BD2) Inhibitor Screening Assay Kit

Catalog # 32523

DESCRIPTION: The *BRD3 (BD2) Inhibitor Screening Assay Kit* is designed to measure the inhibition of BRD3 bromodomain 2 (BD2) from binding to its substrate. The *BRD3 (BD2) Inhibitor Screening Assay Kit* comes in a convenient AlphaLISA[®] format, with biotinylated histone peptide substrate, assay buffer, detection buffer and purified, GST-tagged BRD3 BD2 to perform a total of 384 enzyme reactions. The key to the *BRD3 (BD2) Inhibitor Screening Assay Kit* is the highly specific binding of the BRD3 bromodomain 2 to the acetylated histone substrate. With this kit, only three simple steps on a microtiter plate are required. First, a sample containing BRD3 bromodomain 2 and an inhibitor of choice is incubated with the biotinylated substrate for thirty minutes. Next, acceptor beads are added, then donor beads, followed by reading the Alpha-counts.

COMPONENTS:

Catalog #	Component	Amount	Storage	
31033	BRD3 (306-417), BD2, GST-tag	20 µg	-80 °C	(Avoid freeze/thaw cycles!)
	BET Bromodomain Ligand	400 µl	-80 °C	
	Non-acetylated Ligand	200 µl	-80 °C	
33001	3x BRD Homogeneous Assay Buffer 1	4 ml	-20 °C	
33002	3x BRD Homogeneous Detection Buffer 1	3 ml	-20 °C	

MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

AlphaLISA GSH acceptor beads, 5 mg/ml (PerkinElmer #AL109C)
 AlphaScreen Streptavidin-conjugated donor beads, 5 mg/ml (PerkinElmer #6760002S)
 Optiplate -384 (PerkinElmer #6007290)
 AlphaScreen microplate reader
 Adjustable micropipettor and sterile tips

APPLICATIONS: Useful for the study of bromodomain binding assays, screening inhibitors and selectivity profiling.

CONTRAINDICATIONS: Only limited amounts of DMSO can be included, as it has been shown to disrupt BRD-ligand interaction. Avoid green and blue dyes that absorb light in the AlphaScreen signal emission range (520-620 nm), such as Trypan Blue. Avoid the use of the potent singlet oxygen quenchers such as sodium azide (NaN₃) or metal ions (Fe²⁺, Fe³⁺, Cu²⁺, Zn²⁺ and Ni²⁺). The presence of >1% RPMI 1640 culture medium leads to a signal reduction due to the presence of excess biotin and iron in this medium. MEM, which lacks these components, does not affect AlphaScreen assays.

STABILITY: At least one year from date of receipt when stored as directed.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
 Or you can Email us at: info@bpsbioscience.com
 Please visit our website at: www.bpsbioscience.com

REFERENCE: Muller, S., Filippakopoulos, P., Knapp, S., *Expert Rev. Mol. Med.* 2011 Sep 13;13:e29.

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate. Use slow shaking for all incubations.

Step 1:

- 1) Prepare the master mixture: N wells × (2.5 µl **3x BRD Homogeneous Assay Buffer 1** + 1 µl **BET Bromodomain Ligand** + 1.5 µl **H₂O**).
- 2) Thaw **BRD3 (BD2)** on ice. Upon first thaw, briefly spin tube containing protein to recover full content of the tube. Aliquot both proteins into single use aliquots. Store remaining undiluted protein in aliquots at -80°C immediately. *Note: **BRD3** is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted protein.*
- 3) Dilute **BRD3 (BD2)** in **1x BRD Homogeneous Assay Buffer 1** at 16 ng/µl. Keep diluted proteins on ice until use. Discard any unused diluted protein after use.

Add 5 µl of master mixture to each well designated for the “Positive Control”, “Test Inhibitor”, and “Blank”. For the “Substrate Control”, add 2.5 µl **3x BRD Homogeneous Assay Buffer 1** + 1 µl **Non-acetylated Ligand** + 1.5 µl **H₂O**.

	Blank	Substrate Control	Positive Control	Test Inhibitor
3x BRD Homogeneous Assay Buffer 1	2.5 µl	2.5 µl	2.5 µl	2.5 µl
BET Bromodomain Ligand	1 µl	-	1 µl	1 µl
Non-acetylated Ligand	-	1 µl	-	-
H ₂ O	1.5 µl	1.5 µl	1.5 µl	1.5 µl
Test Inhibitor/Activator	-	-	-	2.5 µl
Inhibitor buffer (no inhibitor)	2.5 µl	2.5 µl	2.5 µl	-
1x BRD Homogeneous Assay Buffer 1	2.5 µl			
BRD3 (BD2) (16 ng/µl)*	-	2.5 µl	2.5 µl	2.5 µl
Total	10 µl	10 µl	10 µl	10 µl

- 4) Add 2.5 µl of **inhibitor solution** to each well designated “Test Inhibitor”. For the “Positive Control”, “Substrate Control” and “Blank”, add 2.5 µl of the same **solution without inhibitor** (inhibitor buffer). *Note: Keep DMSO concentration below 0.5 %.*
- 5) Add 2.5 µl of **1x BRD Homogeneous Assay Buffer 1** to the well designated “Blank”.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
 Or you can Email us at: info@bpsbioscience.com
 Please visit our website at: www.bpsbioscience.com

- 6) Initiate reaction by adding 2.5 μ l of diluted **BRD3 (BD2)** prepared as described above. Incubate at room temperature for 30 minutes.

Step 2:

Note: Protect your samples from direct exposure to light!


- 1) Dilute GSH Acceptor beads (PerkinElmer #AL109C) 250-fold with 1x BRD Homogeneous Detection Buffer 1. Add 10 μ l per well. Shake plate briefly. Incubate at room temperature for 30 minutes.

Step 3:

- 1) Dilute Streptavidin-conjugated donor beads (PE #6760002S) 250-fold with 1x Homogeneous Detection Buffer 1. Add 10 μ l per well. Incubate at room temperature for 15 – 30 minutes.
- 2) Read Alpha-counts.

Due to lot to lot variability in AlphaScreen® bead performance, it may be necessary to optimize assay conditions. For example, slight adjustments to bromodomain or ligand concentrations may improve signal-to-noise ratio.

Example of Assay Results:

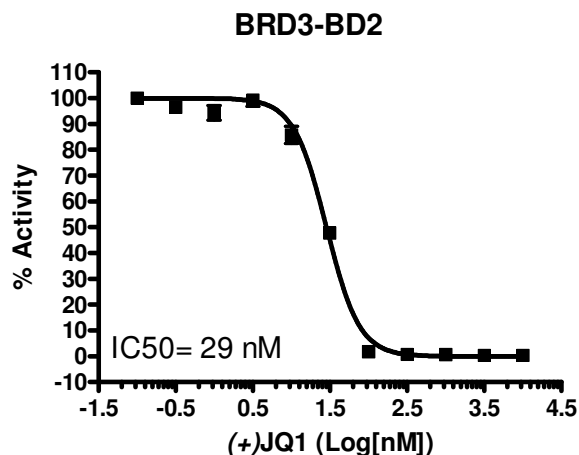

BRD3 (BD2) binding activity, measured using the BRD3 (BD2) Inhibitor Screening Assay Kit, BPS Bioscience, Catalog #32523 and (+)-JQ1 Inhibitor, Catalog #27400. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com.*

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6044 Cornerstone Ct. West, Ste. E
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

RELATED PRODUCTS:

<u>Product Name</u>	<u>Catalog #</u>	<u>Size</u>
(+)-JQ1	27400	10 mg
BRD2 (339 – 459), His-tag	31020	100 µg
BRD2 (65 – 187), GST-tag	31021	100 µg
BRD3 (29 – 145), His-tag	31030	100 µg
BRD3 (29 – 145), GST-tag	31032	100 µg
BRD3 (306 – 417), His-tag	31031	100 µg
BRD3 (306 – 417), GST-tag	31033	100 µg
BRD4 (49 – 170), GST-tag	31040	100 µg
BRD4 (342 – 460), GST-tag	31041	100 µg
BRD4 (49 – 170), His-tag	31042	100 µg
BRD4 (342 – 460), His-tag	31043	100 µg
BRD9 (135 – 242), His-tag	31090	100 µg
BRDT (22 – 138), GST-tag	31108	100 µg
BRDT (22 – 138), His-tag	31101	100 µg
BRDT (257 – 382), His-tag	31100	100 µg
BET Bromodomain Ligand	33000	0.5 mL
BRD3 (BD1) Inhibitor Screening Kit	32513	384 rxns.
BRD4 (BD1) Inhibitor Screening Kit	32514	384 rxns.
BRD4 (BD2) Inhibitor Screening Kit	32524	384 rxns.

AlphaScreen® and AlphaLISA® are registered trademarks of PerkinElmer, Inc.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com