



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6044 Cornerstone Court West, Suite E
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet

Fluorogenic HDAC Substrate (Green fluorescence)

Catalog #: 50038
Lot#: 130903

DESCRIPTION: A fluorogenic, acetylated peptide substrate for HDACs. It is a suitable substrate for many HDACs, except class 2a HDACs.

PURITY: >95% by HPLC.

APPLICATIONS: Study enzyme kinetics, and screen small molecule inhibitors for drug discovery and HTS applications.

SUPPLIED AS: DMSO solution

QUANTITY: 50 μ l @ 5 mM

STORAGE: -20 °C or -70 °C.

REFERENCES:

1. Ito, A., *et al.*, *EMBO J.* 2001;**20**:1331.
2. Barlev, N.A., *et al.*, *Mol Cell.* 2001;**8**:1243.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.
To place your order, please contact us by Phone **1.858.829.3082** Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6044 Cornerstone Court West, Suite E
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

ASSAY PROTOCOL:

Material: Assay buffer (BPS #50031); Assay developer (BPS #50030); HDAC Substrate (BPS #50038)

Step 1: Add all reaction mixture to a low binding NUNC black plate (VWR catalog number 62408-936)

2 μ l of HDAC1 (0.33 μ g/ μ l)
38 μ l of HDAC assay buffer (BPS #50031)
5 μ l of 1 mg/ml BSA
5 μ l of 200 μ M substrate (BPS #50038)

Incubate at 37 °C for 30 min.

Step 2: Stop the reaction

Add 50 μ l of HDAC assay developer (2x) (BPS #50030) and incubate the plate at room temperature for 15 min.

Step 3: Read sample in a microtiter-plate reading fluorimeter capable of excitation at a wavelength in the range of ~485 nm and detection of emitted light in the range of 528 nm. "Blank" value is subtracted from all other values.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.
To place your order, please contact us by Phone **1.858.829.3082** Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com