



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 


Data Sheet

Fluorogenic HDAC6 Assay Kit

Catalog #: 50076

DESCRIPTION: The *Fluorogenic HDAC6 Assay Kit* is a complete assay system designed to measure histone deacetylase 6 (HDAC6) activity for screening and profiling applications. It comes in a convenient 96-well format, with all the reagents necessary for 100 fluorescent HDAC6 activity measurements. In addition, the kit includes purified HDAC6 enzyme and a potent HDAC inhibitor, Trichostatin A for use as a positive and negative control. The *Fluorogenic HDAC6 Assay Kit* is based on a unique fluorogenic substrate and developer combination. This assay method eliminates dealing with the radioactivity, extraction, and chromatography aspects of traditional assays. Using this kit, only two simple steps on a microtiter plate are needed to analyze the HDAC activity level. First, the HDAC fluorometric substrate, containing an acetylated lysine side chain, is incubated with purified HDAC enzyme. The deacetylation sensitizes the substrate so subsequent treatment with the Lysine Developer produces a fluorophore that can then be measured using a fluorescence reader.

HDACs regulate cellular processes by catalyzing the hydrolysis of an acetyl group from acetyllysines in modified proteins. In the HDAC assay, fluorescent-dye molecules are attached to a peptide containing acetyllysine. Attachment to the peptide quenches the fluorescence of the dye. After treatment of the peptide with an HDAC, the reaction is mixed with a development solution that is specific for nonacetylated lysines. If the acetyl group has been removed from the lysine by the HDAC, this solution will release the dye allowing for fluorescence. Fluorescence is therefore directly related to HDAC activity.

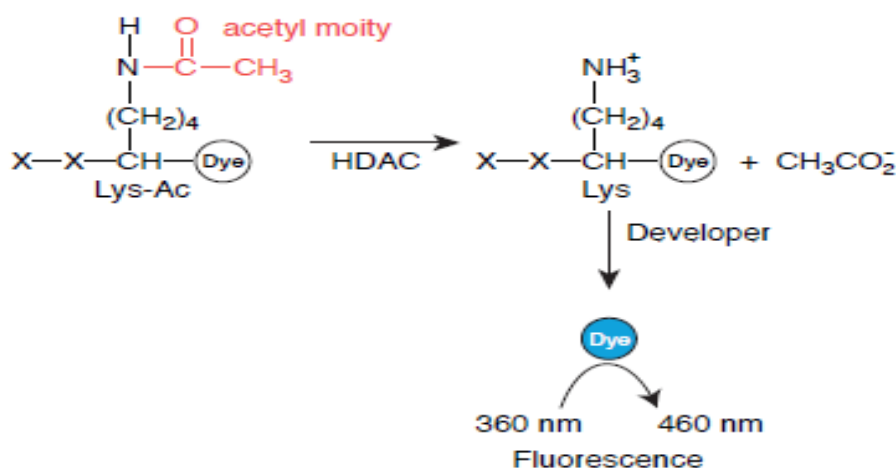

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

COMPONENTS:

Catalog #	Reagent	Amount	Storage	
50006	HDAC6 human recombinant enzyme	9 µg	-80°C	Avoid Freeze/Thaw Cycles!
50037	Fluorogenic HDAC substrate 3 (5 mM)	50 µl	-80°C	
50030	2x HDAC Developer (contains Trichostatin A) (50 µM)	6 ml	-80°C	
	Trichostatin A in DMSO (200 µM)	100 µl	-20°C	
50031	HDAC Assay Buffer	10 ml	-20°C	
79685	Black, low binding NUNC black microtiter plate	1	Room temp.	

APPLICATIONS: Great for studying enzyme kinetics and screening small molecular inhibitors for drug discovery and HTS applications.

MATERIALS REQUIRED BUT NOT SUPPLIED:

0.1% solution (1 mg/ml) of bovine serum albumin (BSA) in water
Fluorimeter capable of excitation at 350-380 nm and detection at 440-460 nm.
Adjustable micropipettor and sterile tips
Rotating or rocker platform

REFERENCE(S):

1. Santo, L., *et al.*, *Blood*. 2012 Mar 15;**119(11)**:2579-89.
2. Bradner, J.E., *et al.*, *Nat Chem Biol*. 2010 Mar;**6(3)**: 238-243.

ASSAY PROTOCOL:

Immediately prior to assay:

- 1) Dilute **Trichostatin A in DMSO (200 µM)** stock 10-fold with **HDAC Assay Buffer** to make a 20 µM solution. (Make only sufficient quantity needed for the assay; store remaining **Trichostatin A in DMSO (200 µM)** stock solution in aliquots at -80°C).
- 2) Dilute **Fluorogenic HDAC substrate 3 (5 mM)** stock 25-fold with **HDAC Assay Buffer** to make a 200 µM solution. (Make only sufficient quantity needed for the assay; store remaining 5 mM stock solution in aliquots at -80°C).
- 3) Dilute **HDAC6 human recombinant enzyme** in **HDAC Assay Buffer** to 7 ng/µl (35 ng/reaction)*. Aliquot any remaining enzyme and store undiluted at -80°C. Keep diluted enzyme on ice. Discard any remaining diluted enzyme after use. *Note: optimal enzyme concentration may vary with the specific activity of the enzyme.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Step 1:

In duplicate, add the reaction mixtures (below) to the microtiter black plate as follows:

- 1) Prepare the master mixture: N wells × (5 μ l **Fluorogenic HDAC substrate 3 (200 μ M)** + 5 μ l BSA (1 mg/ml) + 30 μ l **HDAC Assay Buffer**). Add 40 μ l of master mixture to all wells.
- 2) Add 5 μ l of inhibitor solution of each well designated "Test Inhibitor." For the "Positive Control" and "Blank," add 5 μ l of 10% DMSO in water (inhibitor buffer). Add 5 μ l of diluted **Trichostatin A in DMSO (20 μ M)** to the well designated "Inhibitor Control." Keep final DMSO concentration at or below 1%.
- 3) Add 5 μ l of **HDAC Assay Buffer** to the wells designated "Blank."
- 4) Initiate reaction by adding 5 μ l of diluted **HDAC6 human recombinant enzyme** to the wells designated "Positive Control," "Test Inhibitor," and "Inhibitor Control." Incubate at 37 °C for 30 min.

	"Blank"	Positive Control	Test Inhibitor	Inhibitor Control
HDAC substrate (200 μ M)	5 μ l	5 μ l	5 μ l	5 μ l
BSA (1 mg/ml)	5 μ l	5 μ l	5 μ l	5 μ l
HDAC Assay Buffer	35 μ l	30 μ l	30 μ l	30 μ l
Diluted Trichostatin A (20 μ M)	-	-	-	5 μ l
Test Inhibitor	-	-	5 μ l	-
10% DMSO in water (Inhibitor buffer)	5 μ l	5 μ l	-	-
Diluted HDAC6 (7 ng/ μ l)	-	5 μ l	5 μ l	5 μ l
Total	50 μ l	50 μ l	50 μ l	50 μ l

Step 2:

Add 50 μ l of undiluted **2x HDAC Developer** to each well. Incubate the plate at room temperature for 15 minutes.

Step 3:


Read sample in a microtiter-plate reading fluorimeter capable of excitation at a wavelength in the range of 350-380 nm and detection of emitted light in the range of 440-460 nm. "Blank" value is subtracted from all other values.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Example of Assay Results:

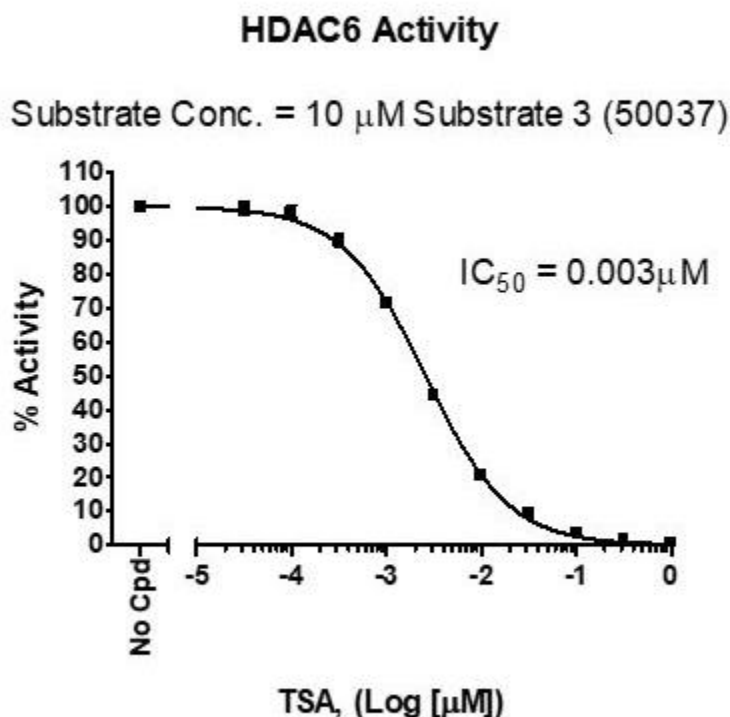

HDAC6 enzyme activity, measured using the *HDAC6 Fluorogenic Assay Kit*, BPS Bioscience Catalog #50076. Fluorescence was measured using a Bio-Tek fluorescent microplate reader. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com.*

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

RELATED PRODUCTS:

<u>Product</u>	<u>Catalog #</u>	<u>Size</u>
HDAC6 (C-Flag)	50056	50 µg
HDAC6 (N-GST)	50006	50 µg
HDAC6 (H216A)	50046	50 µg
HDAC6 (H611A)	50066	50 µg
HDAC1, FLAG-tag, His-tag	50051	50 µg
HDAC2, FLAG-tag	50052	50 µg
HDAC3/NcoR2 Complex, His-tag	50003	50 µg
HDAC4, GST-tag, His-tag	50004	10 µg
HDAC5, Catalytic Domain, His-tag	50005	10 µg
HDAC7, GST-tag	50007	10 µg
HDAC8, His-tag	50008	50 µg
HDAC9, His-tag	50009	10 µg
HDAC10, GST-tag	50010	50 µg
HDAC11, GST-tag	50011	50 µg
HDAC Fluorogenic Assay Kit	50033	96 rxns
HDAC Fluorogenic Assay Kit (Green)	50034	96 rxns
HDAC Class 2a Fluorogenic Assay Kit	50041	96 rxns
HDAC3 Fluorogenic Assay Kit	50073	96 rxns
HDAC8 Assay Kit	50068	96 rxns

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [**info@bpsbioscience.com**](mailto:info@bpsbioscience.com)