



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Ct. West, Ste. B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet
PRMT5 Homogeneous Assay Kit
Catalog #52052
Size: 384 reactions

DESCRIPTION: The *PRMT5 Homogeneous Assay Kit* is designed to measure PRMT5 activity for screening and profiling applications. PRMT5 is a histone methyltransferase that exhibits methylation activity toward H4-R3. The *PRMT5 Homogeneous Assay Kit* comes in a convenient AlphaLISA® format, with biotinylated histone H4 peptide substrate, primary antibody, methylation assay buffer, and purified PRMT5 for 384 enzyme reactions. The key to the *PRMT5 Homogeneous Assay Kit* is a highly specific antibody that recognizes methylated substrate. With this kit, only three simple steps on a microtiter plate are required for methyltransferase detection. First, a sample containing PRMT5 enzyme is incubated with the biotinylated substrate for two hours. Next, acceptor beads and primary antibody are added, then donor beads, followed by reading the Alpha-counts.

COMPONENTS:

Catalog #	Component	Amount	Storage	
51045	PRMT5/MEP50	2x 40 µg	-80°C	<i>Avoid freeze/ thaw cycles!</i>
52120-A	60 µM S-adenosylmethionine	250 µl	-80°C	
52150-3	Primary antibody 4-3 *	100 µl	-80°C	
	Biotinylated histone H4 peptide substrate	500 rxns	-80°C	
	4x PRMT5 assay buffer (add DTT before experiment)	3x1 ml	-20°C	
79389	4x Detection buffer	2 ml	-20°C	

*The detection antibody for this kit was changed in January, 2017.

MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

AlphaLISA® anti-rIgG acceptor beads, 5 mg/ml (PerkinElmer #AL104C)
AlphaScreen® Streptavidin-conjugated donor beads, 5 mg/ml (PerkinElmer #6760002)
Optiplate-384 (PerkinElmer #6007290)
DTT 0.5M

APPLICATIONS: Great for studying enzyme kinetics and HTS applications.

CONTRAINDICATIONS: Green and blue dyes that absorb light in the AlphaScreen® signal emission range (520-620 nm), such as Trypan Blue. Avoid the use of the potent singlet oxygen quenchers such as sodium azide (NaN₃) or metal ions (Fe²⁺, Fe³⁺, Cu²⁺, Zn²⁺ and Ni²⁺). The presence of >1% RPMI 1640 culture medium leads to a signal reduction due to the presence of excess biotin and iron in this medium. MEM, which lacks these components, does not affect AlphaScreen® assays.

STABILITY: At least one year from date of receipt when stored as directed.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Ct. West, Ste. B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

REFERENCE(S): Dillon SC, *et al.* 2005. *Genome Biology* 6:227.

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate.

Step 1:

- 1) Re-suspend tube with **Biotinylated histone H4 peptide substrate** in 500 μ L of distilled water.
- 2) Add 10 μ l **0.5M DTT** (not provided) to 1-ml tube with **4x PRMT5 Assay Buffer**. Prepare **1x PRMT5 Assay Buffer** by adding 1 part of **4x PRMT5 Assay Buffer** to 3 parts water (v/v).
- 3) Prepare the master mixture: N wells x (2 μ l **4x PRMT5 Assay Buffer** + 0.5 μ l **S-adenosylmethionine** (60 μ M) + 1 μ l **Biotinylated substrate** + 1 μ l water). Add 5 μ l to wells designated "Positive Control", "Test Sample", and "Blank". To wells labeled "Substrate Control", add 2 μ l **4x PRMT5 Assay Buffer** + 1 μ l **Biotinylated substrate** + 2 μ l water.
- 4) Add 3 μ l of Inhibitor solution of each well labeled as "Test Inhibitor". For the "Positive Control", "Substrate Control" and "Blank", add 3 μ l of the same solution without inhibitor (Inhibitor buffer).
- 5) Thaw **PRMT5** on ice. Upon first thaw, briefly spin tube containing enzyme to recover full content of the tube. Aliquot **PRMT5** enzyme into single use aliquots. Store remaining undiluted enzyme in aliquots at -80°C. *Note: PRMT5 is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme. Depending on the nature of inhibitor, pre-incubation with the enzyme may be necessary.*
- 6) Dilute **PRMT5** in **1x PRMT5 Assay Buffer** at 100 ng/ μ l. Keep diluted enzyme on ice until use. Discard any unused diluted enzyme after use.

	Positive Control	Test Sample	Substrate Control	Blank
4x PRMT5 Assay Buffer	2 μ l	2 μ l	2 μ l	2 μ l
60 μ M S-adenosylmethionine	0.5 μ l	0.5 μ l	-	0.5 μ l
Biotinylated substrate	1 μ l	1 μ l	1 μ l	1 μ l
H ₂ O	1.5 μ l	1.5 μ l	2 μ l	1.5 μ l
Test Inhibitor/Activator	-	3 μ l	-	-
Inhibitor Buffer (no inhibitor)	3 μ l	-	3 μ l	3 μ l
PRMT5 (100 ng/ μ l)	2 μ l	2 μ l	2 μ l	-
1x PRMT5 Assay Buffer	-	-	-	2
Total	10 μl	10 μl	10 μl	10 μl

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com

- 7) To the wells designated as "Blank", add 2 μ l of 1x PRMT5 Assay buffer.
- 8) Initiate reaction by adding 2 μ l of diluted **PRMT5** enzyme to the wells designated "Positive Control", "Substrate Control", and "Test Sample". Incubate at room temperature for 2 hours.
Protect your samples from direct exposure to light for steps 2 and 3!


Step 2:

- 1) Dilute anti-Rabbit Acceptor beads (PerkinElmer #AL104C) 1:250-fold with **1x Detection Buffer** (made by diluting **4x Detection buffer** 1:4 in distilled water). Add 5 μ l per well. Shake plate briefly.
- 2) Dilute **Primary antibody 4-3** 30-fold with **1x Detection Buffer**. Add 5 μ l per well. Shake plate. Incubate 30-60 min at room temperature.
(Alternatively, dilute anti-Rabbit Acceptor beads (1:500) and Primary antibody 4-3 (1:60) with 1x Detection buffer in one step. Add 10 μ L of acceptor beads/antibody mixture per well.)

Step 3:

- 1) Dilute Streptavidin-conjugated donor beads (PerkinElmer #6760002) 125-fold with **1x Detection Buffer**. Add 10 μ l per well. Incubate for 30 min. at room temperature.
- 2) Read Alpha-counts.

Example of Assay Results:

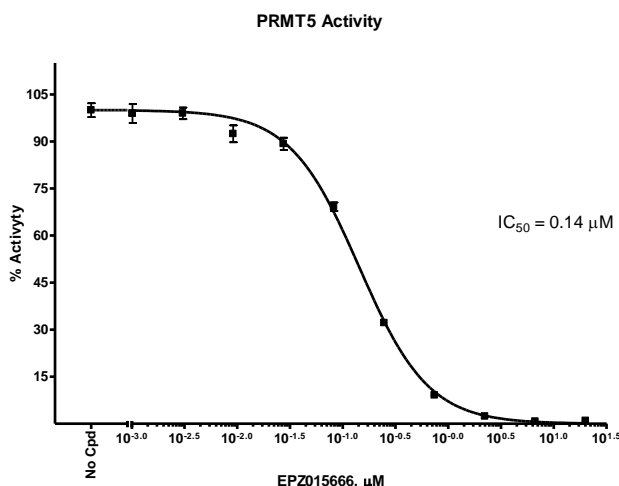

PRMT5 enzyme activity, measured using the *PRMT5 Homogeneous Assay Kit*, BPS Bioscience Cat. #52052. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com*

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Ct. West, Ste. B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

RELATED PRODUCTS:

PRMT1 Homogeneous Assay Kit	#52054	384 reactions
PRMT3 Homogeneous Assay Kit	#52055	384 reactions
PRMT6 Homogeneous Assay Kit	#52056	384 reactions
PRMT8 Homogeneous Assay Kit	#52058	384 reactions
PRMT1 Chemiluminescent Assay Kit	#52004L	96 reactions
PRMT3 Chemiluminescent Assay Kit	#52005L	96 reactions
PRMT4 Chemiluminescent Assay Kit	#52041L	96 reactions
PRMT5 Chemiluminescent Assay Kit	#52002L	96 reactions
PRMT6 Chemiluminescent Assay Kit	#52046	96 reactions
PRMT5 recombinant protein (HEK293)	#51045	20 µg
PRMT5/MEP50 recombinant protein (Sf9)	#51048	20 µg
PRMT1 recombinant protein (<i>E. coli</i>)	#51040	50 µg
PRMT1 recombinant protein (Sf9)	#51041	20 µg
PRMT3 recombinant protein	#51043	50 µg
PRMT4 (CARM 1) recombinant protein	#51047	20 µg
PRMT6 recombinant protein	#51049	20 µg
PRMT7 recombinant protein	#51054	20 µg
PRMT8 recombinant protein	#51052	20 µg
PRMT9 recombinant protein	#51053	20 µg

AlphaScreen® and AlphaLISA® are registered trademarks of PerkinElmer, Inc.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com