



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 


## Data Sheet

### ***PDE3B Assay Kit***

Catalog # 60331  
 Size: 96 reactions

**DESCRIPTION:** Phosphodiesterases (PDEs) play an important role in dynamic regulation of cAMP and cGMP signaling. PDE3B, also known as cGMP-inhibited phosphodiesterase, is involved in mediating the antilipolytic and antiglycogenolytic effects of insulin in adipose and liver tissues. The *PDE3B Assay Kit* is designed for identification of PDE3B inhibitors using fluorescence polarization. The assay is based on the binding of a fluorescent nucleotide monophosphate generated by PDE3B to the binding agent.

Phosphodiesterases catalyze the hydrolysis of the phosphodiester bond in dye-labeled cyclic monophosphates. Beads selectively bind the phosphate group in the nucleotide product. This increases the size of the nucleotide relative to unreacted cyclic monophosphate. In the polarization assay, dye molecules with absorption transition vectors parallel to the linearly-polarized excitation light are selectively excited. Dyes attached to the rapidly-rotating cyclic monophosphates will obtain random orientations and emit light with low polarization. Dyes attached to the slowly-rotating nucleotide-bead complexes will not have time to reorient and therefore will emit highly polarized light.

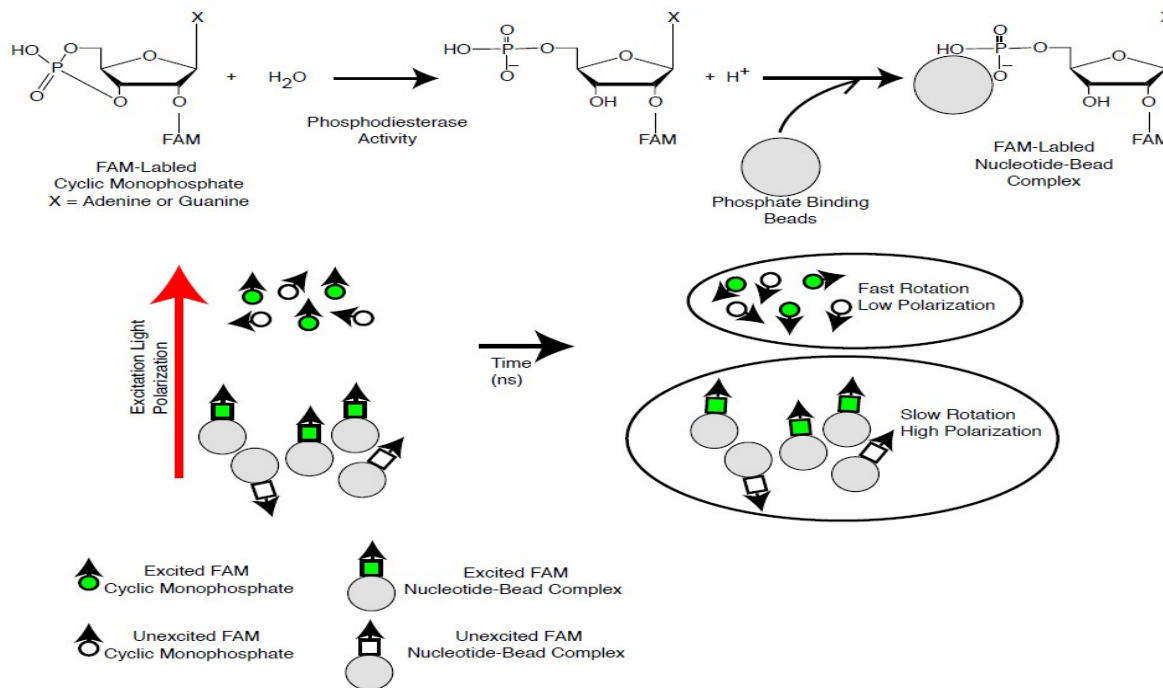

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**  
 Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
 Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

The *PDE3B Assay Kit* comes in a convenient 96-well format, with purified PDE3B enzyme, fluorescently labeled PDE3B substrate (cAMP), binding agent, and PDE assay buffer for 100 enzyme reactions. The key to the *PDE3B Assay Kit* is the specific binding agent. Using this kit, only two simple steps on a microtiter plate are required for PDE3B reactions. First, the fluorescently labeled cAMP is incubated with a sample containing PDE3B for 1 hour. Second, a binding agent is added to the reaction mix to produce a change in fluorescent polarization that can then be measured using a fluorescence reader equipped for the measurement of fluorescence polarization.

#### COMPONENTS:

| Catalog # | Component                            | Amount | Storage    |                                    |
|-----------|--------------------------------------|--------|------------|------------------------------------|
| 60031     | PDE3B recombinant enzyme             | 1 µg   | -80°C      | <b>(Avoid freeze/thaw cycles!)</b> |
| 60200     | FAM-Cyclic-3', 5'-AMP (20 µM)        | 50 µl  | -80°C      |                                    |
| 60393     | PDE assay buffer                     | 25 ml  | -20°C      |                                    |
| 60390     | Binding Agent                        | 100 µl | +4°C       |                                    |
| 60391     | Binding Agent Diluent (cAMP)         | 10 ml  | +4°C       |                                    |
| 79685     | Black, low binding, microtiter plate | 1      | Room temp. |                                    |

#### MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

Fluorescent microplate reader capable of measuring fluorescence polarization

**APPLICATIONS:** Great for studying enzyme kinetics and screening small molecular inhibitors for drug discovery and HTS applications.

**STABILITY:** 6 months from date of receipt when stored as directed.

**REFERENCE:** Chandrasekaran A, *et al.*, *Cell Signal*. 2008; **20(1)**: 139-53.

#### ASSAY PROTOCOL:

**All samples and controls should be tested in duplicate.**

##### Step 1:

- 1) Dilute 20 µM FAM-Cyclic-3', 5'-AMP stock 100-fold with PDE buffer to make a 200 nM solution. Make only sufficient quantity needed for the assay; store remaining 20 µM stock solution in aliquots at -20°C.
- 2) Add 25 µl of FAM-Cyclic-3', 5'-AMP (200 nM) to each well designated "Positive Control", "Test Inhibitor", and "Substrate Control".

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

- 3) Add 20  $\mu$ l of PDE assay buffer to each well designated "Substrate Control" and 45  $\mu$ l of PDE assay buffer to each well designated "Blank".
- 4) Add 5  $\mu$ l of inhibitor solution to each well designated "Test Inhibitor". For the wells labeled "Positive Control", "Substrate Control" and "Blank", add 5  $\mu$ l of the same solution without inhibitor (inhibitor buffer).
- 5) Thaw PDE3B on ice. Upon first thaw, briefly spin tube containing enzyme to recover the full contents of the tube. Aliquot PDE3B enzyme into single-use aliquots. Store remaining undiluted enzyme in aliquots at  $-70^{\circ}\text{C}$  immediately. *Note: PDE3B is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*
- 6) Dilute PDE3B in PDE buffer to 50 pg/ $\mu$ l (1 ng/reaction)\*. Initiate reaction by adding 20  $\mu$ l of diluted PDE3B to the wells designated "Positive Control" and "Test Inhibitor." Discard any remaining diluted enzyme after use. *\*Note: Optimal enzyme concentration may vary with the specific activity of the enzyme.*
- 7) Incubate at room temperature for 1 hour.

|                                 | Positive Control            | Test Inhibitor              | Substrate Control           | "Blank" Negative Control    |
|---------------------------------|-----------------------------|-----------------------------|-----------------------------|-----------------------------|
| FAM-Cyclic-3',5'-AMP (200 nM)   | 25 $\mu$ l                  | 25 $\mu$ l                  | 25 $\mu$ l                  | -                           |
| PDE assay buffer                | -                           | -                           | 20 $\mu$ l                  | 45 $\mu$ l                  |
| Inhibitor (in PDE assay buffer) | -                           | 5 $\mu$ l                   | -                           | -                           |
| Inhibitor Buffer (no inhibitor) | 5 $\mu$ l                   | -                           | 5 $\mu$ l                   | 5 $\mu$ l                   |
| PDE3B (50 pg/ $\mu$ l)          | 20 $\mu$ l                  | 20 $\mu$ l                  | -                           | -                           |
| <b>Total</b>                    | <b>50 <math>\mu</math>l</b> | <b>50 <math>\mu</math>l</b> | <b>50 <math>\mu</math>l</b> | <b>50 <math>\mu</math>l</b> |

## Step 2:

- 1) Mix **binding agent** thoroughly and dilute **binding agent** 1:100 with binding agent diluent.
- 2) Add 100  $\mu$ l diluted binding agent to each microwell. Incubate at room temperature for 1 hour with slow shaking.
- 3) Read the fluorescent polarization of the sample in a microtiter-plate reader equipped for the measurement of fluorescence polarization, capable of excitation at wavelengths ranging from  $485 \pm 5$  nm and detection of emitted light ranging from  $528 \pm 10$  nm. Blank value is subtracted from all other values.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**  
Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

## CALCULATING RESULTS:

### Definition of Fluorescence Polarization

$$P = \frac{I_{\parallel} - I_{\perp}}{I_{\parallel} + I_{\perp}}$$

where  $I_{\parallel}$  = Intensity with polarizers parallel and  $I_{\perp}$  = Intensity with polarizers perpendicular.

The equation above assumes that light is transmitted equally well through both parallel and perpendicular oriented polarizers. In practice, this is generally not true and a correction must be made to measure the absolute polarization state of the molecule. This correction factor is called the "G Factor".

$$\text{FP(measured)} = \frac{(I_{\parallel} - G * I_{\perp})}{(I_{\parallel} + G * I_{\perp})} * 1000$$

The G-factor is instrument-dependent and may vary slightly depending upon instrument and conditions. Please check the manual of your instrument to obtain the information about the establishment of the G-factor.


OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**  
Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

#### EXAMPLE OF ASSAY RESULTS:

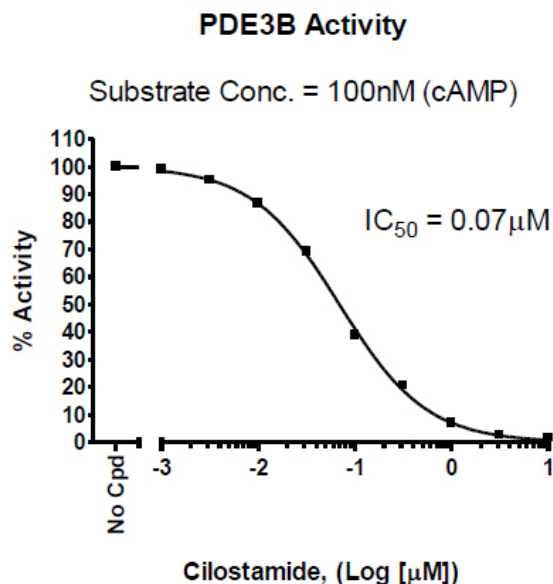

Inhibition of PDE3B activity by cilostamide, measured using the PDE3B Assay Kit, BPS Bioscience # 60331. Fluorescence polarization was measured at 528 nm using a Tecan M1000 fluorescent microplate reader. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at [info@bpsbioscience.com](mailto:info@bpsbioscience.com)*

#### RELATED PRODUCTS :

| <u>Product Name</u> | <u>Catalog #</u> | <u>Size</u> |
|---------------------|------------------|-------------|
| PDE3A               | 60030            | 10 µg       |
| PDE3B               | 60031            | 10 µg       |
| PDE Assay Kit       | 60300            | 96 rxns.    |
| PDE1B Assay Kit     | 60311            | 96 rxns.    |
| PDE3A Assay Kit     | 60330            | 96 rxns.    |
| PDE4A Assay Kit     | 60340            | 96 rxns.    |
| PDE4D Assay Kit     | 60345            | 96 rxns.    |
| PDE10A Assay Kit    | 60400            | 96 rxns.    |
| FAM-cAMP Substrate  | 60200            | 100 nmol    |

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**  
Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)