

# Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten! See the following pages for more information!



# Lieferung & Zahlungsart

siehe unsere Liefer- und Versandbedingungen

# Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

# SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

linkedin.com/company/szaboscandic in





6042 Cornerstone Court W, Ste B San Diego, CA 92121 **Tel:** 1.858.829.3082

Fax: 1.858.481.8694 Email: info@bpsbioscience.com

# Data Sheet Nav1.7 – HEK 293 Cell line Catalog #: 60507

# **Description**

Stable recombinant HEK293 cell line expressing human Nav1.7 (Genbank # Q15858) fused at the C-terminus to Green Fluorescent Protein [Ex. ~395 nm, 475 nm; em ~510 nm] and with C-terminal Streptavidin-Binding Peptide (SBP) fusion. Nav1.7 is a tetrodotoxin-sensitive voltage-gated sodium channel type IX subunit alpha (SCN9A).

## **Background**

Nav1.7 is a voltage-gated sodium ion channel that in humans is encoded by the SCN9A gene. It is usually expressed at high levels in two types of neurons, the nociceptive neurons at dorsal root ganglion and trigeminal ganglion, and sympathetic ganglion neurons, which are part of the autonomic (involuntary) nervous system.

Nav1.7 is present at the endings of pain-sensing nerves, the nociceptors, close to the region where the impulse is initiated. The Nav1.7 channel produces a rapidly activating and inactivating current which is sensitive to the level of tetrodotoxin. Knockout mice that lack Nav1.7 in nociceptors showed reduced response to inflammatory pain [1].

# Sequence

A synthetic codon-optimized DNA sequence encoding human NaV1.7 protein [2] with C-terminal Green Fluorescent Protein (GFP) and C-terminal Streptavidin-Binding Peptide (SBP) [3] tag is stably integrated in tetracycline-inducible HEK293 cells.

## Sequence of the NaV1.7 Tag:

EGGGGSMVSKGEELFTGVVPILVELDGDVNGHKFSVSGEGEGDATYGKLTLKFICTTGKLPVPWPTLVTT LTYGVQCFSRYPDHMKQHDFFKSAMPEGYVQERTIFFKDDGNYKTRAEVKFEGDTLVNRIELKGIDFKED GNILGHKLEYNYNSHNVYIMADKQKNGIKVNFKIRHNIEDGSVQLADHYQQNTPIGDGPVLLPDNHYLST QSALSKDPNEKRDHMVLLEFVTAAGITLGMDELYKVEGLVPRGSGSLVPRGSSAKETAAAKFERQHMDSG ATETSQVAPAGAAAMDEKTTGWRGGHVVEGLAGELEQLRARLEHHPQGQREP

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone 1.858.829.3082 Fax 1.858.481.8694

Or you can Email us at: <a href="mailto:info@bpsbioscience.com">info@bpsbioscience.com</a>

Please visit our website at: <a href="www.bpsbioscience.com">www.bpsbioscience.com</a>



6042 Cornerstone Court W, Ste B San Diego, CA 92121

**Tel:** 1.858.829.3082 **Fax:** 1.858.481.8694

Email: info@bpsbioscience.com

# **Applications**

- Drug compound screening
- Functional assays
- Efficient antigen for mouse immunization

#### **Format**

Each vial contains 2 X 10<sup>6</sup> cells in 1 ml of 10% DMSO

#### Host cell

HEK293 cells, tetracycline-inducible

## **Recommended Storage**

Immediately upon receipt, store in liquid nitrogen.

#### MW

The calculated molecular weight is 242 kDa.

# **Stability**

The cell line has been demonstrated to be stable for at least seven continuous passages. For optimal results, it is recommended to use the cells prior to the 7<sup>th</sup> passage. Upon receipt, amplify the cells in culture and make several frozen aliquots for future use.

## **Propagation Medium and Culture Conditions**

**Thaw Medium 4 (BPS Cat. #60181):** DMEM/F12 (1:1) (Hyclone # SH30271.01) supplemented with 10% FBS (Life Technologies #26140-079), 1% Penicillin/Streptomycin (Hyclone # SV30010)

**Growth Medium 4A (BPS Cat. #79535):** Thaw Medium 4 (BPS Cat. #60181) plus 10  $\mu$ g/ml Blasticidin (Life Technologies # R210-01), and 200  $\mu$ g /ml Zeocin (invivogen # ant-zn-1p).

Cells should be grown at  $37^{\circ}$  with 5% CO<sub>2</sub> using Growth Medium 4A (Thaw Medium 4, Blasticidin, and Zeocin).

**To thaw the cells**, it is recommended to quickly thaw the frozen cells from liquid nitrogen in a 37°C water-bath, transfer to a tube containing 10 ml of Thaw Medium 4 (no Blasticidin and Zeocin), spin down the cells, and resuspend cells in pre-warmed Thaw Medium 4 (no Blasticidin and Zeocin). Transfer resuspended cells to a T25 flask and culture in 37°C CO<sub>2</sub>



6042 Cornerstone Court W, Ste B San Diego, CA 92121

Tel: 1.858.829.3082 Fax: 1.858.481.8694 Email: info@bpsbioscience.com


incubator. At first passage switch to Growth Medium 4A (contains Blasticidin and Zeocin). Cells should be split before they reach complete confluence.

**To passage the cells**, rinse cells with phosphate buffered saline (PBS), detach cells from culture vessel with 0.05% Trypsin/EDTA, add Growth Medium 4A and transfer to a centrifuge tube. Spin down cells, resuspend cells in Growth Medium 4A and seed appropriate aliquots of cell suspension into new culture vessels.

# Induction of the target protein expression

To express Nav1.7, cells are induced with DMEM/F12 50/50 supplemented with 10% FBS, 1% Penicillin Streptomycin, 1  $\mu$ g/ml Doxycycline (Biochemika #44577) and 3 mM Na butyrate (Acros Organics #263190250) for 24 hours prior to cell harvesting or assay.

**Figure 1. Western Blot of the NaV1.7 expressing cells.** Western Blot of HEK293 NaV1.7 cells (lanes 1, 2) stained with anti-Sodium channel Nav1.7, clone N68/6 (Millipore, Cat.No. MABN41) with follow staining with Alkaline Phosphatase conjugated Anti-mouse IgG (Rockland Immunochemicals, Cat.No. 610-1502). M: molecular weight marker.

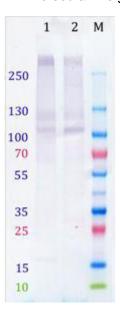





6042 Cornerstone Court W, Ste B San Diego, CA 92121

**Tel:** 1.858.829.3082 **Fax:** 1.858.481.8694

Email: info@bpsbioscience.com

Figure 2. Cell Expression Profile

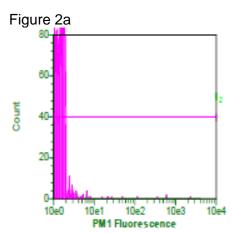

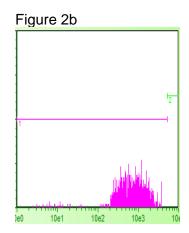

NaV1.7 expression measured by flow cytometry (FACS) using intracellular staining by saponin-permeabilized cells and PE-labeled streptavidin.

Figure 2a: HEK293 Host cells, Figure 2b: NaV1.7 expressing HEK293 cells (#60507) stained with anti-NaV 1.7 clone 68/6 monoclonal antibody and PE-labeled anti-mouse.

#### References

- 1. Catterall, W.A. Cell and Developmental Biology 16: 521–555 (2000).
- 2. Choi, J.S., et al. Neurology 67:1563-1567 (2006).
- 3. Li, Y., el al. Protein Science 20 (1): 140-149 (2011).

## **License Disclosure**

Purchase of this cell line grants you with a 10-year license to use this cell line in your immediate laboratory, for research use only. This license does not permit you to share, distribute, sell, sublicense, or otherwise make the cell line available for use to other laboratories, departments, research institutions, hospitals, universities, or biotech companies. The license does not permit use of this cell line in humans or for therapeutic or drug use. The license does not permit modification of the cell line in any way. Inappropriate use or distribution of this cell line will result in revocation of the license and result in an immediate cease of sales and distribution of BPS products to your laboratory. BPS does not warrant the suitability of the cell line for any particular use, and does not accept any liability in connection with the handling or use of the cell line. Modifications of this cell line, transfer to another facility, or commercial use of the cells may require a separate license and additional fees; contact sales@bpsbioscience.com for details. Publications using this cell line should reference BPS Bioscience, Inc., San Diego.