



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: info@bpsbioscience.com

**Data Sheet**  
**NAMPT Inhibitor Screening Assay Kit**  
Catalog: 71276  
Size: 96 reactions

**BACKGROUND:** NAMPT catalyzes the formation of nicotinamide mononucleotide from nicotinamide and 5-phosphoribosyl-1-pyrophosphate (PRPP). It is the rate limiting component in the mammalian NAD biosynthesis pathway. The protein is thought to be involved in many important biological processes, including metabolism, stress response and aging.

**DESCRIPTION:** The *NAMPT Inhibitor Screening Assay Kit* is designed to measure NAMPT activity for screening and profiling applications. The NAMPT assay kit comes in a convenient 96-well format, with purified recombinant NAMPT enzyme, NAMPT assay buffer, NAMPT dilution buffer, ATP, Nicotinamide, PRPP, and Ethanol sufficient for 96 enzyme reactions.

**COMPONENTS:**

Catalog #	Reagent	Amount	Storage	
91004	NAMPT	50 µg	-80°C	<b>Avoid multiple freeze/thaw cycles!</b>
	NAMPT dilution buffer	3 ml	-80°C	
	4x NAMPT assay buffer	2 x 750 µl	-80°C	
79686	ATP (400 µM)	250 µl	-80°C	
	Nicotinamide (400 µM)	250 µl	-80°C	
	PRPP (800 µM)	250 µl	-80°C	
	Ethanol (30%)	1 ml	-80°C	
79685	Low binding black microtiter 96-well plate	1	Room Temp.	

**MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:**

Fluorescent microplate reader  
Adjustable micropipettor and sterile tips  
Aluminum foil

**APPLICATIONS:** Great for studying enzyme kinetics and screening small molecular inhibitors for drug discovery and HTS applications.

**STABILITY:** Up to 6 months from date of receipt, when stored as recommended.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
**Tel:** 1.858.202.1401  
**Fax:** 1.858.481.8694  
**Email:** [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

**REFERENCE:** Ramsey, K.M., et al. (2009) *Science* **324**: 651-654  
Nakahata, Y., et al. (2009) *Science* **324**: 654-657.

**ASSAY PROTOCOL:**

***All samples and controls should be tested in duplicate.***

1. Thaw **4x NAMPT assay buffer**, **ATP**, **Nicotinamide**, **Phosphoribosyl pyrophosphate (PRPP)**, and **30% Ethanol** on ice.
2. Thaw **NAMPT dilution buffer** and **NAMPT** enzyme on ice. Upon first thaw, briefly spin tube containing enzyme to recover full contents of the tube. Calculate the amount of **NAMPT** required for the assay and dilute enzyme to 20-50 ng/ $\mu$ l with **NAMPT dilution buffer**. Aliquot remaining **NAMPT** enzyme into single-use aliquots. Store remaining undiluted enzyme in aliquots at  $-80^{\circ}\text{C}$ . *NAMPT enzyme is sensitive to freeze/thaw cycles. Avoid multiple freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*
3. Add 10  $\mu$ l of diluted **NAMPT** enzyme to the wells designated "Positive Control" and "Test Inhibitor". To the wells designated as "Blank," add 10  $\mu$ l of **NAMPT dilution buffer**.
4. Add 5  $\mu$ l of Inhibitor solution of each well labeled as "Test Inhibitor." For the wells labeled "Positive Control" and "Blank," add 5  $\mu$ l of the same solution without inhibitor (Inhibitor buffer). *Note: Final DMSO concentration must be  $\leq 1\%$ . Higher DMSO levels can significantly decrease the enzyme activity. For example, to test an inhibitor dissolved in 100% DMSO at 10  $\mu$ M, dilute 1 mM inhibitor with water to make a 100  $\mu$ M inhibitor in 10% DMSO (aq). Then, add 5  $\mu$ l of the 100  $\mu$ M solution into the 50  $\mu$ l assay to make a 1% DMSO concentration in the final reaction mixture.*
5. Preincubate **NAMPT** with the inhibitors for 30 min at room temperature (slow shaking).
6. During enzyme preincubation with the inhibitors, prepare the **Master Mixture** (35  $\mu$ l per well): N wells x (12.5  $\mu$ l **4x NAMPT assay buffer** + 2.5  $\mu$ l **ATP** (400  $\mu$ M) + 2.5  $\mu$ l **Nicotinamide** (400  $\mu$ M) + 2.5  $\mu$ l **PRPP** (800  $\mu$ M) + 2.5  $\mu$ l 30% **Ethanol** + 12.5  $\mu$ l **water**). Add 35  $\mu$ l to every well. Store remaining components at  $-80^{\circ}\text{C}$  in single use aliquots.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

	Positive Control	Test Inhibitor	Blank
4x NAMPT assay buffer	12.5 $\mu$ l	12.5 $\mu$ l	12.5 $\mu$ l
ATP (400 $\mu$ M)	2.5 $\mu$ l	2.5 $\mu$ l	2.5 $\mu$ l
Nicotinamide (400 $\mu$ M)	2.5 $\mu$ l	2.5 $\mu$ l	2.5 $\mu$ l
PRPP (800 $\mu$ M)	2.5 $\mu$ l	2.5 $\mu$ l	2.5 $\mu$ l
30% Ethanol	2.5 $\mu$ l	2.5 $\mu$ l	2.5 $\mu$ l
Water	12.5 $\mu$ l	12.5 $\mu$ l	12.5 $\mu$ l
Test Inhibitor	-	5 $\mu$ l	-
NAMPT dilution buffer in 10% DMSO (Inhibitor buffer)	5 $\mu$ l	-	5 $\mu$ l
NAMPT dilution buffer	-	-	10 $\mu$ l
NAMPT (20-50 ng/ $\mu$ l)	10 $\mu$ l	10 $\mu$ l	-
<b>Total</b>	<b>50 <math>\mu</math>l</b>	<b>50 <math>\mu</math>l</b>	<b>50 <math>\mu</math>l</b>

7. Initiate the reaction by adding 35  $\mu$ l of **Master Mixture** to all the wells. Incubate at 30°C for two hours.
8. After the reaction, read sample in a microtiter-plate reading fluorimeter capable of excitation at a wavelength of ~340 nm and detection of emitted light of ~460 nm. "Blank" value is subtracted from all other values.


OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)

**Example of Assay Results:**

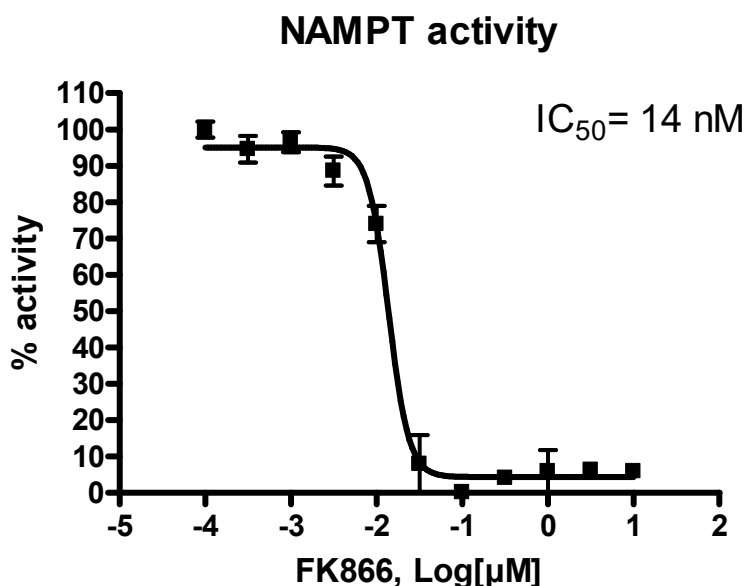

NAMPT inhibition by FK866, measured using the NAMPT Inhibitor Screening Assay Kit, BPS Bioscience Cat.71276. The compound was pre-incubated with the NAMPT enzyme for 30 min at room temperature before the reaction was initiated with the addition of master mix. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at [info@bpsbioscience.com](mailto:info@bpsbioscience.com)*

**RELATED PRODUCTS:**

<u>Product Name</u>	<u>Catalog #</u>	<u>Size</u>
NAMPT (PBEF1), GST tag	91004	50 $\mu$ g
NAMPT (PBEF1)	71098	50 $\mu$ g
NMNAT, His-tag	71090	100 $\mu$ g
CD38, His-Tag (Human), HiP™	71227	100 $\mu$ g
CD38 Inhibitor Screening Assay Kit	71275	96 rxns
5'-Nucleotidase/CD73, His-tag	71184	50 $\mu$ g
CD73 Inhibitor Screening Assay Kit	72055	96 rxns
CD73 Inhibitor Screening Assay Kit	72058	384 rxns
Adenosine Deaminase (ADA), His-tag	70016	100 $\mu$ g
NAD <sup>+</sup> , Biotin-Labeled	80610	500 $\mu$ l

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)