



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: info@bpsbioscience.com

## Data Sheet

### ***PD-L1:B7-1[Biotinylated] Inhibitor Screening Assay Kit***

**Catalog # 72026**  
**Size: 96 reactions**

**BACKGROUND:** There have been a number of studies defining a role for PD-L1 in the regulation of immune tolerance and cancer immunity through binding to T cell-expressed PD-1. Interestingly, PD-L1 has also been shown to interact with B7-1, also known as CD80. Like PD-1:PD-L1 interaction, B7-1:PD-L1 signaling also leads to inhibition of T cell activation, proliferation, and cytokine production, suggesting that this pathway may also serve as a target in cancer immunotherapy and inflammatory diseases.

**DESCRIPTION:** The *PD-L1:B7-1[Biotinylated] Inhibitor Screening Assay Kit* is designed for screening and profiling inhibitors of the PD-L1:B7-1 interaction. This kit comes in a convenient 96-well format, with biotin-labeled B7-1 (CD80), purified PD-L1, streptavidin-labeled HRP, and assay buffer for 100 binding reactions. The key to this kit is the high sensitivity of detection of biotin-labeled B7-1 by streptavidin-HRP. Only a few simple steps on a microtiter plate are required for the assay. First, PD-L1 is coated on a 96-well plate. Next, B7-1-biotin is incubated with PD-L1 on the plate. Finally, the plate is treated with streptavidin-HRP followed by addition of an HRP substrate to produce chemiluminescence, which can then be measured using a chemiluminescence reader.

#### **COMPONENTS:**

Catalog #	Component	Amount	Storage	
71104	PD-L1, Fc fusion,	25 µg	-80°C	<b>(Avoid freeze/thaw cycles!)</b>
71114	B7-1 (CD80), Fc fusion, Biotin-labeled	15 µg	-80°C	
	Streptavidin-HRP	15 µl	+4°C	
79311	3x Immuno Buffer 1	50 ml	-20°C	
	Blocking Buffer	50 ml	+4°C	
	HRP chemiluminescent substrate A (transparent bottle)	6 ml	+4°C	
	HRP chemiluminescent substrate B (brown bottle)	6 ml	+4°C	
	White 96-well microplate	1	+4°C	

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
**Tel:** 1.858.202.1401  
**Fax:** 1.858.481.8694  
**Email:** info@bpsbioscience.com

**MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:**

PBS (Phosphate buffered saline)  
Luminometer or fluorescent microplate reader capable of reading chemiluminescence  
Rotating or rocker platform

**APPLICATIONS:** This kit is useful for screening for inhibitors of PD-L1 binding to B7-1.

**STABILITY:** One year from date of receipt when stored as directed.

**REFERENCES:** 1. Butte, M.J., *et al.*, *Immunity*. 2007; **27(1)**: 111-122.  
2. Butte, M.J. *et al.*, *Mol. Immunol.* 2008; **45**: 3567-3572.

**ASSAY PROTOCOL:**

All samples and controls should be tested in duplicate.

**Coating the plate with PD-L1:**

- 1) Thaw **PD-L1** on ice. Upon first thaw, briefly spin tube containing **PD-L1** to recover the full contents of the tube. Aliquot into single use aliquots. Immediately store remaining PD-L1 in aliquots at -80°C. *Note: PD-L1 is very sensitive to freeze/thaw cycles. Avoid multiple freeze/thaw cycles.*
- 2) Dilute **PD-L1** to 5 µg/ml in PBS.
- 3) Add 50 µl of diluted **PD-L1** solution to each well and incubate overnight at 4°C. Leave a couple of wells empty (uncoated), for use with the "Ligand Control" (see below).
- 4) Dilute **3x Immuno Buffer 1** to **1x Immuno Buffer 1** in water.
- 5) Decant to remove supernatant. Wash the plate 3 times with 100 µl **1x Immuno Buffer 1**. Tap plate onto clean paper towels to remove liquid.
- 6) Block wells by adding 100 µl of **Blocking Buffer** to each well. Incubate for 1 hour at room temperature. Remove supernatant as described in step 4.

**Step 1:**

- 1) Prepare the master mixture: N wells × (10 µl **3x Immuno Buffer 1** + 15 µl H<sub>2</sub>O).
- 2) Add 25 µl of master mixture to each well. Use uncoated wells for the "Ligand Control".

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: info@bpsbioscience.com

- 3) Add 5  $\mu$ l of inhibitor solution to each well designated "Test Inhibitor". For the "Positive Control", "Substrate Control" and "Blank", add 5  $\mu$ l of the same solution without inhibitor (inhibitor buffer). Incubate at room temperature for one hour.

	Blank	Ligand Control	Positive Control	Test Inhibitor
3x Immuno Buffer 1	10 $\mu$ l	10 $\mu$ l	10 $\mu$ l	10 $\mu$ l
H <sub>2</sub> O	15 $\mu$ l	15 $\mu$ l	15 $\mu$ l	15 $\mu$ l
Test Inhibitor/Activator	-	-	-	5 $\mu$ l
Inhibitor buffer (no inhibitor)	5 $\mu$ l	5 $\mu$ l	5 $\mu$ l	-
1x Immuno Buffer 1	20 $\mu$ l	-	-	-
B7-1-biotin (7.5 $\mu$ g/ml)	-	20 $\mu$ l	20 $\mu$ l	20 $\mu$ l
<b>Total</b>	<b>50 <math>\mu</math>l</b>	<b>50 <math>\mu</math>l</b>	<b>50 <math>\mu</math>l</b>	<b>50 <math>\mu</math>l</b>

- 4) Thaw **B7-1-biotin** on ice. Upon first thaw, briefly spin tube containing protein to recover full contents of the tube. Aliquot **B7-1-biotin** into single use aliquots. Immediately store remaining undiluted protein in aliquots at -80°C. *Note: B7-1-biotin is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted protein.*
- 5) Dilute **B7-1-biotin** in **1x Immuno Buffer 1** at 7.5  $\mu$ g/ml. Keep diluted protein on ice until use. Discard any unused diluted protein after use.
- 6) Add 20  $\mu$ l of **1x Immuno Buffer 1** to the well designated "Blank".
- 7) Initiate reaction by adding 20  $\mu$ l of diluted **B7-1-biotin** (see Step 1-5) to wells labeled "Positive Control", "Ligand Control" and "Test Inhibitor". Incubate at room temperature for two hours.
- 8) Decant to remove supernatant. Wash the plate 3 times with 100  $\mu$ l/well **1x Immuno Buffer 1**. Tap plate onto clean paper towels to remove liquid.
- 9) Block wells by adding 100  $\mu$ l of **Blocking Buffer** to each well. Incubate for 10 minutes at room temperature. Tap plate onto clean paper towels to remove liquid.

## Step 2:

- 1) Dilute **Streptavidin-HRP** 1000-fold with **Blocking Buffer**.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
**Tel:** 1.858.202.1401  
**Fax:** 1.858.481.8694  
**Email:** [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

- 2) Add 100  $\mu$ l to each well. Incubate for 1 hour at room temperature with slow shaking.
- 3) Wash plate three times with **1x Immuno Buffer 1**. Tap onto clean paper towels to remove liquid.
- 4) Block wells by adding 100  $\mu$ l of **Blocking Buffer** to each well. Incubate for 10 minutes at room temperature. Decant to remove supernatant. Tap plate onto clean paper towels to remove liquid.
- 5) Just before use, mix on ice 50  $\mu$ l **HRP Chemiluminescent Substrate A** and 50  $\mu$ l **HRP Chemiluminescent Substrate B** per well of the reaction, then add 100  $\mu$ l to each well. Discard any unused chemiluminescent reagent after use.
- 6) Immediately read sample in a luminometer or microtiter-plate capable of reading chemiluminescence. "Blank" value is subtracted from all readings.

#### **Reading Chemiluminescence:**

Chemiluminescence is the emission of light (luminescence) which results from a chemical reaction. The detection of chemiluminescence requires no wavelength selection because the method used is emission photometry and is not emission spectrophotometry.

To properly read chemiluminescence, make sure the plate reader is set for LUMINESCENCE mode. Typical integration time is 1 second, delay after plate movement is 100 msec. Do not use a filter when measuring light emission. Typical settings for the Synergy 2 BioTek plate reader are: use the "hole" position on the filter wheel; Optics position: Top; Read type: endpoint. Sensitivity may be adjusted based on the luminescence of a control assay without binding partner (typically we set this value as 100).

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**



Or you can Email us at: [\*\*info@bpsbioscience.com\*\*](mailto:info@bpsbioscience.com)

Please visit our website at: [\*\*www.bpsbioscience.com\*\*](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
 San Diego, CA 92121  
 Tel: 1.858.202.1401  
 Fax: 1.858.481.8694  
 Email: info@bpsbioscience.com

**Example of Assay Results:**

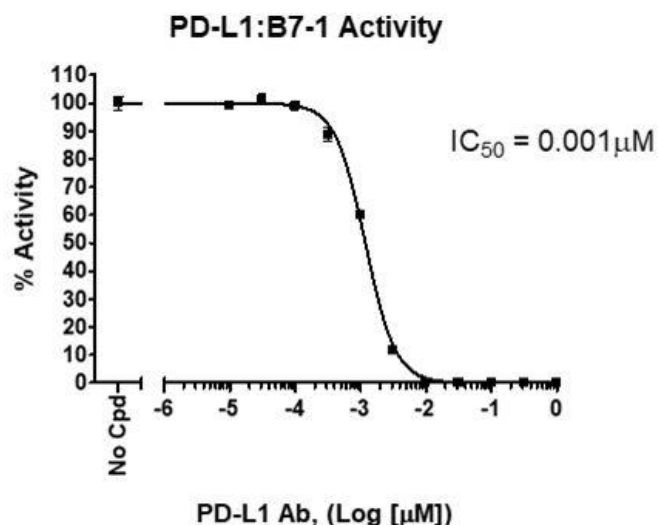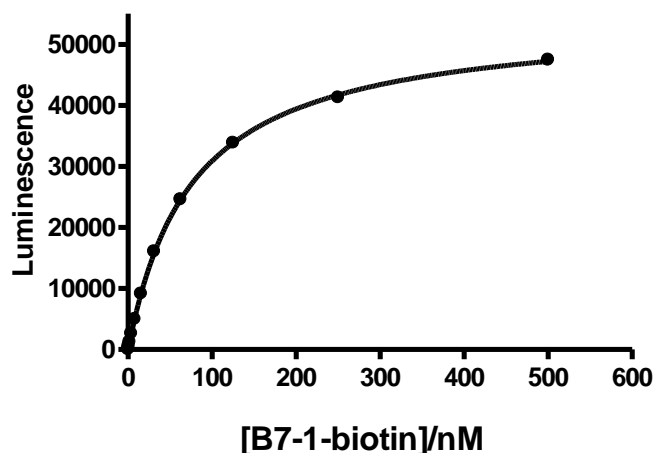

*PD-L1:B7-1-[Biotinylated]* binding (left) and inhibition (right), measured using the *PD-L1:B7-1-[Biotinylated]* Inhibitor Screening Assay Kit, BPS Bioscience, Catalog #72026. Luminescence was measured using a Bio-Tek fluorescent microplate reader. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at [info@bpsbioscience.com](mailto:info@bpsbioscience.com).*

**RELATED PRODUCTS:**

<u>Product Name</u>	<u>Catalog #</u>	<u>Size</u>
B7-1	71125	100 µg
B7-1, Biotin labeled	71114	50 µg
B7-2	71150	100 µg
CD28	71113	200 µg
PD-1	71106	100 µg
PD-1, Biotin labeled	71109	50 µg
PD-L1	71104	100 µg
PD-L1, Biotin-labeled	71105	50 µg
PD-L2	71107	100 µg
PD-L2, Biotin-labeled	71108	50 µg
CTLA4:B7-1[Biotinylated] Inhibitor Screening Assay Kit	72009	96 rxns
CD28:B7-1[Biotinylated] Inhibitor Screening Assay Kit	72007	96 rxns
PD-1:PD-L1[Biotinylated] Inhibitor Screening Assay Kit	72003	96 rxns
PD-1:PD-L2[Biotinylated] Inhibitor Screening Assay Kit	72004	96 rxns

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
 San Diego, CA 92121  
**Tel:** 1.858.202.1401  
**Fax:** 1.858.481.8694  
**Email:** info@bpsbioscience.com

### TROUBLESHOOTING GUIDE

Problem	Possible Cause	Solution
Luminescence signal of positive control reaction is weak	PD-L1 or B7-1 has lost binding capacity	Protein loses binding capacity upon repeated freeze/thaw cycles. Use fresh B7-1-biotin, (BPS Bioscience #71114) and fresh PD-L1 (BPS Bioscience #71104). Store proteins in single-use aliquots. Increase time of protein incubation. Increase protein concentration.
	Incorrect settings on instruments	Refer to instrument instructions for settings to increase sensitivity of light detection.
	Chemiluminescent reagents mixed too soon	Chemiluminescent solution should be used within 15 minutes of mixing. Ensure both reagents are properly mixed.
Luminescent signal is erratic or varies widely among wells	Inaccurate pipetting/technique	Run duplicates of all reactions. Use a multichannel pipettor. Use master mixes to minimize errors.
	Bubbles in wells	Pipette slowly to avoid bubble formation. Tap plate lightly to disperse bubbles; be careful not to splash between wells.
Background (signal to noise ratio) is high	Insufficient washes	Increase number of washes. Increase wash volume. Increase Tween-20 concentration to 0.1% in PBST.
	Sample solvent is inhibiting the protein	Run negative control assay including solvent. Maintain DMSO level at <1% Increase time of protein incubation.
	Results are outside the linear range of the assay	Use different concentrations of B7-1-biotin (BPS Bioscience #71114) to create a standard curve.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)