



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Ct. West, Ste. B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet
CD226:CD155 Homogeneous Assay Kit
Catalog: #72052
Size: 384 reactions

BACKGROUND: CD226 is an activating receptor expressed on the surface of natural killer (NK) cells, monocytes, platelets, and T cells. Its natural ligands are CD112 and CD155 which are expressed by antigen presenting cells (APC). Binding of CD226 to either CD112 or CD155 triggers NK cell effector function and can promote Th1 cell differentiation in T cells. CD226 competes for binding with CD112 and CD155 with the co-inhibitory receptor TIGIT. Agonistic CD226 monoclonal antibodies are being explored as potential therapeutics in settings where immune activation is beneficial, such as cancer.

DESCRIPTION: The *CD226:CD155 Homogeneous Assay Kit* is designed to measure the inhibition of CD226 binding to CD155 (PVR). The *CD226:CD155 Homogeneous Assay Kit* comes in a convenient AlphaLISA[®] format with purified biotinylated CD226, His-tagged CD155, and assay buffer to perform a total of 384 reactions. With this kit, only three simple steps on a microtiter plate are required. First, a sample containing CD226 and an inhibitor of choice is incubated with the CD155 for 60 minutes. Next, acceptor beads are added, then donor beads, followed by reading the Alpha-counts.

COMPONENTS:

| Catalog # | Component | Amount | Storage | |
|-----------|--------------------|---------|---------|---|
| 71253 | CD226-Fc-biotin | 2x 4 µg | -80°C | (Avoid freeze/ thaw cycles!) |
| 71181 | CD155-His | 2x 5 µg | -80°C | |
| 79311 | 3x Immuno Buffer 1 | 4 ml | -20°C | |

MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

AlphaLISA Ni Chelate Acceptor beads, 5 mg/ml (PerkinElmer #AL108C)
AlphaScreen Streptavidin-conjugated Donor beads, 5 mg/ml (PerkinElmer #6760002S)
Optiplate-384 (PerkinElmer #6007290)
AlphaScreen microplate reader
Adjustable micropipettor and sterile tips

APPLICATIONS: Useful for screening for inhibitors of CD226 binding to CD155

CONTRAINDICATIONS: Only limited amounts of DMSO can be included, as it has been shown to disrupt CD226:CD155 interaction. Avoid green and blue dyes that absorb light in the AlphaScreen signal emission range (520-620 nm), such as Trypan Blue. Avoid the use of the potent singlet oxygen quenchers such as sodium azide (NaN₃) or metal ions (Fe²⁺, Fe³⁺, Cu²⁺, Zn²⁺ and Ni²⁺). The presence of >1% RPMI 1640 culture medium leads to a signal reduction due

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Ct. West, Ste. B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

to the presence of excess biotin and iron in this medium. MEM, which lacks these components, does not affect AlphaScreen assays.

STABILITY: At least one year from date of receipt when stored as directed.

REFERENCES: 1. Bottino, C. *et al.*, *J. Exp. Med.* 2003; **198(4)**: 557-567.
2. Pende, D. *et al.*, *Blood.* 2005; **105(5)**: 2066-2073.

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate. Use slow shaking for all incubations.

Step 1:

- 1) Thaw **CD155-His** on ice. Upon first thaw, briefly spin tube containing protein to recover full contents of the tube. Aliquot the protein into single use aliquots. Store remaining undiluted protein in aliquots at -80°C immediately. *Note: CD155-His is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted protein.*
- 2) Dilute one part **3x Immuno Buffer 1** with 2 parts of distilled water (3-fold dilution) to make **1x Immuno Buffer 1**. Make only a sufficient quantity needed for the assay; store remaining stock solution in aliquots at -20°C.
- 3) Dilute **CD155-His** in **1x Immuno Buffer 1** to 4 ng/μl. Keep diluted protein on ice until ready to use. Discard any remaining unused diluted protein after use.
- 4) Prepare the master mixture: N wells × (2 μl **3x Immuno Buffer 1** + 2 μl diluted **CD155-His** + 2 μl distilled water). Add 6 μl of master mixture to every well.

| | Blank | Positive Control | Test Inhibitor |
|---------------------------------|--------------|------------------|----------------|
| 3x Immuno Buffer 1 | 2 μl | 2 μl | 2 μl |
| CD155-His (4 ng/μl) | 2 μl | 2 μl | 2 μl |
| Distilled water | 2 μl | 2 μl | 2 μl |
| Test Inhibitor | - | - | 2 μl |
| Inhibitor buffer (no inhibitor) | 2 μl | 2 μl | - |
| 1x Immuno Buffer 1 | 2 μl | | |
| CD226-biotin (3 ng/μl) | - | 2 μl | 2 μl |
| Total | 10 μl | 10 μl | 10 μl |

- 5) Add 2 μl of inhibitor solution to each well designated “Test Inhibitor”. For the “Positive Control” and “Blank”, add 2 μl of the same solution without inhibitor (inhibitor buffer). *Note: If possible, keep final DMSO concentration below 0.5%.*

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Ct. West, Ste. B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

- 6) Add 2 μ l of **1x Immuno Buffer 1** to the well designated "Blank".
- 7) Thaw **CD226-biotin** on ice. Upon first thaw, briefly spin tube containing protein to recover full contents of the tube. Aliquot the protein into single use aliquots. Store remaining undiluted protein in aliquots at -80°C immediately. *Note: **CD226-biotin** is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted protein.*
- 8) Dilute **CD226-biotin** in **1x Immuno Buffer 1** to 3 ng/ μ l. Keep diluted proteins on ice until use. Discard any remaining unused diluted protein after use.
- 9) Initiate reaction by adding 2 μ l of diluted **CD226-biotin** prepared as described above to each well designated "Positive Control" and "Test Inhibitor". Incubate at room temperature for 60 minutes.

Step 2:

Note: Protect your samples from direct exposure to light!

- 1) Dilute Ni Chelate Acceptor beads (PerkinElmer #AL108C) 250-fold with **1x Immuno Buffer 1**. Add 10 μ l per well. Shake plate briefly. Incubate at room temperature for 30 minutes.

Step 3:

Note: Protect your samples from direct exposure to light!


- 1) Dilute Streptavidin-conjugated donor beads (PE #6760002S) 125-fold with **1x Immuno Buffer 1**. Add 10 μ l per well. Incubate at room temperature for 30 minutes.
- 2) Read Alpha-counts.

Due to lot to lot variability in AlphaScreen[®] bead performance, it may be necessary to optimize assay conditions. For example, slight adjustments to CD226-biotin or CD155-His concentrations may improve signal-to-noise ratio.

Example of Assay Results:

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com

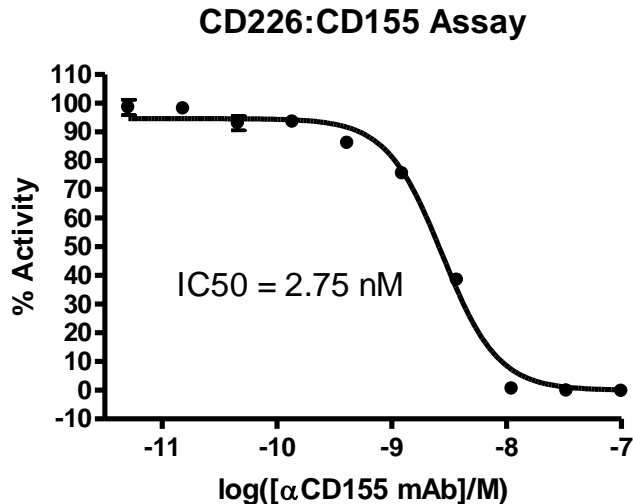

CD226:CD155 inhibition, measured using the CD226:CD155 Inhibitor Screening Assay Kit, BPS Bioscience, Catalog #72052 and a CD155 antibody. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com.*

RELATED PRODUCTS:

| <u>Product Name</u> | <u>Catalog #</u> | <u>Size</u> |
|--|------------------|-------------|
| Human TIGIT, Fc fusion | #71186 | 100 µg |
| Human CD112, His-tag | #71197 | 100 µg |
| Human CD112, His-tag, Biotin-labeled | #71234 | 50 µg |
| Human CD155 (PVR), His-tag | #71181 | 100 µg |
| Mouse CD155 (PVR), His-tag | #71167 | 100 µg |
| Mouse CD155 (PVR), His-tag, Biotin-labeled | #71168 | 50 µg |
| TIGIT:CD112 Homogeneous Assay Kit | #72030 | 384 rxns. |
| TIGIT:CD155 Homogeneous Assay Kit | #72029 | 384 rxns. |
| CD226:CD112 Homogeneous Assay Kit | #72051 | 384 rxns. |

AlphaScreen® and AlphaLISA® are registered trademarks of PerkinElmer, Inc.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
 Or you can Email us at: info@bpsbioscience.com
 Please visit our website at: www.bpsbioscience.com