



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court West, Ste. B
San Diego Ca 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet
TMPRSS2 Fluorogenic Assay Kit
Catalog # 78083
Size: 96 reactions

BACKGROUND: Transmembrane protease, serine 2 is an enzyme that belongs to the serine protease family. It has been implicated as a target in prostate cancer. TMPRSS2 facilitates SARS-CoV-2 particle entry into host cells via S protein priming, and its inhibition blocks virus fusion with angiotensin-converting enzyme 2 (ACE2). This, in turn, restricts SARS-CoV-2 viral entry, making TMPRSS2 an important therapeutic target.

DESCRIPTION: The *TMPRSS2 Fluorogenic Assay Kit* is provided in a convenient 96-well format, with purified TMPRSS2, TMPRSS2 Fluorogenic Substrate, and TMPRSS2 assay buffer for 96 enzymatic reactions. The key to the *TMPRSS2 Fluorogenic Assay Kit* is the fluorogenic substrate. Using this kit, only one simple step on a microtiter plate is required for TMPRSS2 reactions. A sample containing TMPRSS2 is incubated in a reaction mixture with the fluorogenic substrate and fluorescence ($\lambda_{ex} = 383 \text{ nm}$, $\lambda_{em} = 455 \text{ nm}$) is measured using a plate reader. Camostat is supplied as a protease inhibitor control.

COMPONENTS:

Catalog #	Component	Amount	Storage	
	TMPRSS2	15 μg	-80°C	Avoid freeze/thaw cycles!
78047	TMPRSS2 Fluorogenic Substrate (5 mM)	10 μl	-20°C	
78048	1X TMPRSS2 Assay Buffer	5 ml	-20°C	
78049	Camostat	500 μg	-80°C	
79685	96-well black plate	1	Room temp.	

MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

Fluorescent microplate reader capable of reading fluorescence at $\lambda_{ex} = 383 \pm 15 \text{ nm}$, $\lambda_{em} = 455 \pm 15 \text{ nm}$

Adjustable micropipettor and tips

Rotating or rocker platform

APPLICATIONS: Great for studying enzyme kinetics and screening small molecular inhibitors for drug discovery and HTS applications.

STABILITY: One year from date of receipt when stored as directed.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
San Diego Ca 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

REFERENCES:

1. Tomlins, S. A., *et al.* 2008. "Role of the TMPRSS2-ERG gene fusion in prostate cancer." *Neoplasia* **10 (2)**: 177-188.
2. Hoffmann, M., *et al.*, 2020. "SARS-CoV-2 Cell Entry Depends on ACE2 and TMPRSS2 and Is Blocked by a Clinically Proven Protease Inhibitor". *Cell* **181**: 271–280.

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate.

Step 1:

- 1) Thaw **TMPRSS2** on ice. Upon first thaw, briefly spin tube containing enzyme to recover full content of the tube. Aliquot **TMPRSS2** into single use aliquots. Store remaining undiluted enzyme in aliquots at -80°C. *Note: TMPRSS2 is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*
- 2) Dilute **TMPRSS2** in **1X TMPRSS2 Assay Buffer** at 5 ng/μl (150 ng/reaction). Keep diluted enzyme on ice until use. Discard any unused diluted enzyme after use.
- 3) Re-constitute the 500 μg of **Camostat** with 20 μl of water (50 mM). Dilute **Camostat** 1000-fold with assay buffer (50 μM). Add 10 μl to the wells labeled "Inhibitor Control." The final concentration of **Camostat** in the assay will be 10 μM.
- 4) Dispense 10 μl of Inhibitor solution to each well labeled as "Test Inhibitor." For the "Positive Control" and "Blank," add 10 μl of 5% DMSO in buffer (Inhibitor buffer). Maintain the same level of DMSO in the controls as the test sample(s).

Note: Final DMSO concentration must be ≤2%. Higher DMSO level can significantly decrease the enzyme activity. For example, to test an inhibitor at 10 μM that is dissolved in 100% DMSO, dilute 1 mM inhibitor with buffer to make 50 μM inhibitor in 5% DMSO (aqueous). Then, use 10 μl of the 50 μM solution for the 50 μl assay.

	Positive Control	Inhibitor Control	Test Inhibitor	Blank
Camostat	-	10 μl	-	-
Test Inhibitor	-	-	10 μl	-
Inhibitor buffer (5% DMSO in buffer)	10 μl	-	-	10 μl
1x TMPRSS2 Assay Buffer	-	-	-	30 μl
TMPRSS2 (5 ng/μl)	30 μl	30 μl	30 μl	-
TMPRSS2 Substrate (50 μM)	10 μl	10 μl	10 μl	10 μl
Total	50 μl	50 μl	50 μl	50 μl

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
San Diego Ca 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

- 5) Add 30 μ l of diluted **TMPRSS2** (5 ng/ μ l) to all wells except for "Blank". For "Blank," add 30 μ l of **1x TMPRSS2 Assay buffer**. Incubate plate for 30 minutes at room temperature.
- 6) Begin thawing **TMPRSS2 Fluorogenic Substrate (5 mM)** at room temperature.
Note: Protect this substrate from direct exposure to light!

Step 2:

Note: Protect TMPRSS2 Fluorogenic Substrate from direct exposure to light!

- 7) Dilute **TMPRSS2 Fluorogenic Substrate (5 mM)** 100-fold with **1x TMPRSS2 Assay buffer** to 50 μ M.
- 8) Initiate reactions by adding 10 μ l of **TMPRSS2 Fluorogenic Substrate (50 μ M)** to all wells. Final concentration of **TMPRSS2 Fluorogenic Substrate** is 10 μ M.
- 9) Protect plate from light by covering with aluminum foil. Incubate plate for ten minutes at room temperature.

Step 3:


- 10) Read fluorescence at $\lambda_{\text{ex}} = 383 (\pm 15)$ nm and $\lambda_{\text{em}} = 455 (\pm 15)$ nm. Alternatively, kinetic measurements can also be monitored for up to 30 minutes at 2-5 minute intervals. "Blank" value is subtracted from all measurements. *Note: Make sure the wells are not exposed to direct light.*

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com

Example of Assay Results:

TMPRSS2 Activity

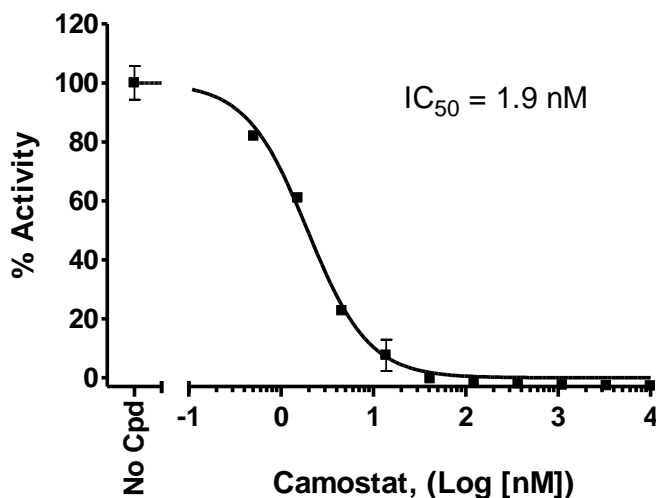

Inhibition of TMPRSS2 by Camostat (BPS Bioscience, #78049), measured using the *TMPRSS2 Fluorogenic Assay Kit* (BPS Bioscience #78083). *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com*

Related Products

<u>Product</u>	<u>Cat. #</u>	<u>Size</u>
Furin Protease Assay Kit	78040	96 rxns.
3CL Protease, Untagged (SARS-CoV-2) Assay Kit	78042-1	96 rxns.
3CL Protease, Untagged (SARS-CoV-2) Assay Kit	78042-2	384 rxns.
3CL Protease (SARS-CoV-2), no tag	100823-1	50 µg
3CL Protease (Mpro) (SARS-CoV-2)	100755	50 µg
3CL Protease (SARS-CoV-2) Assay Kit	79955	96 rxns.
3CL Protease, His-tag (SARS-CoV-1)	100807	100 µg
3CL Protease (SARS-CoV-1) Assay Kit	78015	96 rxns

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com