



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

DCAF11 Intrachain TR-FRET Assay Kit

Description

The DCAF11 Intrachain TR-FRET Assay Kit is a sensitive high-throughput screening (HTS) TR-FRET Assay Kit designed to measure DCAF11 auto-ubiquitination activity in a homogeneous 384-reaction format. It utilizes Europium-labeled Ubiquitin (donor) and Cy5-labeled Ubiquitin (acceptor) to complete the TR-FRET pairing. This assay measures poly-ubiquitination since both the TR-FRET donor and acceptor are incorporated into poly-ubiquitin chains formed on DCAF11. As a homogeneous assay, it requires no time-consuming washing steps, making it especially suitable for HTS applications as well as real-time analyses.

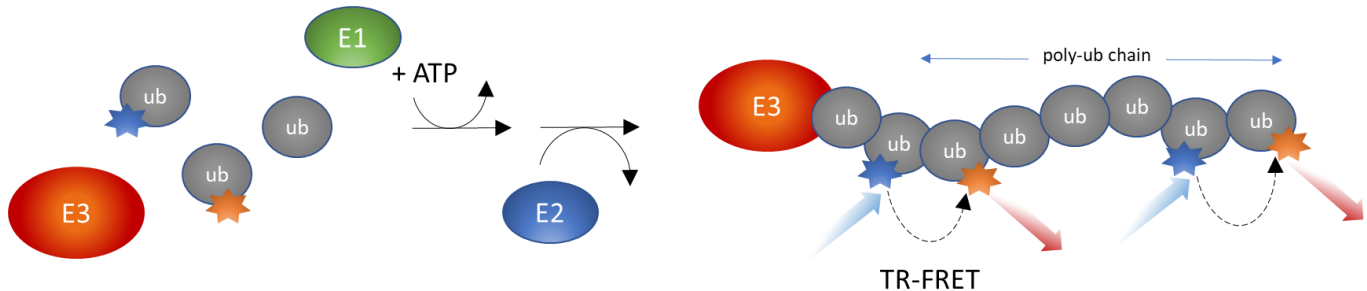


Figure 1. DCAF11 Intrachain TR-FRET Assay Kit Schematic.

Background

Covalent conjugation to ubiquitin (Ub) is one of the major post-translational modifications that regulates protein stability, function, and localization. Ubiquitination is the concerted action of three enzymes: a Ub-activating enzyme (E1), a Ub-conjugating enzyme (E2), and a Ub ligase (E3). The specificity and efficiency of ubiquitination are largely determined by the E3 enzyme, which directs the last step of the Ub-conjugating cascade by binding to both an E2~Ub conjugate and a substrate protein. This step ensures the transfer of Ub from E2~Ub to the substrate, leading to its mono- or poly-ubiquitination.

DDB1 and CUL4-associated Factor 11 (DCAF11) is a protein that associates with CUL4A or CUL4B to form a complex with E3 ligase activity. DCAF11 interacts with and ubiquitinates nuclear factor E2-related factor 2 (Nrf2), which is known to regulate cellular stress response genes, including those associated with oxidative stress. Additionally, DCAF11 activity is repressed by transcriptional factor EB (TFEB), allowing for stabilization of proteins such as Nrf2. Because Nrf2 and TFEB are involved in misfolded protein clearance, and oxidative stress is known to contribute to diseases such as Alzheimer's, Huntington's, and Parkinson's disease, DCAF11 is a potential target for the treatment of multiple neurodegenerative disorders. Like most E3 ligases, DCAF11 ubiquitinates itself.

Application(s)

- Screen molecules that inhibit DCAF11 Ubiquitin ligase activity in drug discovery HTS applications
- Determine Inhibitor IC₅₀
- Perform DCAF11 real-time kinetics.

Supplied Materials

Catalog #	Name	Amount	Storage	
80301	UBE1 (UBA1), FLAG-Tag (E1)*	50 µg	-80°C	Avoid multiple freeze/thaw cycles
80314	UbcH5b, His-Tag (E2)*	60 µg	-80°C	
101495	DCAF11/Rbx1/CUL4B/DDB1 complex, (E3)*	62 µg	-80°C	
78307	TRF Ubiquitin Mix (200x)	40 µl	-80°C	
	ATP (4 mM)	2 x 1 ml	-80°C	
	CBL assay buffer 2	2 x 10 ml	-80°C	
79969	White, nonbinding Corning, low volume 384-well microtiter plate		Room Temp	

* The initial concentration of enzyme is lot-specific and will be indicated on the tube containing the protein.

Materials Required but Not Supplied

- Fluorescent microplate reader capable of measuring Time Resolved Fluorescence Resonance Energy Transfer (TR-FRET)
- Adjustable micropipettor and sterile tips
- Rotating or rocker platform

Storage Conditions

This assay kit will perform optimally for up to 6 months from date of receipt when the materials are stored as directed. **Avoid multiple freeze/thaw cycles!**

Safety

This product is for research purposes only and not for human or therapeutic use. This product should be considered hazardous and is harmful by inhalation, in contact with skin, eyes, clothing, and if swallowed. If contact occurs, wash thoroughly.

Contraindications

The DCAF11 Intrachain TR-FRET Assay Kit is compatible with up to 1% final DMSO concentration. We recommend preparing the inhibitor in no higher than 5% DMSO solution in buffer and using 4 µl per well.

Assay Protocol

- All samples and controls should be performed in triplicates
 - The assay should include a “Blank”, a “Positive control”, and a “Negative control”
1. Thaw **UBE1**, **UBCH5b**, **DCAF11**, **Ubiquitin Mix**, **Assay Buffer**, and **ATP** on ice. Briefly spin the tubes to recover their full contents.
 2. Prepare a 5x Ubiquitin Mix in Assay Buffer by making a 40-fold dilution of the stock 200x TRF Ubiquitin Mix.

Example: Add 1 volume of stock Ubiquitin Mix to 39 volumes of Assay Buffer.

3. Calculate the amount of protein required for the assay and dilute enough for the assay. Refer to the protocol below to calculate how much of each protein is needed.

4. Prepare the appropriate amounts of diluted proteins. The concentration of each protein is lot-specific and is indicated on the tube. Verify the initial concentration and dilute accordingly.
 - a. Dilute UBE1 in Assay Buffer to 96 ng/ μ l (800 nM - the final concentration in the reaction will be 40 nM)
 - b. Dilute UBCH5b in Assay Buffer to 144 ng/ μ l (8 μ M - the final concentration in the reaction will be 400 nM)
 - c. Dilute the DCAF11 complex in Assay Buffer to 31 ng/ μ l (100 nM – the final concentration in the reaction will be 25 nM)

Note: Keep all diluted proteins on ice until use. Do not freeze and re-use diluted proteins.

5. Aliquot the remainder of each protein, the assay buffer, and the ATP into 2-4 single-use aliquots as may be necessary and immediately store at -80°C .

Note: UBE1, UBCH5b, DCAF11, TRF Ubiquitin Mix, and Assay Buffer are sensitive to freeze/thaw cycles. Avoid multiple freeze-thaw cycles.

6. Prepare the Test Inhibitor (4 μ l/well): for a titration, prepare serial dilutions at concentrations 5-fold higher than the desired final concentrations. The final volume of the reaction is 20 μ l.

Without DMSO

- a) If the Test Inhibitor is water-soluble, prepare serial dilutions in the Assay Buffer, 5-fold more concentrated than the desired final concentrations. For the positive and negative controls, use 1x Assay Buffer (Diluent Solution).

Or

With DMSO

- a) If the Test inhibitor is soluble in DMSO, prepare the test inhibitor at a concentration 100-fold higher than the highest desired concentration in DMSO, then dilute the inhibitor 20-fold in Assay Buffer to prepare the highest concentration of the 5-fold intermediate dilutions. The concentration of DMSO is now 5%.
 - b) Prepare serial dilutions of the Test Inhibitor at 5-fold the desired final concentrations using 5% DMSO in Assay Buffer to keep the concentration of DMSO constant.
 - c) For positive and negative controls, prepare 5% DMSO in water (vol/vol) so that all wells contain the same amount of DMSO (Diluent Solution).
The final concentration of DMSO should not exceed 1%.
7. To the wells designated as “Blank”, add 4 μ l of **5x Ubiquitin Mix** + 1 μ l of **UBE1** + 1 μ l of **UBCH5b** + 4 μ l of **diluent solution** (for example DMSO 5% in assay buffer) + 5 μ l of **Assay Buffer**.

Component	Blank
Ubiquitin Mix (5x)	4 μ l
UBE1	1 μ l
UBCH5b	1 μ l
DCAF11	-
Test Inhibitor	-
Diluent solution* (no inhibitor)	4 μ l
Assay Buffer	5 μ l
ATP (4 mM)	5 μ l
Total	20 μl

*The diluent solution contains the assay buffer with the same concentration of solvent (e.g., DMSO) as the Test Inhibitor solution.

8. Prepare a Master Mix using diluted reagents prepared in steps 2 and 3:

N wells \times (4 μ l **5x Ubiquitin Mix** + 1 μ l **UBE1** + 1 μ l **UBCH5b** + 5 μ l **DCAF11**).

9. Add 11 μ l of Master Mix to each well designated “Negative Control”, “Positive Control”, and “Test Sample”.
10. Add 4 μ l of Test Inhibitor solution to each well designated “Test Inhibitor”. For “Positive Control” and “Negative Control”, add 4 μ l of the diluent solution without inhibitor.
11. Initiate the reaction by adding 5 μ l of **ATP** to the wells labeled “Positive Control,” “Test Inhibitor,” and “Blank.” Add 5 μ l of **Assay Buffer** to the well designated “Negative Control.” Cover the plate with a plate sealer. Incubate the reaction at room temperature for two hours or at 30°C for one hour.

Component	Test Inhibitor	Negative Control	Positive Control
Master Mix	11 μ l	11 μ l	11 μ l
Test Inhibitor	4 μ l	-	-
Diluent solution* (no inhibitor)	-	4 μ l	4 μ l
Assay Buffer	-	5 μ l	-
ATP (4 mM)	5 μ l	-	5 μ l
Total	20 μl	20 μl	20 μl

*The diluent solution contains the assay buffer with the same concentration of solvent (e.g. DMSO) as the test Inhibitor solution.

12. Read the fluorescent intensity in a microtiter-plate reader capable of measuring TR-FRET. “Blank” value is subtracted from all other values.

Instrument Settings

Eu Donor Emission	Time Resolved	Dye-acceptor Emission	Time Resolved
Excitation Wavelength	317±20 nm	Excitation Wavelength	317±20 nm
Emission Wavelength	620±10 nm	Emission Wavelength	665±10 nm
Lag Time	60 µs	Lag Time	60 µs
Integration Time	500 µs	Integration Time	500 µs

Calculating Results:

Two sequential measurements should be conducted. Eu donor emission should be measured at 620 nm followed by dye-acceptor emission at 665 nm. Data analysis is performed using the TR-FRET ratio (665 nm emission/620 nm emission).

When percentage activity is calculated, the FRET value from the Blank (it is expected that Blank and Negative Control represent similar value) can be set as zero percent activity and the FRET value from the positive control can be set as one hundred percent activity.

$$\% \text{ Activity} = \frac{\text{FRET}_s - \text{FRET}_{\text{blank}}}{\text{FRET}_p - \text{FRET}_{\text{blank}}} \times 100\%$$

Where FRET_s = Sample FRET, FRET_{blank} = Blank FRET, and FRET_p = Positive control FRET.

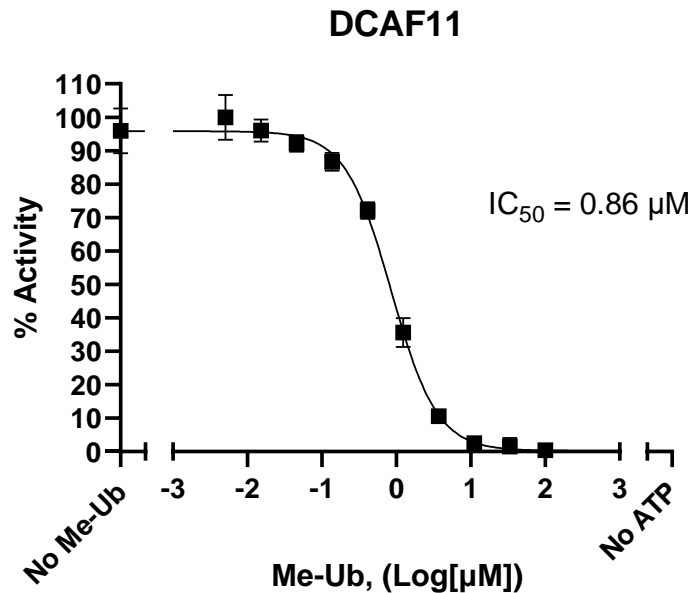
Example Results

Figure 1: Inhibition of DCAF11 auto-ubiquitination.

DCAF11 auto-ubiquitination was measured in presence of increasing concentrations of Methylated Ubiquitin using the DCAF11 Intrachain TR-FRET Assay Kit, BPS Bioscience #78542.

Data shown is representative. For lot-specific information, please contact BPS Bioscience, Inc. at support@bpsbioscience.com.

Troubleshooting Guide

Visit bpsbioscience.com/assay-kits-faq for detailed troubleshooting instructions. For all further questions, please email support@bpsbioscience.com

Related Products

<i>Products</i>	<i>Catalog #</i>	<i>Size</i>
Cereblon Intrachain TR-FRET Assay Kit	78301	384 reactions
MDM2 Intrachain TR-FRET Assay Kit	78302	384 reactions
SMURF1 Intrachain TR-FRET Assay Kit	78303	384 reactions
VHL Intrachain TR-FRET Assay Kit	78305	384 reactions
XIAP Intrachain TR-FRET Assay Kit	78306	384 reactions
DCAF15 Intrachain TR-FRET Assay Kit	78543	384 reactions