



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Description

The Chemi-Verse™ PARP7 Assay Kit is an ELISA-type chemiluminescent assay designed to measure PARP7 enzymatic activity for screening and profiling applications. PARP7 catalyzes the NAD-dependent ADP-ribosylation of histones. This 96-well format assay kit contains sufficient amounts of purified PARP7 enzyme (amino-acids 400-657), histone mixture, and PARP assay buffer for 100 enzyme reactions. The Chemi-Verse™ PARP7 Assay Kit takes advantage of a highly specific ADP-ribose binding reagent. First, histone proteins are coated on a 96-well plate. Next, an NAD⁺ substrate is incubated with the PARP7 enzyme in an optimized assay buffer. Finally, the plate is treated with the ADP-ribose specific binding reagent and secondary HRP-conjugated antibody, followed by addition of the ELISA ECL substrate to produce chemiluminescence that can be measured using a chemiluminescence reader.

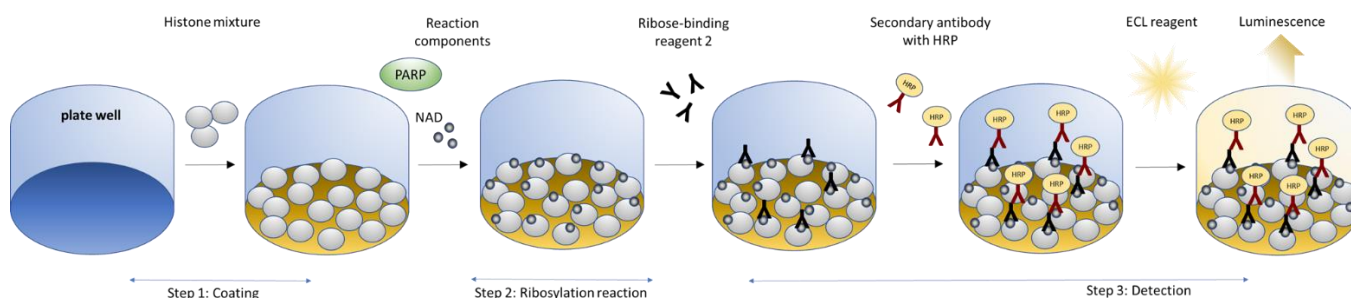


Figure 1. Chemi-Verse™ PARP7 chemiluminescent assay schematic.

Applications

Study enzyme kinetics and screening small molecule inhibitors for drug discovery and HTS applications.

Supplied Materials

| Catalog # | Name | Amount | Storage | |
|-----------|--|-----------|-----------|--|
| 80527 | PARP7, FLAG-Tag (Amino-acids 400-657)* | 2 µg | -80°C | Avoid multiple freeze/thaw cycles |
| 52029 | 5x Histone Mixture | 1 ml | -80°C | |
| | NAD ⁺ (50 µM) | 250 µl | -80°C | |
| | PP-16-2 Assay Buffer | 1 ml | -20°C | |
| 79743 | Blocking Buffer 3 | 2 x 25 ml | +4°C | |
| | Ribose-Binding Reagent 2-D | 10 µl | -80°C | |
| 52131H | Secondary HRP--Antibody 2 | 10 µl | -80°C | |
| 79670 | ELISA ECL Substrate A (translucent bottle) | 6 ml | Room Temp | |
| | ELISA ECL Substrate B (brown bottle) | 6 ml | Room Temp | |
| 79837 | 96-well module plate | | Room Temp | |

*The concentration of the protein is lot-specific and will be indicated on the tube.

Materials Required but Not Supplied

| Name | Ordering Information |
|--|----------------------|
| DTT (10 mM in water, prepared fresh) | |
| 1x PBS (phosphate buffer saline) | |
| PBST buffer (1x PBS, containing 0.05% Tween-20) | |
| Luminometer or plate reader capable of reading chemiluminescence | |
| Adjustable micropipettor and sterile tips | |
| Rotating or rocker platform | |

Storage Conditions

This assay kit will perform optimally for up to 6 months from date of receipt when the materials are stored as directed. **Avoid multiple freeze/thaw cycles!**

Safety

This product is for research purposes only and not for human or therapeutic use. This product should be considered hazardous and is harmful by inhalation, in contact with skin, eyes, clothing, and if swallowed. If contact occurs, wash thoroughly.

Contraindications

The Chemi-Verse™ PARP7 Chemiluminescent Assay Kit is compatible with up to 1% final DMSO concentration. We recommend preparing the inhibitor in no higher than 10% DMSO solution in buffer and using 5 µl per well.

Assay Protocol

- All samples and controls should be performed in duplicates.
- The assay should include a “Blank” and a “Positive control.”

Step 1: Coat a 96-well module (VWR catalog no. 62409-300) with 50 µl per well of histone solution

- 1) Dilute 5x histone mixture 1:5 with PBS to make 1x histone mixture.
- 2) Add 50 µl of histone mixture to each well and incubate at 4°C overnight.
- 3) Wash the plate three times using 200 µl/well of PBST buffer (1x PBS containing 0.05% Tween 20).
- 4) Tap the plate onto a clean paper towel to remove the liquid.
- 5) Block the wells by adding 200 µl of Blocking buffer 3 to every well. Incubate at room temperature for at least 90 minutes.
- 6) Wash the plate three times with 200 µl/well of PBST buffer.

- 7) Tap the plate onto a clean paper towel to remove the liquid.

Step 2: Ribosylation reaction

- 1) Prepare a fresh solution of 10 mM DTT in water.
- 2) Prepare the Master Mix (25 µl/well): N wells x (2.5 µl of PP-16-2 buffer + 2.5 µl of NAD + 17.5 µl of water + 2.5 µl of 10 mM fresh DTT).

Note: The concentration of DTT in the Master Mix is 1 mM.

- 3) Add 25 µl of Master Mix to every well.
- 4) Prepare 1x Assay Buffer with DTT. Dilute the stock of PP-16-2 buffer 10-fold by adding 1 volume of PP-16-2 buffer to 1 volume of 10 mM DTT and 8 volumes of water.

Note: The concentration of DTT in the 1x Assay Buffer is 1 mM

- 5) Prepare the Test Inhibitor (5 µl/well): for a titration, prepare serial dilutions at concentrations 10-fold higher than the desired final concentrations. The final volume of the reaction is 50 µl.

Without DMSO

- 5.1. If the Test Inhibitor is water-soluble, prepare serial dilutions in the 1x Assay Buffer, 10-fold more concentrated than the desired final concentrations.

With DMSO

- 5.2. If the Test inhibitor is soluble in DMSO, prepare the test inhibitor at 100-fold the highest desired concentration in DMSO, then dilute the inhibitor 10-fold in 1x Assay Buffer to prepare the highest concentration of the 10-fold intermediate dilutions. The concentration of DMSO is now 10%.
 - a. Using 1x Assay Buffer in 10% DMSO, prepare serial dilutions of the Test Inhibitor at 10-fold the desired final concentrations to keep the concentration of DMSO constant.
 - b. For positive and negative controls, prepare 10% DMSO in 1x Assay Buffer (vol/vol) so that all wells contain the same amount of DMSO (Diluent Solution).

Note: The final concentration of DMSO in the assay should not exceed 1%.

- 6) Add 5 µl of Test Inhibitor to each well labeled as "Test Inhibitor."
 - a. For the "Positive Control" and "Blank," add 5 µl of the diluent solution used to dilute the inhibitor, but without inhibitor (Diluent Solution).

- 7) Thaw PARP7 enzyme on ice. Briefly spin the tube containing the enzyme to recover the full content of the tube. Calculate the amount of PARP7 required for the assay and dilute enzyme to **0.78 ng/μl** with 1x Assay Buffer containing DTT. The final concentration of PARP7 will be 10 nM. Aliquot the remaining undiluted PARP7 enzyme into aliquots and store at -80°C. Do not re-use these aliquots more than once.

*Note: PARP7 enzyme is very sensitive to freeze/thaw cycles. Avoid multiple freeze/thaw cycles. **Do not re-use the diluted enzyme.***

- 8) Initiate the reaction by adding 20 μl of diluted PARP7 enzyme to the wells designated "Positive Control" and "Test Inhibitor."
- To the wells designated as "Blank," add 20 μl of 1x assay buffer with DTT.
 - Incubate at room temperature for 1 hour.

| | Blank | Positive Control | Test Inhibitor |
|--------------------------|--------------|-------------------------|-----------------------|
| Master Mix | 25 μl | 25 μl | 25 μl |
| Test Inhibitor | - | - | 5 μl |
| Diluent Solution | 5 μl | 5 μl | - |
| 1x Assay Buffer with DTT | 20 μl | - | - |
| PARP7 (0.78 ng/μl) | - | 20 μl | 20 μl |
| Total | 50 μl | 50 μl | 50 μl |

- 9) Wash the plate three times with 200 μl of PBST buffer and tap the plate onto a clean paper towel to remove the excess liquid.

Step 3: Detection

- Dilute the Ribose-binding reagent 1,000-fold in Blocking Buffer 3.
- Add 50 μl of diluted Ribose-binding reagent to each well. Incubate for 45 minutes at room temperature.
- Wash three times with 200 μl of PBST buffer and tap the plate onto a clean paper towel.
- Dilute the Secondary HRP-Antibody 5,000-fold in Blocking buffer 3.
- Add 50 μl of diluted Secondary HRP-Antibody to each well. Incubate for 30 minutes at room temperature.
- Wash three times with 200 μl of PBST buffer and tap the plate onto a clean paper towel.
- Just before use, mix 1 volume of ELISA ECL Substrate A and 1 volume of ELISA ECL Substrate B. Add 100 μl per well.
- Immediately read the plate in a luminometer or microtiter-plate reader capable of reading chemiluminescence. The "Blank" value is subtracted from all other values.

Reading Chemiluminescence

Chemiluminescence is the emission of light (luminescence) which results from a chemical reaction. The detection of chemiluminescence requires no wavelength selection because the method used is emission photometry and is not emission spectrophotometry. To properly read chemiluminescence, make sure the plate reader is set for LUMINESCENCE mode. Typical integration time is 1 second, delay after plate movement is 100 msec. Do not use a filter when measuring light emission. Typical settings for the Synergy 2 BioTek plate reader are: use the “hole” position on the filter wheel; Optics position: Top; Read type: endpoint. Sensitivity may be adjusted based on the luminescence of a control assay without enzyme (typically we set this value as 100).

Validation Data

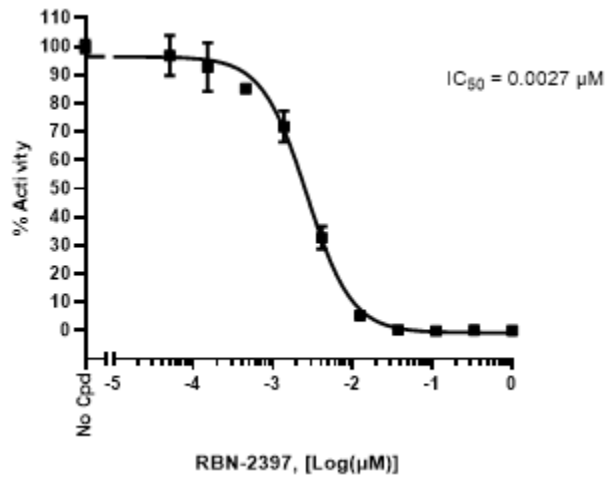


Figure 2: Inhibition of PARP7 activity by RBN-2397.

PARP7 activity was measured in the presence of increasing concentrations of RBN-2397 (MedKool Biosciences #462554) using the Chemi-Verse™ PARP7 Assay Kit (BPS Bioscience #78814). Luminescence was measured using a Bio-Tek microplate reader.

Data shown is representative, for lot-specific information, please contact BPS Bioscience, Inc. at support@bpsbioscience.com.

Troubleshooting Guide

Visit bpsbioscience.com/assay-kits-faq for detailed troubleshooting instructions. For all further questions, please email support@bpsbioscience.com

Related Products

| <i>Products</i> | <i>Catalog #</i> | <i>Size</i> |
|---|------------------|--------------|
| PARP1 Chemiluminescent Assay Kit | 80551 | 96 reactions |
| PARP2 Chemiluminescent Assay Kit | 80552 | 96 reactions |
| PARP3 Chemiluminescent Assay Kit | 80553 | 96 reactions |
| PARP6 Chemiluminescent Assay Kit | 80556 | 96 reactions |
| PARP7 Chemiluminescent Assay Kit | 79729 | 96 reactions |
| PARP10 Chemiluminescent Assay Kit | 80560 | 96 reactions |
| PARP11 Chemiluminescent Assay Kit | 80561 | 96 reactions |
| PARP12 Chemiluminescent Assay Kit | 78504 | 96 reactions |
| PARP14 Chemiluminescent Assay Kit | 80568 | 96 reactions |
| PARP15 Chemiluminescent Assay Kit | 80567 | 96 reactions |
| PARP15FL Chemiluminescent Assay Kit | 78596 | 96 reactions |
| TNKS1 (PARP5A) Chemiluminescent Assay Kit | 78405 | 96 reactions |
| TNKS2 (PARP5B) Chemiluminescent Assay Kit | 78406 | 96 reactions |
| PARP1 Enzyme | 80501 | 20 µg |
| PARP2 Enzyme | 80502 | 10 µg |
| PARP3 Enzyme | 80503 | 10 µg |
| PARP6 Enzyme | 80506 | 10 µg |
| PARP7 Enzyme | 80527 | 10 µg |
| PARP10 Enzyme | 80522 | 10 µg |
| PARP11 Enzyme | 80511 | 10 µg |
| PARP12 Enzyme | 80513 | 10 µg |
| PARP14 Enzyme | 80514 | 10 µg |
| PARP15 Enzyme | 80517 | 10 µg |
| TNKS1 Enzyme | 80504 | 10 µg |
| TNKS2 Enzyme | 80515 | 10 µg |