



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court West, Ste. B
San Diego CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet
METTL3/METTL14 Complex
Chemiluminescent Assay Kit
Catalog #79614
Size: 96 reactions

DESCRIPTION: The *METTL3/METTL14 Complex Chemiluminescent Assay Kit* is designed to measure METTL3/METTL14 complex activity for screening and profiling applications. The *METTL3/METTL14 Complex Chemiluminescent Assay Kit* comes in a convenient format, with 96-well plate precoated with a specific substrate, the antibody against methylated adenosine, a secondary HRP-labeled antibody, S-adenosylmethionine, methyltransferase assay buffer, and purified *METTL3/METTL14 complex* enzyme for 96 enzyme reactions. The key to the *METTL3/METTL14 Complex Chemiluminescent Activity Assay Kit* is a highly specific antibody that recognizes N6-methylated adenosine. With this kit, only three simple steps are required for methyltransferase detection. First, S-adenosylmethionine is incubated with a sample containing assay buffer and methyltransferase enzyme. Next, primary antibody is added. Finally, the plate is treated with an HRP-labeled secondary antibody followed by addition of the HRP substrate to produce chemiluminescence that can then be measured using a chemiluminescence reader.

COMPONENTS:

Catalog #	Component	Amount	Storage	
100105	METTL3/METTL14 human enzyme	50 µg	-80°C	(Avoid freeze/thaw cycles!)
	20 µM S-adenosylmethionine	250 µl	-80°C	
52140Z4	Primary antibody 29	100 µl	-80°C	
52131H	Secondary HRP-labeled antibody 2	10 µl	-80°C	
52193Z	4x METTL assay buffer*	3 ml	-20°C	
52100B	Blocking buffer	50 ml	+4°C	
	HRP chemiluminescent substrate A (translucent bottle)	6 ml	+4°C	
	HRP chemiluminescent substrate B (brown bottle)	6 ml	+4°C	
	White microplate precoated with the RNA substrate*	1	+4°C	

***Important: Use RNase-free conditions!**

***Add 30 µl of 0.5 M DTT.**

MATERIALS REQUIRED BUT NOT SUPPLIED:

TBST buffer (1 x Tris-buffered saline (TBS), pH 8.0, containing 0.05% Tween-20)

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
San Diego CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Dithiothreitol (0.5 M DTT)
Luminometer or fluorescent microplate reader capable of reading chemiluminescence
Adjustable micropipettor and sterile tips
Rotating or rocker platform

APPLICATIONS: Great for studying enzyme kinetics and HTS applications.

CONTRAINDICATIONS: DMSO >1%, strong acids or bases, ionic detergents, high salt

STABILITY: One year from date of receipt when stored as directed.

REFERENCE: Liu, J., *et al. Nat Chem Biol.* 2014; **10(2)**: 93–95.

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate.

Step 1:

- 1) Rehydrate the microwells by adding 150 μ l of TBST buffer (1x TBS, pH 8.0, containing 0.05% Tween-20) to every well. Incubate 15 minutes at room temperature. Tap the strip onto clean paper towels to remove liquid.
- 2) Thaw **S-adenosylmethionine** on ice. Upon first thaw, briefly spin tube containing S-adenosylmethionine to recover full content of the tube. Aliquot S-adenosylmethionine into single use aliquots. Store remaining S-adenosylmethionine in aliquots at -80°C immediately. *Note: **S-adenosylmethionine** is sensitive to freeze/thaw cycles. Avoid multiple freeze-thaw cycles.*
- 3) Add 30 μ l of 0.5 M DTT to 4x METTL assay buffer. Prepare the master mixture: N wells \times (7.5 μ l **4x METTL assay buffer** + 2.5 μ l 20 μ M **S-adenosylmethionine** + 15 μ l distilled **H₂O**).
- 4) Add 25 μ l of master mixture to each well designated for the "Positive Control," "Test Inhibitor," and "Blank." For the "Substrate Control," add 7.5 μ l **4x METTL assay buffer** + 17.5 μ l **H₂O**.
- 5) Thaw **METTL3/METTL14 complex enzyme** on ice. Upon first thaw, briefly spin tube containing enzyme to recover full content of the tube. Aliquot **METTL3/METTL14 complex enzyme** into single use aliquots. Store remaining undiluted enzyme in aliquots at -80°C immediately. *Note: **METTL3/METTL14 complex***

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
San Diego CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

enzyme is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.

- 6) Dilute **METTL3/METTL14 complex enzyme** in **1x METTL assay buffer** to 25 ng/ μ l (500 ng/20 μ l). Keep diluted enzyme on ice until use. Discard any unused diluted enzyme after use.

	Blank	Substrate Control	Positive Control	Test Inhibitor
4x METTL assay buffer	7.5 μ l	7.5 μ l	7.5 μ l	7.5 μ l
20 μ M S-adenosylmethionine	2.5 μ l	–	2.5 μ l	2.5 μ l
H ₂ O	15 μ l	17.5 μ l	15 μ l	15 μ l
Test Inhibitor/Activator	–	–	–	5 μ l
10% DMSO in water (Inhibitor buffer)	5 μ l	5 μ l	5 μ l	–
1x METTL assay buffer	20 μ l	–	–	–
METTL3/METTL14 Complex (25 ng/ μ l)	–	20 μ l	20 μ l	20 μ l
Total	50 μl	50 μl	50 μl	50 μl

- 7) Add 5 μ l of inhibitor solution of each well designated “Test Inhibitor.” For the “Positive Control,” “Substrate Control,” and “Blank,” add 5 μ l of 10% DMSO in water (**Inhibitor buffer**).
- 8) Add 20 μ l of **1x METTL buffer** to the well designated “Blank.”
- 9) Initiate reaction by adding 20 μ l of diluted **METTL3/METTL14 complex** (prepared as described above) to the wells labeled “Test Inhibitor,” “Positive Control,” and “Substrate Control.” Incubate overnight at room temperature on a rotating platform. Seal the wells if necessary.
- 10) Wash the wells three times with 200 μ l TBST buffer. Blot dry onto clean paper towels.
- 11) Add 100 μ l of **Blocking buffer** to every well. Shake on a rotating platform for 10 min. Remove supernatant as above.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
San Diego CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Step 2:

- 1) Dilute **Primary antibody 29** 100-fold with **Blocking buffer**.
- 2) Add 100 μ l per well. Incubate 1 hour at room temperature with slow shaking.
- 3) Wash plate three times with 200 μ l TBST buffer and incubate in **Blocking buffer** as in steps 1-10 and 1-11.

Step 3:

- 1) Dilute **Secondary HRP-labeled antibody 2** 1,000-fold with **Blocking buffer**.
- 2) Add 100 μ l per well. Incubate for 30 minutes at room temperature with slow shaking.
- 3) Wash plate three times with TBST buffer and incubate in **Blocking buffer** as in steps 1-10 and 1-11.
- 4) Just before use, mix on ice 50 μ l **HRP chemiluminescent substrate A** and 50 μ l **HRP chemiluminescent substrate B** and add 100 μ l per well. Discard any unused chemiluminescent reagent after use.
- 5) Immediately read sample in a luminometer or microtiter-plate reader capable of reading chemiluminescence. "Blank" value is subtracted from all readings.

Reading Chemiluminescence:

Chemiluminescence is the emission of light (luminescence) which results from a chemical reaction. The detection of chemiluminescence requires no wavelength selection because the method used is emission photometry and is not emission spectrophotometry.


To properly read chemiluminescence, make sure the plate reader is set for LUMINESCENCE mode. Typical integration time is 1 second, delay after plate movement is 100 msec. Do not use a filter when measuring light emission. Typical settings for the Synergy 2 BioTek plate reader are: use the "hole" position on the filter wheel; Optics position: Top; Read type: endpoint. Sensitivity may be adjusted based on the luminescence of a control assay without enzyme (typically we set this value as 100).

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com

Example of Assay Results:

METTL3/METTL14 Activity

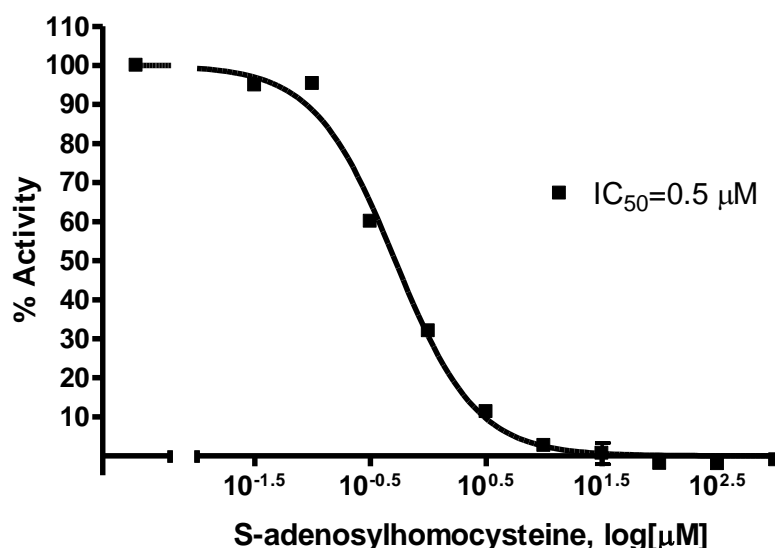

METTL3/METTL14 complex enzyme activity, measured using the *METTL3/METTL14 Complex Chemiluminescent Assay Kit*, BPS Bioscience Catalog #79614 Luminescence was measured using a Bio-Tek fluorescent microplate reader. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com.*

RELATED PRODUCTS

<u>Product Name</u>	<u>Catalog #</u>	<u>Size</u>
NSD2 enzyme (catalytic)	#51026	50 μg
NSD2 (782-end)/ReBP11 enzyme	#51025	20 μg
NSD1 enzyme	#51024	50 μg
NSD3 (1021-1322) enzyme	#51036	50 μg
SETD2 enzyme	#53019	50 μg
SETD2 Chemiluminescent Assay Kit	#52060	96 reactions
Chaetocin	#27221	1 mg
4x HMT Assay Buffer 7	#52193	30 mL
FBXL10(KDM2B, JHDM1B) enzyme	#50120	20 μg

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
 Or you can Email us at: info@bpsbioscience.com
 Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
San Diego CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

FBXL11(KDM2A) enzyme

#50102

20 µg

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court West, Ste. B
 San Diego CA 92121
 Tel: 1.858.829.3082
 Fax: 1.858.481.8694
 Email: info@bpsbioscience.com

TROUBLESHOOTING GUIDE

Problem	Possible Cause	Solution
Luminescence signal of positive control reaction is weak	METTL3/METTL14 complex enzyme has lost activity	Enzyme loses activity upon repeated freeze/thaw cycles. Use fresh enzyme (METTL3, BPS Bioscience #100105). Store enzyme in single-use aliquots. Increase time of enzyme incubation. Increase enzyme concentration.
	Antibody reaction is insufficient	Increase time for antibody incubation. Avoid freeze/thaw cycles of antibody.
	Incorrect settings on instruments	Refer to instrument instructions for settings to increase sensitivity of light detection.
	Chemiluminescent reagents mixed too soon	Chemiluminescent solution should be used within 15 minutes of mixing. Ensure both reagents are properly mixed.
Luminescent signal is erratic or varies widely among wells	Inaccurate pipetting/technique	Run duplicates of all reactions. Use a multichannel pipettor. Use master mixes to minimize errors.
	Bubbles in wells	Pipette slowly to avoid bubble formation. Tap strip lightly to disperse bubbles; be careful not to splash between wells.
Background (signal to noise ratio) is high	Insufficient washes	Increase number of washes. Increase wash volume. Increase Tween-20 concentration to 0.1% in TBST.
	Sample solvent is inhibiting the enzyme	Run negative control assay including solvent. Maintain DMSO level at <1%. Increase time of enzyme incubation.
	Results are outside the linear range of the assay	Use different concentrations of enzyme (METTL3, BPS Bioscience #100105) to create a standard curve.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082**, Fax **1.858.481.8694**
 Or you can Email us at: info@bpsbioscience.com
 Please visit our website at: www.bpsbioscience.com