



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

CSL Reporter – HEK293 Cell line Notch Signaling Pathway Catalog #: 79754

Background

The Notch signaling pathway controls cell fate decisions in vertebrate and invertebrate tissues. Notch signaling is triggered through the binding of a transmembrane ligand to Notch transmembrane receptor (NOTCH1/ NOTCH2/NOTCH3/NOTCH4) on a neighboring cell. This results in proteolytic cleavage of the Notch receptor, releasing the constitutively active intracellular domain of NOTCH (NICD). NICD translocates to the nucleus and associates with transcription factors CSL (CBF1/RBPJk/Suppressor of Hairless/Lag-1) and coactivator Mastermind to turn on transcription of Notch-responsive genes.

Description

The Notch CSL Reporter – HEK293 cell line contains the firefly luciferase gene under the control of Notch-response elements (CSL responsive elements) stably integrated into HEK293 cells. Transfection of this cell line with a Notch expression vector and activation of the Notch pathway leads to expression of luciferase reporter. This cell line is validated with a Notch1DeIE expression vector, which is constitutively processed by γ -secretase, leading to the release of NICD and the expression of luciferase.

Application

- Monitor Notch signaling pathway activity.
- Screen for inhibitors of the Notch signaling pathway.
- Test synthetic Notch receptors (synNotch) or other genetic constructs that conditionally activate the CSL promoter

Format

Each vial contains $\sim 1.5 \times 10^6$ cells in 1 ml of FBS with 10% DMSO.

Storage

Immediately upon receipt, store cells in liquid nitrogen.

General Culture Conditions

Thaw Medium 1 (BPS Bioscience #60187): MEM medium (Hyclone #SH30024.01) supplemented with 10% FBS (Thermo Fisher, #26140079), 1% non-essential amino acids (Hyclone #SH30238.01), 1 mM Na pyruvate (Hyclone #SH30239.01), 1% Penicillin/Streptomycin (Hyclone #SV30010.01).

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Growth Medium 1B (BPS Bioscience #79531): Thaw Medium 1, 400 µg/ml Geneticin® (G418 Sulfate; Thermo Fisher, #11811031).

Cells should be grown at 37°C with 5% CO₂ in Growth Medium 1B. If using different medium components, it may be necessary to adjust the percentage of CO₂ in the incubator depending on the NaHCO₃ level in the basal medium.

Cell Thawing: Thaw the frozen cells in a 37°C water-bath for 40 seconds, until freezing medium is partially thawed but with some ice remaining. Transfer to a tube containing 10 ml of Thaw Medium 1 (**no geneticin**), and spin down cells at 200 x g for 5 minutes. Remove supernatant and resuspend cells in pre-warmed Thaw Medium 1 (**no geneticin**). Transfer resuspended cells to a T25 flask and culture in a 5% CO₂ incubator at 37°C overnight. The next day, replace the medium with fresh Thaw Medium 1 (**no geneticin**) and continue growing culture in a CO₂ incubator at 37°C until the cells are ready to be split. At first passage, switch to Growth Medium 1B (**contains geneticin**). Cells should be split before they reach complete confluence.

To passage the cells, rinse cells with phosphate buffered saline (PBS), and detach cells from the culture vessel with 0.05% Trypsin/EDTA. Add Growth Medium 1B and transfer to a tube. Spin down the cells, then resuspend cells and seed appropriate aliquots of cell suspension into new culture vessels. Subcultivation ration: 1:5 to 1:7 twice a week. *Note: This cell line will grow more slowly than wildtype HEK293 cells.*

To freeze down the cells, rinse cells with phosphate buffered saline (PBS), and detach cells from culture vessel with 0.05% Trypsin/EDTA. Add Growth Medium 1B, transfer to a tube, spin down cells, and resuspend in freezing medium (10% DMSO + 90% FBS). Dispense 1 ml of cell aliquots into cryogenic vials. Place vials in an insulated container for slow cooling (-0.5°C to -1°C per min) and freeze at -80°C overnight. The next day, transfer the cells to liquid nitrogen for long term storage.

Functional Validation and Assay Performance

The following assays are designed for a 96-well format. To perform the assay in different tissue culture formats, the cell number and reagent volume should be scaled appropriately.

Materials Required but Not Supplied

- Assay Medium: Thaw Medium 1 (BPS Bioscience #60187)
- Growth Medium 1B (BPS Bioscience #79531)
- Lipofectamine 2000 (Invitrogen #11668027) or other transfection reagent
- DAPT (Selleckchem #S2215): inhibitor of Notch pathway (γ-secretase inhibitor). Prepare 10 mM of stock solution in DMSO.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

- 96-well tissue culture plate or 96-well tissue culture-treated white clear-bottom assay plate (Corning #3610)
- ONE-Step™ Luciferase Assay System (BPS Bioscience #60690)
- Luminometer

Mycoplasma testing

The cell line has been screened using the MycoAlert Mycoplasma Detection kit (Lonza #LT07-318) to confirm the absence of Mycoplasma species.

Inhibition of Notch reporter activity by inhibitor of Notch signaling pathway in Notch CSL Reporter – HEK293 cells

1. Harvest CSL Reporter – HEK293 cells from culture in Growth Medium 1B and seed cells at a density of 35,000 cells per well into white clear-bottom 96-well microplate in 90 μ l Thaw Medium 1. Grow cells overnight at 37°C in a 5% CO₂ incubator. Leave a couple wells empty for use as the control for background luminescence.
2. Using Lipofectamine 2000 or other transfection reagent, transfect cells according to manufacturer's protocol with 50,ng per well of Notch1DeIE expression vector, or a similar genetic construct of interest. Grow cells overnight at 37°C in a 5% CO₂ incubator.
3. Remove medium from the cells, and replace with 90 μ l Thaw Medium 1. Add 10 μ l of diluted inhibitor (DAPT) or vehicle control to the test wells. The final concentration of DMSO in assay medium should be below 0.3%.
Add 100 μ l of Thaw Medium 1 with DMSO to cell-free control wells (for determining background luminescence).
Set up each treatment in at least triplicate.
3. Incubate the plate at 37°C in a 5% CO₂ incubator for 24 - 48 hours.
4. Perform luciferase assay using the ONE-Step™ Luciferase Assay System according to the protocol provided: Mix component B with component A (1:100 ratio), and add 100 μ l master mix per well. Rock at room temperature for ~15 minutes, and measure luminescence using a luminometer. *If using other luciferase reagents from other vendors, follow the manufacturer's assay protocol.*
5. Data Analysis: Obtain the background-subtracted luminescence by subtracting the average background luminescence (cell-free control wells) from the luminescence reading of all wells.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.


To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com

Figure 1. Response of Notch CSL Reporter – HEK293 cells to Notch pathway inhibitor.

A. DAPT Blocked Notch Reporter Activity. CSL-HEK293 cells transfected with Notch1DeIE, Notch2DeIE, Notch3DeIE, Notch4DeIE, or empty vector were treated with 10 μ M DAPT or DMSO control. DAPT inhibited γ -secretase processing of Notch, resulting in decreased luciferase expression.

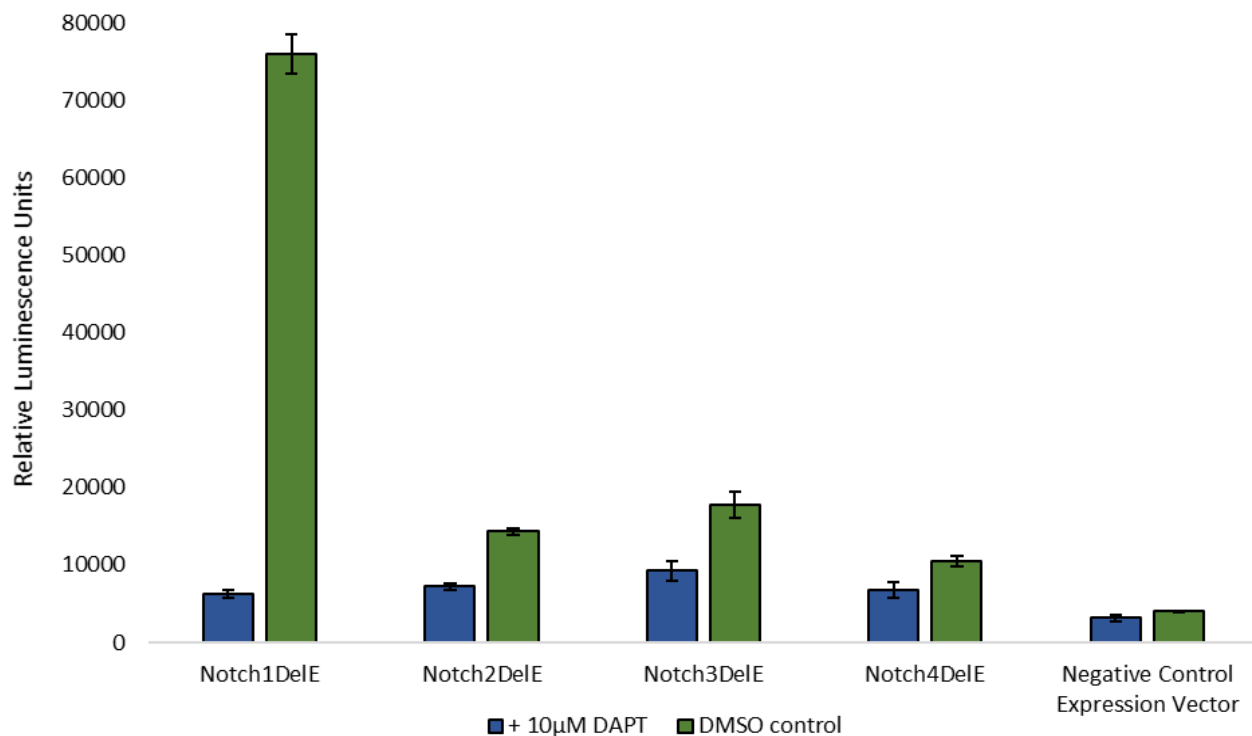

References

Lu, F.M., *et al.* (1996) Constitutively active human Notch1 binds to the transcription factor CBF1 and stimulates transcription through a promoter containing a CBF1-responsive element. *Proc. Natl. Acad. Sci. USA* **93(11)**: 5663-5667.

Kanungo, J., *et al.* (2008) The Notch signaling inhibitor DAPT down-regulates cdk5 activity and modulates the distribution of neuronal cytoskeletal proteins. *J. Neurochem.* **106**: 2236-2248.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

License Disclosure

Purchase of this cell line grants you with a 10-year license to use this cell line in your immediate laboratory, for research use only. This license does not permit you to share, distribute, sell, sublicense, or otherwise make the cell line available for use to other laboratories, departments, research institutions, hospitals, universities, or biotech companies. The license does not permit use of this cell line in humans or for therapeutic or drug use. The license does not permit modification of the cell line in any way. Inappropriate use or distribution of this cell line will result in revocation of the license and result in an immediate cease of sales and distribution of BPS products to your laboratory. BPS does not warrant the suitability of the cell line for any particular use, and does not accept any liability in connection with the handling or use of the cell line. Modifications of this cell line, transfer to another facility, or commercial use of the cells may require a separate license and additional fees; contact sales@bpsbioscience.com for details. Publications using this cell line should reference BPS Bioscience, Inc., San Diego.

Related Products

| <u>Product name</u> | <u>Catalog#</u> | <u>Size</u> |
|---|------------------------|--------------------|
| Notch1 / CSL Reporter – HEK293 Cell line | 60652 | 2 vials |
| Human Notch1 Pathway Reporter Kit | 79503 | 500 rxns |
| Mouse Notch1 Pathway Reporter Kit | 60509 | 500 rxns |
| Mouse NICD Notch1 Pathway Reporter Kit | 79504 | 500 rxns |
| TCF/LEF Reporter Kit (Wnt / β -catenin signaling Pathway) | 60500 | 500 rxns |
| SRE Reporter Kit (MAPK/ERK signaling Pathway) | 60511 | 500 rxns |
| CRE/CREB Kit (cAMP/PKA signaling Pathway) | 60611 | 500 rxns |
| Dual Luciferase (Firefly-Renilla) Assay System | 60683-1 | 10 mL |
| Dual Luciferase (Firefly-Renilla) Assay System | 60683-2 | 100 mL |
| Wnt Signaling Pathway TCF/LEF (Luc) Reporter HEK293 | 60501 | 2 vials |
| Hedgehog Signaling Pathway Gli Reporter-NIH3T3 Cell Line | 60409 | 2 vials |
| ERK Signaling Pathway SRE Reporter-HEK293 Cell Line | 60406 | 2 vials |
| JNK Signaling Pathway AP1 Reporter-HEK293 Cell Line | 60405 | 2 vials |
| JAK/STAT Signaling Pathway ISRE Reporter-HEK293 | 60409 | 2 vials |
| ONE-Step™ Luciferase Assay System | 60690-1 | 10 ml |
| ONE-Step™ Luciferase Assay System | 60690-2 | 100 ml |
| Thaw Medium 1 | 60187-1 | 100 ml |
| Thaw Medium 1 | 60187-2 | 500 ml |
| Growth Medium 1B | 79531 | 500 ml |

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com