



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: info@bpsbioscience.com

**Data sheet**  
**SARS-CoV-2 IgG Detection Kit**  
**(Colorimetric Anti-Spike RBD IgG detection)**  
Catalog #79985  
Size: 96 reactions

**DESCRIPTION:** The SARS-CoV-2 IgG detection kit is designed for qualitative detection of human IgG antibodies in serum collected from individuals suspected of prior infection with the virus that causes COVID-19. This fast and simple ELISA uses the receptor binding domain (RBD) of the SARS-CoV-2 Spike protein (BPS Bioscience #100696) to identify IgG antibodies that indicate a previous infection with SARS-CoV-2. The SARS-CoV-2 IgG (Colorimetric Anti-Spike RBD IgG detection) Detection Kit comes in a convenient 96-well format, with purified SARS-CoV-2 RBD protein and an HRP-conjugated anti-human Fc antibody for 100 binding reactions.

**BACKGROUND:** The pandemic coronavirus disease 2019 (COVID-19) is caused by Severe Acute Respiratory Syndrome Coronavirus 2 (SARS-CoV-2). As a first step of the viral replication strategy, the virus attaches to the host cell surface before entering the cell. The Spike protein recognizes and binds to the Angiotensin-Converting Enzyme 2 (ACE2) receptor found on the surface of type I and II pneumocytes, endothelial cells, and ciliated bronchial epithelial cells. Drugs targeting the interaction between the Spike protein of SARS-CoV-2 and ACE2 may offer some protection against the viral infection.

**COMPONENTS:**

| Catalog # | Component                              | Amount | Storage   |                                    |
|-----------|--|--------|-----------|------------------------------------|
| 100696    | Spike S1 RBD, Avi-His tag (SARS-CoV-2) | 5 µg   | -80°C     | Avoid multiple freeze/thaw cycles! |
|           | anti-human Fc-HRP conjugate (1 mg/ml)  | 5 µl   | +4°C      |                                    |
|           | Colorimetric HRP substrate             | 10 ml  | +4°C      |                                    |
|           | Transparent 96-well microplate         | 1      | Room Temp |                                    |

**APPLICATIONS:** This kit is useful for detection of anti SARS-CoV-2 Spike RBD IgG in human serum

**STABILITY:** Up to 6 months from date of receipt, when stored as recommended.

**REFERENCES:** Long, Q.-X., *et al.* 2020. Antibody responses to SARS-CoV-2 in patients with COVID-19. *Nat. Med.* (in press). <https://doi.org/10.1038/s41591-020-0897-1>

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

**MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:**

PBS (Phosphate buffered saline)  
PBST (Phosphate buffered saline containing 0.05% Tween-20)  
Dry Milk (Fisher #115668 or compatible)  
1N HCl (aqueous)  
Rotating or rocker platform  
UV/Vis spectrophotometer microplate reader capable of reading absorbance at 450 nm

**ASSAY PROTOCOL:**

All samples and controls should be tested in duplicate.

**Coating the plate with SARS-CoV-2 Spike S1 RBD protein:**

- 1) Thaw **Spike S1 RBD protein** on ice. Upon first thaw, briefly spin tube containing **Spike S1 RBD protein** to recover the full contents of the tube. Aliquot into single use aliquots. Immediately store remaining **Spike S1 RBD protein** in aliquots at -80°C. Note: **Spike S1 RBD protein** is very sensitive to freeze/thaw cycles. Avoid multiple freeze/thaw cycles.
- 2) Dilute **Spike S1 RBD protein** to 1 µg/ml in PBS.
- 3) Add 50 µl of diluted **Spike S1 RBD protein** solution to each well and incubate at 4°C overnight.
- 4) After the overnight coating, discard the solution and wash the plate once with 200 µl PBST. Tap plate onto clean paper towels to remove liquid.
- 5) Block wells by adding 200 µl blocking buffer (PBST containing 5% dry milk) to each well. Incubate for 1 hour at room temperature with slow shaking. Remove the blocking solution and wash the plate twice with 200 µl PBST. Tap plate onto clean paper towels to remove liquid.

**Serum sample dilution, IgG capture and detection**

- 1) Dilute the serum sample 1:50 in the blocking buffer (PBST containing 5% dry milk) and add 50 µl diluted sample to the test well.
- 2) For the "Negative Control", add 50 µl of diluted serum that was collected be from a non-Covid19 patient. For the "Blank", add 50 µl of the blocking buffer only. Incubate at room temperature for one hour with slow shaking.
- 3) After one hour, discard the solution and wash the plate three times with 200 µl PBST. Tap plate onto clean paper towels to remove liquid.
- 4) Dilute **anti-human Fc-HRP conjugate** in the blocking buffer (PBST containing 5% dry milk) at 1:2000. Add 50 µl of diluted anti-human Fc-HRP conjugate to each well, and incubate at room temperature for 30 minutes.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)


Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: info@bpsbioscience.com

- 5) After 30 minutes, discard the solution and wash the plate three times with 200  $\mu$ l PBST. Tap plate onto clean paper towels to remove liquid.
- 6) Prepare the stop solution (1N HCl in water). Note: 2N H<sub>2</sub>SO<sub>4</sub> or other compatible acidic solutions can be substituted.
- 7) Add 100  $\mu$ l of the **Colorimetric HRP substrate** to each well and incubate the plate at room temperature until blue color is developed in the positive control well. This usually takes ~ 1 minute but could take less than a minute depending on the IgG level in the sample serum. The optimal incubation time may vary, and should be determined empirically by the user. It is recommended that the reaction be stopped when the 'Negative Control' well is lower than ~ 0.1 absorbance at 450 nm.
- 8) Once blue color has developed in the positive well, add 100  $\mu$ l stop solution prepared in step 6. The positive well should become yellow.
- 9) Read the absorbance at 450 nm using a UV/Vis spectrophotometer microplate reader. "Blank" value is subtracted from all readings. The blank wells should exhibit an absorbance of ~ 0.1 or less at 450 nm as noted above.

**Example of assay results:**

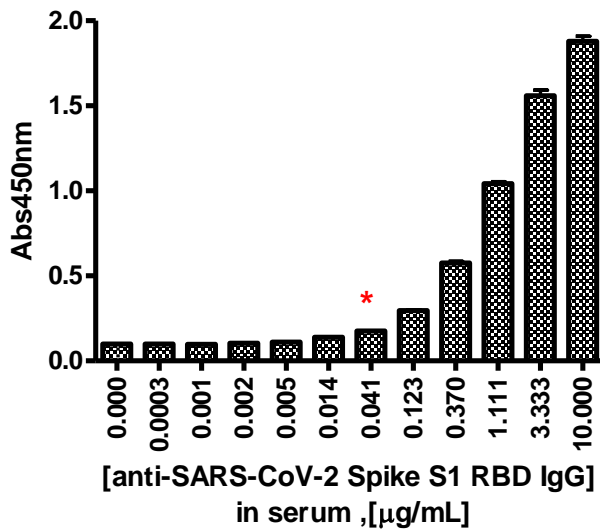

Anti-SARS-CoV-2 Spike S1 RBD IgG in serum was detected by using **SARS-CoV-2 IgG Detection Kit (Colorimetric Anti-Spike RBD IgG detection)**. Positive control human anti-Spike (IgG1) was added to COVID-19 negative human serum at indicated concentrations, and diluted in the blocking buffer as described in the protocol. ~ 50 ng/ml of IgG in serum can be detected with confidence (absorbance should be at least 1.5 fold higher than the negative serum; e.g. Abs<sub>Negative Control</sub> = 0.1, Abs<sub>41ng/ml</sub> = 0.17, Abs<sub>123ng/ml</sub> = 0.3)

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
San Diego, CA 92121  
Tel: 1.858.202.1401  
Fax: 1.858.481.8694  
Email: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

**RELATED PRODUCTS:**

| <b><u>Product Name</u></b>                                    | <b><u>Catalog#</u></b> | <b><u>Size</u></b> |
|---|------------------------|--------------------|
| SARS-CoV-2 Spike:ACE2 Inhibitor Screening Assay Kit           | 79931                  | 96 reactions       |
| ACE2:SARS-CoV-2 Spike Inhibitor Screening Assay Kit           | 79936                  | 96 reactions       |
| ACE2:SARS-CoV-2 Spike S1-Biotin Inhibitor Screening Assay Kit | 79945                  | 96 reactions       |
| SARS-CoV-2 Spike S1-Biotin:ACE2 TR-FRET Assay Kit             | 79949                  | 96 reactions       |
| Spike S1, Fc Fusion, Avi-tag (SARS-CoV-2)                     | 100678                 | 100 µg/1 mg        |
| Spike S1, Fc fusion, Avi-tag, Biotin-Labeled (SARS-CoV-2)     | 100679                 | 25 µg/50 µg        |
| Spike S1 RBD, His-tag (SARS-CoV-2)                            | 100687                 | 50 µg/100 µg       |
| Spike S1, Fc fusion (SARS-CoV-2)                              | 100688                 | 20 µg/50 µg        |
| Spike S1 RBD, Fc fusion (SARS-CoV-2)                          | 100699                 | 50 µg/100 µg       |
| ACE2 Inhibitor Screening Assay Kit                            | 79923                  | 96 reactions       |
| ACE2, His-tag   | 11003                  | 20 µg/100 µg       |
| ACE2, His-Avi-Tag, Biotin-labeled HiP™                        | 100665                 | 20 µg/50 µg        |
| ACE2, Fc Fusion (Monkey)                                      | 100701                 | 50 µg/1 mg         |
| ACE2, His-tag (Monkey)  | 100702                 | 50 µg/1 mg         |

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Court W, Ste B  
 San Diego, CA 92121  
 Tel: 1.858.202.1401  
 Fax: 1.858.481.8694  
 Email: info@bpsbioscience.com

**TROUBLESHOOTING GUIDE:**

| Problem   | Possible cause   | Solution   |
|---|--|--|
| Colorimetric signal of positive control reaction is weak    | Spike trimer or serum sample has lost activity           | Increase time of incubation with serum. Increase serum concentration. Proteins lose activity upon repeated freeze/thaw cycles. Store S1 trimer and anti-Fc HRP conjugate in single-use aliquots. |
|   | Incorrect settings on instruments                        | Refer to instrument instructions for settings to increase sensitivity of light detection.  |
|   | Colorimetric HRP substrate was not incubated long enough | Increase the amount of time that the colorimetric HRP substrate is incubated in the wells. Avoid azides.   |
| Colorimetric signal is erratic or varies widely among wells | Inaccurate pipetting/technique                           | Run duplicates of all reactions. Use a multichannel pipettor. Use master mixes to minimize errors.   |
|   | Bubbles in wells   | Pipette slowly to avoid bubble formation. Tap plate lightly to disperse bubbles; be careful not to splash between wells.   |
|   | Signal is out of range of detection (too high)           | Decrease the amount of time that the colorimetric HRP substrate is incubated in the wells  |
| Background (signal to noise ratio) is high                  | Insufficient washes                                      | Increase number of washes. Increase wash volume.   |
|   | Sample solvent is inhibiting the enzyme                  | Run negative control assay including solvent. Maintain DMSO level at <1% Increase time of enzyme incubation.   |
|   | Results are outside the linear range of the assay        | Use different concentrations of proteins to create a standard curve  |

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)