



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet ***Renin Assay Kit*** **Catalog # 80211**

DESCRIPTION: Renin is a protease that is synthesized by the kidneys. Renin catalyzes the hydrolysis of angiotensinogen to angiotensin I, which is converted to angiotensin II, a highly potent vasoconstrictor. Renin inhibitors have been used therapeutically in the treatment of hypertension (high blood pressure). The *Renin Assay Kit* is designed to measure Renin activity for screening and profiling applications. It comes in a convenient 96-well format, with purified, activated renin, renin substrate, and assay buffer for 100 enzyme reactions. Aliskiren Hydrochloride is also included as a control for renin inhibition. The key to the Renin Assay Kit is the specific, fluorogenic substrate. Using this kit, only one simple step on a microtiter plate is required for renin reactions. The fluorometric substrate is incubated with a sample containing renin enzyme to produce a fluorophore that can then be measured using a fluorescence reader.

COMPONENTS:

Catalog #	Component	Amount	Storage	
80200	Activated Renin	2 µg	-80°C	(Avoid freeze/thaw cycles!)
	Renin Assay Buffer	25 ml	-20°C	
	Renin Substrate	30 µl @ 1 mM	-80°C	
27303	Aliskiren Hydrochloride	20 µl @ 100µM in DMSO	-20°C	
	Black 96-well plate		RT	

MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

Fluorescent microplate reader

APPLICATIONS: Great for studying enzyme kinetics, screening small molecular inhibitors for drug discovery and HTS applications.

STABILITY: 6 months from date of receipt when stored as directed.

REFERENCE(S):

1. Paschalidou, K. *et al.*, *Biochem. J.* 2004; **382**: 1031.
2. Gossas, T., *et al.*, *Naunyn Schmiedebergs Arch Pharmacol.* 2012; **385(2)**:219-24.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate.

Step 1:

- 1) Dilute the Renin Substrate 1:100 using the Renin Assay Buffer. For example, add 30 μ l of Renin Substrate to 2,970 μ l of Renin Assay Buffer.
- 2) Add 25 μ l of diluted Renin Substrate to each well designated as "Positive Control", "Test Inhibitor", "Inhibitor Control", and "Blank".
- 3) Dilute Aliskiren HCl (100 μ M stock) 10-fold with Renin Assay Buffer to make a 10 μ M solution. Make a series of dilutions of the 10 μ M solution with 10% DMSO (prepared in Renin Assay Buffer) for IC₅₀ assay. You can use the same method to prepare the test inhibitor.
- 4) Add 5 μ l of the prepared test inhibitor solution to each well designated "Test Inhibitor". For the "Positive Control" and "Blank", add 5 μ l of the same solution without the inhibitor (Inhibitor buffer). For the "Inhibitor Control", add 5 μ l of the Aliskiren HCl solution.
- 5) Thaw the human recombinant Renin on ice. Upon the first thaw, briefly spin the tube containing the enzyme to recover the full contents of the tube. Aliquot the Renin enzyme into single-use aliquots. Store the remaining undiluted enzyme in aliquots at - 80°C immediately. *Note: Renin is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*

	Blank	Positive Control	Inhibitor Control	Test Inhibitor
Renin Substrate (10 μ M)	25 μ l	25 μ l	25 μ l	25 μ l
Test Inhibitor				5 μ l
Renin Assay Buffer	20 μ l			
Aliskiren HCl			5 μ l	
Inhibitor Buffer (no inhibitor)	5 μ l	5 μ l		
Human recombinant Renin		20 μ l	20 μ l	20 μ l
Total	50 μl	50 μl	50 μl	50 μl

- 6) Dilute the human recombinant Renin to 0.1 ng/ μ l using the Renin Assay Buffer. Add 20 μ l of diluted Renin to each well designated "Positive Control", "Test Inhibitor", and "Inhibitor Control". For the "Blank", add 20 μ l of Renin Assay Buffer.

- 7) Incubate the plate at room temperature for 30 minutes.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com




6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.202.1401
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Step 2:

Read the sample in a microtiter-plate fluorimeter that is capable of excitation at wavelengths 490 nm and detection of emitted light at 520 nm. The "Blank" value is subtracted from all other values.

Example of Assay Results:

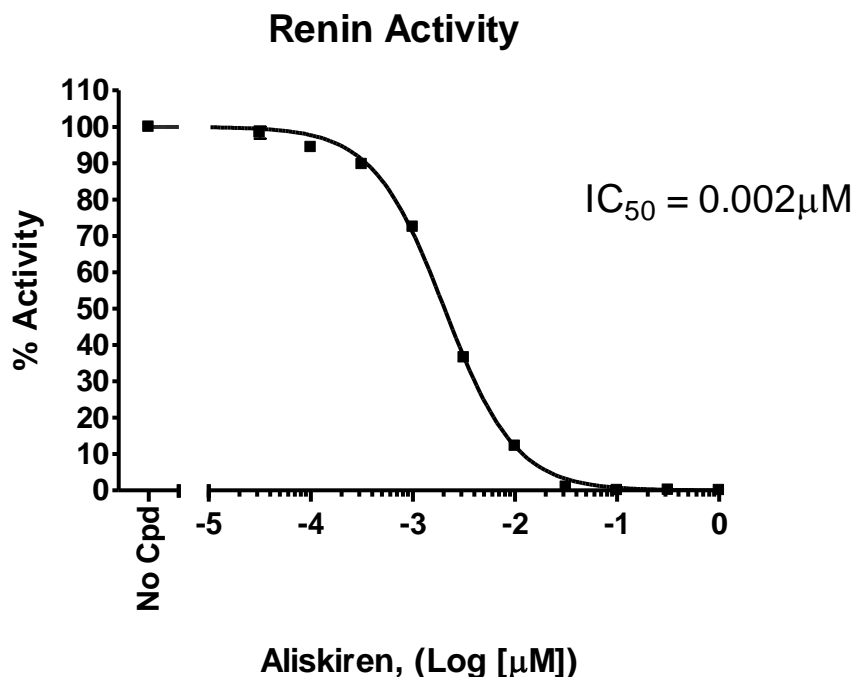

Inhibition of Activated Renin (BPS Bioscience, #80200) by Aliskiren Hydrochloride (BPS Bioscience, #27303), measured using the *Renin Assay Kit* (BPS Bioscience, #80211). Fluorescence intensity was measured using a Tecan fluorescent microplate reader. *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com*

RELATED PRODUCTS

<u>Product Name</u>	<u>Catalog#</u>	<u>Size</u>
Activated Renin	80200	100 µg
Aliskiren Hydrochloride	27303	1 mg

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.
To place your order, please contact us by Phone **1.858.202.1401** Fax **1.858.481.8694**
Or you can Email us at: info@bpsbioscience.com
Please visit our website at: www.bpsbioscience.com