



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

## Description

The Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 Cell Line is a HEK293 cell line stably expressing the full-length human muscarinic acetylcholine receptor M1 (CHRM1/M1R/M1; accession number: AF498915) and a firefly luciferase reporter under the control of NFAT (nuclear factor of activated T cells) response element.

This cell line has been validated for its responses to the muscarinic acetylcholine receptor agonists Oxotremorine M and Carbamoylcholine chloride (carbachol). Luciferase activity increases in proportion to mAChR activation.

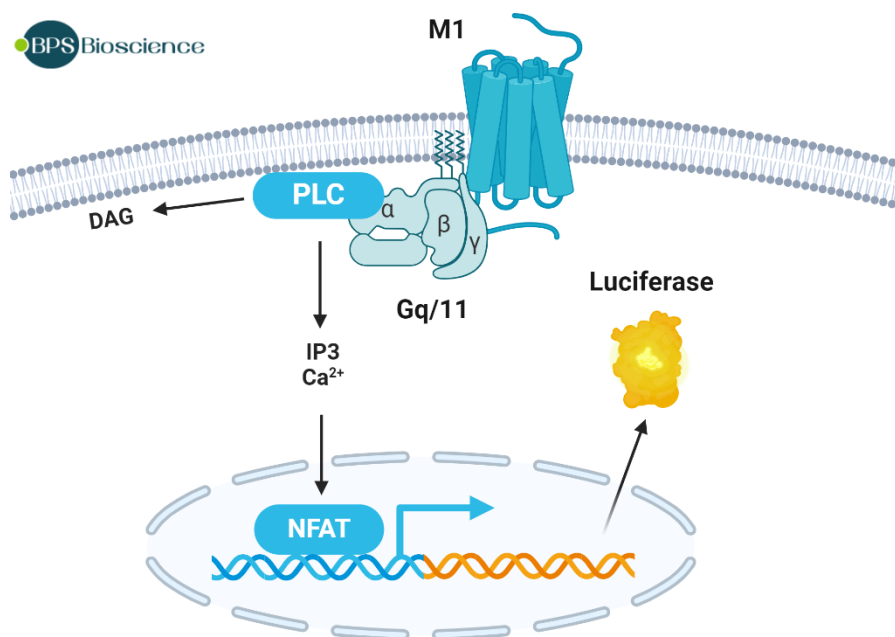


Figure 1: Diagram illustrating the mechanism of action of Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 Cell Line.

## Background

Muscarinic acetylcholine receptors (mAChRs) are relatively abundant and mediate many of the diverse actions of acetylcholine in the CNS (central nervous system), as well as throughout non-nervous tissues innervated by the parasympathetic nervous system. They are involved in regulating a large number of cognitive, behavioral, motor, and autonomic functions. There are five muscarinic receptor subtypes referred to as M1, M2, M3, M4 and M5 which belong to the GPCR (G-protein-coupled receptor) superfamily. M1, M3 and M5 receptor subtypes couple efficiently through Gq/11, to activate phospholipase C (PLC), which initiates the phosphatidylinositol (IP3) turnover response and leads to the release of intracellular calcium ions (Ca<sup>2+</sup>). The M2 and M4 receptor subtypes preferentially couple to Gi/o G-proteins and inhibit adenylyl cyclase (AC) activity, which leads to a decrease in the level of cAMP. M2 and M4 receptors also activate G protein-coupled potassium channel by  $\beta\gamma$ -dimer dissociated from the active Gi/o G-protein, which leads to hyperpolarization of the plasma membrane of excitable cells. Disruption of muscarinic signaling contributes to several pathophysiological conditions in the CNS and in the periphery. Thus, muscarinic agonists have a wide therapeutic potential in the treatment of neurodegenerative and neuropsychiatric disorders and conditions, e.g. Alzheimer's disease, schizophrenia, pain, ischemia or heart failure. To target these diseases, selective modulation of individual muscarinic receptor subtypes is necessary to avoid undesired side effects. The orthosteric binding site of all subtypes of muscarinic receptors is virtually the

same and no affinity-based selective agonists of muscarinic receptors have been discovered so far. The development of biased agonists can be the right approach to achieve selective modulation of individual subtypes of muscarinic receptors.

### Application

- Screen for activators or inhibitors in M1 receptor-related research and drug discovery.
- Counter screen functionally selective agonists or biased agonists for M3 and M5 receptors.
- Characterize mAChR M1 antibodies/antagonists and agonists in binding assays.

### Materials Provided

| Components              | Format  |
|-------------------------|---|
| 2 vials of frozen cells | Each vial contains >1 x 10 <sup>6</sup> cells in 1 ml of Cell Freezing Medium (BPS Bioscience #79796) |

### Host Cell

HEK293, epithelial-like cells, adherent

### Mycoplasma Testing

The cell line has been screened to confirm the absence of Mycoplasma species.

### Materials Required but Not Supplied



These materials are not supplied with this cell line but are necessary for cell culture and cellular assays. BPS Bioscience reagents systems are validated and optimized for use with this cell line and are highly recommended for best results. Media components are provided in the Media Formulations section.

#### Materials Required for Cell Culture

| Name             | Ordering Information                  |
|------------------|---------------------------------------|
| Thaw Medium 1    | <a href="#">BPS Bioscience #60187</a> |
| Growth Medium 1G | <a href="#">BPS Bioscience #79544</a> |

#### Materials Required for Cellular Assays

| Name                              | Ordering Information                  |
|-----------------------------------|---------------------------------------|
| Thaw Medium 1                     | <a href="#">BPS Bioscience #60187</a> |
| Growth Medium 1G                  | <a href="#">BPS Bioscience #79544</a> |
| MEM medium                        | Hyclone #SH30024.01                   |
| Carbamoylcholine chloride         | TOCRIS #2810                          |
| Oxotremorine M                    | TOCRIS #1067                          |
| ONE-Step™ Luciferase Assay System | <a href="#">BPS Bioscience #60690</a> |
| Luminometer                       |                                       |

### Storage Conditions



Cells are shipped in dry ice and should immediately be thawed or stored in liquid nitrogen upon receipt. Do not use a -80°C freezer for long term storage. Contact technical support at support@bpsbioscience.com if the cells are not frozen in dry ice upon arrival.

### Media Formulations

For best results, the use of validated and optimized media from BPS Bioscience is *highly recommended*. Other preparations or formulations of media may result in suboptimal performance.



Note: Thaw Media do *not* contain selective antibiotics. However, Growth Media *do* contain selective antibiotics, which are used to maintain selective pressure on the cell population expressing the gene of interest.

Cells should be grown at 37°C with 5% CO<sub>2</sub>. BPS Bioscience's cell lines are stable for at least 10 passages when grown under proper conditions.

#### Media Required for Cell Culture

*Thaw Medium 1 (BPS Bioscience #60187):*

MEM medium supplemented with 10% FBS, 1% non-essential amino acids, 1 mM Na pyruvate, 1% Penicillin/Streptomycin.

*Growth Medium 1G (BPS Bioscience #79544):*

MEM medium supplemented with 10% FBS, 1% non-essential amino acids, 1 mM Na pyruvate, 1% Penicillin/Streptomycin plus 400 µg/ml of Geneticin and 50 µg/ml of Hygromycin B.

### Functional Validation

#### A. Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 Cell Line Agonist Response Assay

- The following assay was designed for a 96-well format. To perform the assay in different tissue culture formats, the cell number and reagent volume should be scaled appropriately.
  - All conditions should be performed in triplicate.
  - Assay should include "Stimulated", "Unstimulated" and "Background Luminescence" conditions.
1. Harvest Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 cells from culture in Growth Medium 1G and seed cells at a density of ~35,000 cells per well into a white clear-bottom 96-well plate in 90 µl of Thaw Medium 1. Leave a few empty wells to use as "Background Luminescence" control.
  2. Incubate the plate at 37°C in a CO<sub>2</sub> incubator for 16 to 24 hours.
  3. Prepare a three-fold serial dilution at 10x the final concentrations of Oxotremorine M and Carbamoylcholine chloride agonists separately in Thaw Medium 1 (10 µl/well). For an EC<sub>50</sub> dose curve, we recommend a range of approximately 0.002 to 50 µM, final concentration.

4. Add 10  $\mu$ l diluted Oxotremorine M or Carbamoylcholine to the “Stimulated” wells.
5. Add 10  $\mu$ l Thaw Medium 1 to “Unstimulated” wells.
6. Add 100  $\mu$ l Thaw Medium 1 to “Background Luminescence” wells (for determining background luminescence).
7. Incubate the plate at 37°C in a CO<sub>2</sub> incubator for about 5 hours.
8. Add 100  $\mu$ l of ONE-Step™ Luciferase reagent per well and rock at Room Temperature (RT) for 15-30 minutes.
9. Measure luminescence using a luminometer.
10. Data Analysis: Subtract the average background luminescence (cell-free control wells) from the luminescence reading of all wells. The fold induction of NFAT luciferase reporter expression is the background-subtracted luminescence of treated well divided by the average background-subtracted luminescence of control wells.

$$\text{Fold induction} = \frac{(\text{luminescence of stimulated cells} - \text{avg background})}{(\text{avg luminescence of unstimulated cells} - \text{avg background})}$$

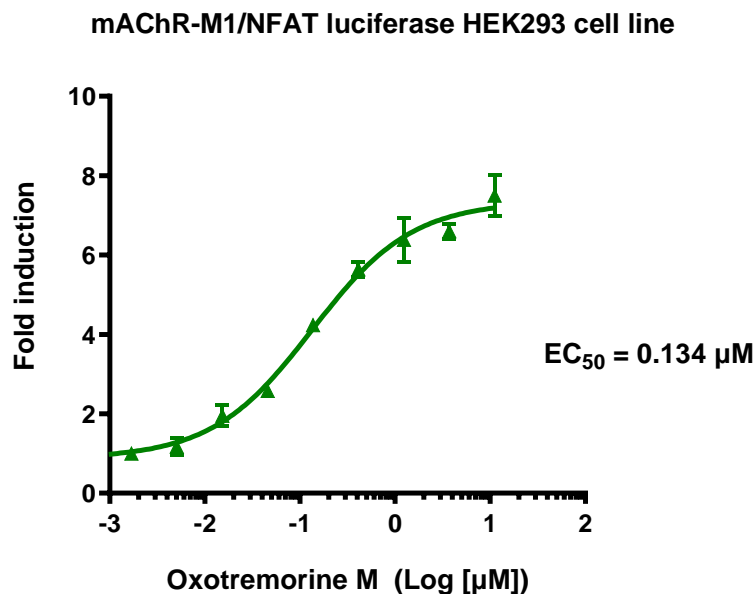


Figure 2. Dose response of Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 Cell Line to Oxotremorine M.

Oxotremorine M agonist was diluted and added to Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 cells for about 5 hours at 37°C in a cell culture incubator. Luciferase activity was measured using ONE-Step™ Luciferase Assay System (#60690).

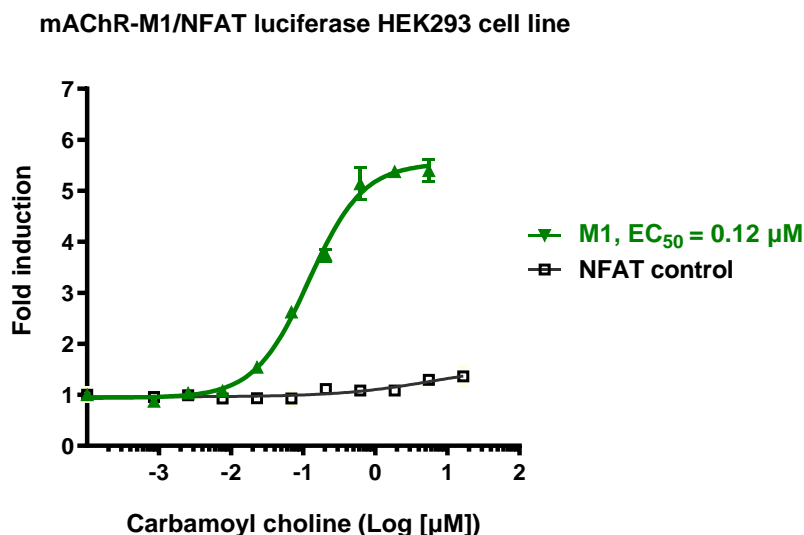


Figure 3. Dose response of Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 Cell Line to carbamoyl choline.

Carbamoylcholine chloride agonist was diluted and added to Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 cells for about 5 hours at 37°C in a cell culture incubator. Luciferase activity was measured using ONE-Step™ Luciferase Assay System (#60690).

Data shown is representative. For lot-specific information, please contact BPS Bioscience, Inc. at [support@bpsbioscience.com](mailto:support@bpsbioscience.com).

### Sequence

Human mAChR M1(CHRM1) sequence (accession number: AF498915)

```
MNTSAPPAVSPNITVLAPGKGPWQVAFIGITTGLLSLATVTGNLLVLISFKVNTLKTVNNNYFLLSLACADLIIGTFSMNLTYTLLM
GHWALGTLACDLWLALDYVASNASVMNLLISFDYFVTRPLSYRAKRTPRRAALMIGLAWLVSFVLWAPAILFWQYLVGERT
VLAGQCYIQFLSQPIITFGTAMAAFYLPVTVMCTLYWRIYRETENRARELAALQGSETPGKGGGSSSSSERSQPGAEGSPETPPGR
CCRCCRAPRLLQAYSWKEEEEEDEGSMESLTSSEGEPEGSEVVIKMPMVDPEAQPTKQPPRSSPNTVKRPTKKGRDRAGKQK
PRGKEQLAKRKTFSVLVKEKKAARTLSAILLAFILTWTPYNIMVLVSTFCKDCVPETLWELGYWLCYVNSTINPMCYALCNKAFRDFT
RLLLLCRWDKRRWRKIPKRPGSVHRTPSRQC
```

### License Disclosure

Visit [bpsbioscience.com/license](https://bpsbioscience.com/license) for the label license and other key information about this product.

### Troubleshooting Guide

Muscarinic Acetylcholine Receptor (mAChR) M1/NFAT Luciferase Reporter HEK293 Cell Line

Visit [bpsbioscience.com/cell-line-faq](https://bpsbioscience.com/cell-line-faq) for detailed troubleshooting instructions. For all further questions, please email [support@bpsbioscience.com](mailto:support@bpsbioscience.com).

**Related Products**

| <i>Products</i>  | <i>Catalog #</i> | <i>Size</i> |
|--|------------------|-------------|
| NFAT Reporter– HEK293 Cell Line (PKC/Ca <sup>2+</sup> Pathway)                         | 79298            | 2 vials     |
| Muscarinic Acetylcholine Receptor (mAChR) M5/NFAT Luciferase Reporter HEK293 Cell Line | 82729            | 2 vials     |
| Muscarinic Acetylcholine Receptor (mAChR) M3/NFAT Luciferase Reporter HEK293 Cell Line | 82728            | 2 vials     |

*Version 011625*