



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Product Information

Resazurin Cell Viability Assay Kit

Kit Contents

30025-1 (2500 assays): 25 mL resazurin solution
 30025-2 (10,000 assays): 100 mL resazurin solution

Storage and Handling

Store at 4°C, protected from light. Product is stable for at least 6 months from date of receipt when stored as recommended.

Spectral Properties

After reduction of resazurin to resorufin (at neutral pH):
 Absorbance/excitation: 571 nm
 Emission: 585 nm

Product Description

Resazurin Fluorometric Cell Viability Assay Kit offers a simple, rapid, reliable, sensitive, safe and cost-effective measurement of cell viability. The assay is homogenous and requires no cell lysis or washing. Biotium's resazurin assay performs at least as well as other commercial resazurin-based cell proliferation assay kits with the trademark name AlamarBlue®.

Cell growth creates a reduced environment while inhibition of growth maintains an oxidized environment. In response to chemical reduction of growth medium resulting from cell growth, resazurin (which is purple and non-fluorescent) is reduced to form the red fluorescent dye resorufin (1-3). Reduction of resazurin can be monitored by measuring fluorescence or absorbance. The fluorescent and colorimetric signal generated from the assay is proportional to the number of living cells in the sample. The resazurin assay is as sensitive as [³H] thymidine assay for detecting cell proliferation (1). Depending on the cell type, the resazurin assay can be used to detect as few as 40 cells with reproducible and sensitive signal.

References

- Ahmed SA, Gogal RM Jr, Walsh JE. (1994). A new rapid and simple non-radioactive assay to monitor and determine the proliferation of lymphocytes: an alternative to [³H]thymidine incorporation assay. *J Immunol Methods*. 170(2):211-24.
- Shahan TA, Siegel PD, Sorenson WG, Kuschner WG, Lewis DM. (1994). A sensitive new bioassay for tumor necrosis factor. *J Immunol Methods*. 175(2):181-7.
- Nociari MM, Shalev A, Benias P, Russo C. (1998). A novel one-step, highly sensitive fluorometric assay to evaluate cell-mediated cytotoxicity. *J Immunol Methods*. 213(2):157-67.

Please visit our website at www.biotium.com for information on our life science research products, including cell viability, proliferation, and apoptosis assays, non-mutagenic GelRed™ & GelGreen™ nucleic acid gel stains, environmentally friendly EvaGreen® qPCR master mixes, fluorescent CF™ dye antibody conjugates and antibody labeling kits, and more probes and kits for life science research.

AlamarBlue is a registered trademark of Trek Diagnostics Systems, Inc. Materials from Biotium are sold for research use only, and are not intended for food, drug, household, or cosmetic use.

Resazurin Cell Viability Assay Kit
 PSF006

Assay Protocol

Note: Resorufin can be further reduced to hydroresorufin (colorless and non-fluorescent). At higher cell densities or with prolonged development times, the assay signal can initially increase, then decrease after all resazurin is converted into resorufin, which then begins to be further reduced to hydroresorufin. Therefore, it is important to conduct a cell number titration (standard curve) to identify the optimal plating density and development time that generates signal that increases proportionally with cell number.

- Plate cells in 96-well tissue culture plates in 100 uL/well. For a standard curve, plate a series of cell dilutions in the range of 40-20,000 cells per well for adherent cells, and 2,000 to 500,000 cells per well for suspension cells. For fluorescence-based detection, include a well with 100 uL of cell culture medium without cells to use as a background control.
- After cells have reached the desired density, add 10 uL resazurin solution to the medium in each well, and mix thoroughly.
- Incubate the plate for between 1 hour and 24 hours at 37°C.

Note: Signal from the same plate can be read at multiple time points to determine the optimal incubation time for your cell type and density.

- For colorimetric detection, measure absorbance at 570 nm and 600 nm using an absorbance microplate reader. For fluorescence-based detection, measure fluorescence with excitation/emission at 570/585 nm using a fluorescence microplate reader.

Note: Fluorescence-based detection is more sensitive and has broader dynamic range than colorimetric detection.

Note: The excitation and emission spectra of resorufin are fairly broad, excitation filters between 530–570 nm and emission filters between 580-620 nm can be used.

- For the colorimetric detection method, subtract background absorbance at 600 nm from resorufin absorbance at 570 nm. For fluorescence-based detection, subtract fluorescence at 585 nm from the background control (culture medium without cells) from each cell sample.
- Plot cell plating density vs. background-subtracted absorbance or fluorescence for your cell number titration to determine the optimal assay conditions for your cell line.

Related Products

| Cat. no. | Product |
|-------------|---|
| 30026 | Calcein AM Cell Viability Assay Kit |
| 30006 | MTT Cell Viability Assay Kit |
| 30007 | XTT Cell Viability Assay Kit |
| 30020 | ATP-Glo™ Bioluminometric Cell Viability Assay Kit |
| 30068 | ViaFluor™ 405-SE Cell Proliferation Kit |
| 30066 | Apoptotic, Necrotic, and Healthy Cells Quantitation Kit Plus |
| 30060 | CF™488A Annexin V and 7-AAD Apoptosis Kit |
| 30061 | CF™488A Annexin V and PI Apoptosis Kit |
| 32002-32009 | Live-or-Dye™ Fixable Viability Staining Kits (choose from 8 fluorescent dye colors) |
| 30029 | NucView™ 488 Caspase-3 Assay Kit for live cells |
| 30067 | Dual Apoptosis Assay with NucView™ 488 and CF™594 Annexin V |
| 30072 | NucView™ 488 and RedDot™ 2 Apoptosis and Necrosis Kit |
| 10405 | NucView™ 405 Caspase-3 Substrate |
| 30019 | MCB Glutathione Detection Kit |