



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

CGRP Affinity Sorbent

ORDERING INFORMATION

Catalogue Number: A19482

Stability: ≥ 2 years

Storage: +4°C

Shipping: ice pack

• **PRODUCT DESCRIPTION**

The following is the purification protocol for CGRP sorbent provided by SPI-BIO.


- A. Pour the gel in an empty column (around 10 x 1 cm).
- B. Let the storage solution elute entirely (the gel should become dry but not for a long time).
- C. Add around 20 mL of plasma on top of the column with the outlet of the column closed.
- D. Let it incubate overnight at +4°C with gentle agitation.
- E. Elute the plasma that is now free of CGRP and ready to be used to dilute standards and samples in the assay.
- F. So as to regenerate the gel (eliminate the bound CGRP), wash the column with around 20 mL of 0.1 M HCl pH 1.5 and then with around 50 mL of 0.1 M Phosphate buffer pH 7.4.
- G. You can then store the sorbent in 3 to 4 mL of EIA buffer (buffer of the kit) at +4°C till the next use. It is stable at +4°C for many months.

Related Products

1. Human CGRP EIA kit ref. #A05481.
2. Rat CGRP EIA kit ref. #A05482.

**For research laboratory
use only.**

**Not for human
diagnostic use.**

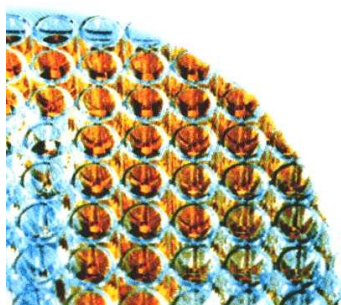

November 05

Should you wish further information, do not hesitate to contact us.