

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten! See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere Liefer- und Versandbedingungen

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

linkedin.com/company/szaboscandic in



Product Information

Cell-Based Assay Soluble Epoxide Hydrolase Substrate

Item No. 600095

Contents: This vial contains 100 µl of 10 mM Epoxy Fluor 7 in DMSO

Stability: ≥1 year at -20°C

Laboratory Procedures

Prior to assaying, dilute 10 µl of Soluble Epoxide Hydrolase (sEH) Substrate with 10 ml of the diluted Assay Buffer. Add 100 µl of this substrate solution to each well except the wells containing 6-methoxy-2-Naphthaldehyde Standards.

CAUTIONS

For laboratory research use only. This reagent is not for human or veterinary use. This product has been tested and formulated to work exclusively with the sEH Cell-Based Assay Kit from Cayman (Item No. 600090). This product may not perform as described if used with other assay reagents or protocols.

Related Product

Soluble Epoxide Hydrolase Cell-Based Assay Kit - Item No. 600090

WARNING: This product is for laboratory research only: not for administration to humans. Not for human or veterinary DIAGNOSTIC OR THERAPEUTIC USE.

MATERIAL SAFETY DATA

This material should be considered hazardous until information to the contrary becomes available. Do not ingest, swallow, or inhale. Do not get in eyes, on skin, or on clothing. Wash thoroughly after handling. This information contains some, but not all, of the information required for the safe and proper use of this material. Before use, the user must review the complete Material Safety Data Sheet, which has been sent via email to your institution.

Cayman Chemical Company makes **no warranty or guarantee** of any kind, whether written or oral, expressed or implied, including without limitation, any warranty of fitness for a particular purpose, suitability and merchantability, which extends beyond the description of the chemicals hereof. Cayman **warrants only** to the original customer that the material will <u>meet our specifications</u>

tools not linest out specifications.

Said refund or replacement is conditioned on Buyer giving written notice to Cayman within thirty (30) days after arrival of the material at its destination. Failure of Buyer to give said notice within thirty (30) days shall constitute a waiver by Buyer of all claims hereunder with respect to said material.

For further details, please refer to our Warranty and Limitation of Remedy located on our website and in our catalog.

Copyright Cayman Chemical Company, 01/26/2011

www.caymanchem.com

custserv@caymanchem.com

Cayman Chemical

Mailing address 1180 E. Ellsworth Road Ann Arbor, MI 48108 USA **Phone**

(800) 364-9897

(734) 971-3335

(734) 971-3640

WARRANTY AND LIMITATION OF REMEDY

art the time of delivery.

Cayman will carry out its delivery obligations with due care and skill. Thus, in no event will Cayman have any obligation or liability, whether in tort (including negligence) or in contract, for any direct, incidental or consequential damages, even if Cayman is informed about their possible existence.

This limitation of liability does not apply in the case of intentional acts or negligence of Cayman, districtors or its employees.

Buyer's exclusive remedy and Cayman's sole liability hereunder shall be limited to a refund of the purchase price, or at Cayman's option, the replacement, at no cost to Buyer, of all material that does not meet our specifications.

Said refund or replacement is conditioned on Buyer giving written notice to Cayman within thirty (30) days after a trival of the material at its destination. Failure of Buyer to give said notice within