

# Produktinformation



Forschungsprodukte & Biochemikalien
Zellkultur & Verbrauchsmaterial
Diagnostik & molekulare Diagnostik
Laborgeräte & Service

Weitere Information auf den folgenden Seiten! See the following pages for more information!



Lieferung & Zahlungsart siehe unsere Liefer- und Versandbedingungen

# Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

# SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien T. +43(0)1 489 3961-0 F. +43(0)1 489 3961-7 <u>mail@szabo-scandic.com</u> www.szabo-scandic.com

# **Product Information**



### Cell-Based Assay Monochlorobimane

Item No. 600361

**Contents:** This vial contains 100 µl of monochlorobimane in DMSO. Stability: ≥1 year at -20°C

#### Laboratory Procedures

To prepare a 1 ml of Substrate Solution, add 50 µl of the Cell-Based Assay Monochlorobimane to 1 ml of Cell-Based Assay Buffer (prepared from Cell-Based Assay Buffer Tablet, Item No. 10009322). If you are not using all the Monochlorobimane Substrate at one time, we recommend that you make small aliquots for long-term storage. The Monochlorobimane Substrate is stable for six months at -20°C.

#### CAUTIONS

For laboratory research use only. This reagent is not for human or veterinary use. This product has been tested and formulated to work exclusively with the Glutathione Cell-Based Detection Kit (Blue Fluorescence) from Cayman (Item No. 600360). This product may not perform as described if used with other assay reagents or protocols.

## **Related Product**

DIAGNOSTIC OR THERAPEUTIC USE.

Sheet, which has been sent via email to your institution. WARRANTY AND LIMITATION OF REMEDY

MATERIAL SAFETY DATA

at the time of delivery.

Glutathione Cell-Based Detection Kit (Blue Fluorescence) - Item No. 600360

#### **Cayman Chemical**

#### **Mailing address**

1180 E. Ellsworth Road Ann Arbor, MI 48108 USA

Phone (800) 364-9897 (734) 971-3335

Fax (734) 971-3640

#### E-Mail

custserv@caymanchem.com

#### Web www.caymanchem.com

any direct, incidental or consequential damages, even if Cayman is informed about their possible existence. This limitation of liability does not apply in the case of intentional acts or negligence of Cayman, its directors or its employees. Buyer's **exclusive remedy** and Cayman's sole liability hereunder shall be limited to a <u>refund</u> of the purchase price, or at Cayman's option, the <u>replacement</u>, at no cost to Buyer, of all material that does not meet our specifications.

Cayman Chemical Company makes no warranty or guarantee of any kind, whether written or oral, expressed or implied, including without limitation, any warranty of fitness for a particular pose, suitability and merchantability, which extends beyond the description of the chemicals hereof. Cayman warrants only to the original customer that the material will meet our specifications

2 unic on tearers, and the second second

WARNING: This product is for laboratory research only: not for administration to humans. Not for human or veterinary

This material should be considered hazardous until information to the contrary becomes available. Do not ingest, swallow, or inhale. Do not get in eyes, on skin, or on clothing. Wash thoroughly after handling. This information contains some, but not all, of the information required for the safe and proper use of this material. Before use, the user must review the complete Material Safety Data

a does not meet our specifications. Said refund or replacement is conditioned on Buyer giving written notice to Cayman within thirty (30) days after arrival of the material at its destination. Failure of Buyer to give said notice within thirty (30) days shall constitute a waiver by Buyer of all claims hereunder with respect to said material. For further details, please refer to our Warranty and Limitation of Remedy located on our website and in our catalog. Copyright Cayman Chemical Company, 01/03/2011