



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Technically
Speaking

CEDARLANE[®] 
www.cedarlanelabs.com

Conveniently Delivering You Today's Innovations
for the Science of Tomorrow™

Human IgG Sensitized Glutaraldehyde Stabilized Sheep RBCs

CLD68LS-1

Lot: 1292A

DESCRIPTION:

Cedarlane's **CLD68LS-1** are Human IgG Sensitized Glutaraldehyde-stabilized freshly prepared sheep Red Blood Cells. Glutaraldehyde stabilized red blood cells are produced from freshly washed sheep RBCs by brief exposure to a glutaraldehyde-buffer solution and exhaustive washing in saline. This treatment largely preserves the native antigenicity of the erythrocytes while rendering the cells generally resistant to lysis by osmotic shock, freeze-thawing or immune hemolysis. This reagent can be used directly in hemagglutination procedures, or may be coupled with various proteins for hemagglutination testing.

ATTACHMENT OF HUMAN IGG TO RBC'S:

The above lot of sensitized sheep RBCs clump when mixed with rabbit anti-human IgG, and incubated at 37°C. No clumping of the above lot of sensitized sheep RBCs is observed if rabbit anti-human albumin is used in place of anti-human IgG.

PRESENTATION:

10 ml of Human IgG Sensitized Glutaraldehyde-Stabilized sheep RBCs in saline with 0.1% NaN₃ (1%).

STORAGE:

Store at 4°C. DO NOT FREEZE.

APPLICATIONS:

This product is suitable for use in hemagglutination assays.

Optimum conditions for use of this product in analytical procedures should be established in each laboratory

Laboratory Reagent For Research Use Only

MW 02/09/15

Visit our website for your local distributor.

CEDARLANE[®] 

www.cedarlanelabs.com

An ISO 9001:2000 and ISO 13485:2003
registered company.

In CANADA: Toll Free: 1-800-268-5058

4410 Paletta Court, Burlington, ON L7L 5R2 ph: (289) 288-0001, fax: (289) 288-0020
e-mail: general@cedarlanelabs.com

In the USA: Toll Free: 1-800-721-1644

1210 Turrentine Street, Burlington, NC 27215 ph: (336) 513-5135, fax: (336) 513-5138
e-mail: service@cedarlanelabs.com