SZABO SCANDIC

Part of Europa Biosite

Produktinformation

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mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Technically
Speaking

CEDARLANE[®] 
www.cedarlanelabs.com

Conveniently Delivering You Today's Innovations
for the Science of Tomorrow™

Glutaraldehyde Stabilized, Tanned Sheep RBCs

CLD68LT

Lot: 1657

DESCRIPTION:

Cedarlane's **CLD68LT** are glutaraldehyde-stabilized, tanned freshly prepared sheep red blood cells. Glutaraldehyde stabilized, tanned sheep RBCs are produced from freshly washed sheep RBCs by brief exposure to a glutaraldehyde-buffer solution and exhaustive washing in saline. This is followed by brief exposure to a tannic acid-buffer solution and washing with saline. This treatment largely preserves the native antigenicity of the erythrocytes while rendering the cells generally resistant to lysis by osmotic shock, freeze-thawing or immune hemolysis. It also allows for coupling of purified serum proteins directly to the red blood cell surface. This reagent can be used directly in hemagglutination procedures, or may be coupled with various proteins for hemagglutination testing.

PRESENTATION:

10 mL of glutaraldehyde-stabilized, tanned sheep RBCs in saline with 0.1% NaN₃ (10%).

STORAGE:

Store at 4°C. DO NOT FREEZE.

APPLICATIONS:

This product is suitable for use in hemagglutination assays.

Optimum conditions for use of this product in analytical procedures should be established in each laboratory

Laboratory Reagent For Research Use Only

JK 03/03/15

Visit our website for your local distributor.

CEDARLANE[®] 

www.cedarlanelabs.com

An ISO 9001:2000 and ISO 13485:2003
registered company.

In CANADA: Toll Free: 1-800-268-5058

4410 Paletta Court, Burlington, ON L7L 5R2 ph: (289) 288-0001, fax: (289) 288-0020
e-mail: general@cedarlanelabs.com

In the USA: Toll Free: 1-800-721-1644

1210 Turrentine Street, Burlington, NC 27215 ph: (336) 513-5135, fax: (336) 513-5138
e-mail: service@cedarlanelabs.com