



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 


Go Alcohol-free and Eliminate Ongoing Costs with IPA-free Corning® CoolCell Freezing Systems

CORNING



Many labs use isopropanol (IPA) containers for cryogenic freezing, but they require costly alcohol replacements every 5 uses, can be cumbersome to handle, and may have inconsistent freezing rates.

Corning CoolCell is a reusable system that uniformly freezes your cells without the mess and cost of IPA. With CoolCell containers, you can depend on high reproducibility and high cell viability, to ensure you preserve the most cells possible for your research.

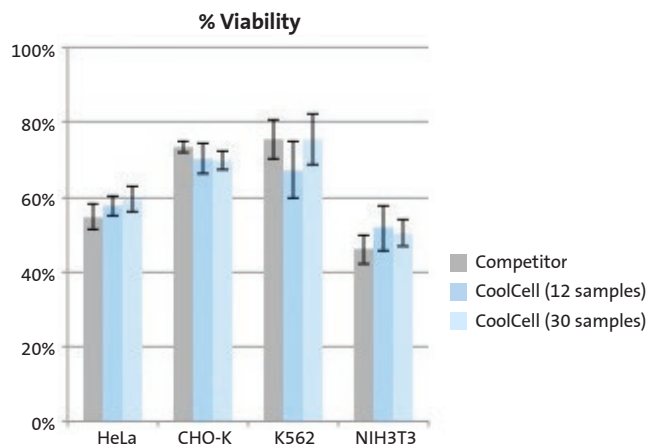

HeLa, CHO-K, K562, NIH3T3. Corning CoolCell 12-well, CoolCell FTS30 30-well, or competitor freezing containers were used to freeze all four cell lines. Identical transfection efficiencies and viabilities were observed after thawing.

Ensure temperature consistency and minimize contamination risk with Corning® sample cooling systems



To support consistent, reproducible, and standardized sample handling, Corning now offers solutions for sample cooling that address the pitfalls of temperature regulation and reduce contamination risk and variability which originates from tubes or wells placed directly into ice.

Corning CoolRack® thermo-conductive modules enable consistent, reproducible, standardized temperature control. These solutions may be used with Corning ice buckets and pans or if you prefer to go completely ice free, you can pair it with Corning CoolBox™ ice-free cooling systems.

GOOD



Corning Ice Bucket

BETTER



Corning CoolRack with Ice Bucket

BEST



Corning CoolRack with Corning CoolBox

For more specific information on claims, visit the Certificates page at www.corning.com/lifesciences.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

For additional product or technical information, visit www.corning.com/lifesciences or call 800.492.1110. Customers outside the United States, call +1.978.442.2200 or contact your local Corning sales office.

Corning Incorporated Life Sciences

836 North St.
Building 300, Suite 3401
Tewksbury, MA 01876
t 800.492.1110
t 978.442.2200
f 978.442.2476

www.corning.com/lifesciences

Worldwide Support Offices

ASIA/PACIFIC
Australia/New Zealand
t 61 427286832

China
t 86 21 3338 4338
f 86 21 3338 4300

India
t 91 124 4604000
f 91 124 4604099

Japan
t 81 3-3586 1996
f 81 3-3586 1291

Korea
t 82 2-796-9500
f 82 2-796-9300

Singapore
t 65 6572-9740
f 65 6861-2913

Taiwan
t 886 2-2716-0338
f 886 2-2516-7500

EUROPE

France
t 0800 916 882
f 0800 918 636

Germany
t 0800 101 1153
f 0800 101 2427

The Netherlands
t 31 20 655 79 28
f 31 20 659 76 73

United Kingdom
t 0800 376 8660
f 0800 279 1117

All Other European Countries

t 31 (0) 20 659 60 51
f 31 (0) 20 659 76 73

LATIN AMERICA
grupoLA@corning.com

Brasil
t (55-11) 3089-7400

Mexico
t (52-81) 8158-8400

CORNING | **FALCON** | **AXYGEN** | **GOSELIN** | **PYREX**

For a listing of trademarks, visit www.corning.com/clstrademarks.
All other trademarks are the property of their respective owners.