



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

**AMO (Amoxicillin) ELISA Kit**

Catalog No: E-FS-SP001

96T/96T\*3

This manual must be read attentively and completely before using this product.

If you have any problems, please contact our Technical Service Center for help.

Toll-free: 1-888-852-8623 Tel: 1-832-243-6086 Fax: 1-832-243-6017

Email: [techsupport@elabscience.com](mailto:techsupport@elabscience.com)

Website: [www.elabscience.com](http://www.elabscience.com)

Please kindly provide us the lot number (on the outside of the box) of the kit for more efficient service.

## Test principle

This kit uses Competitive-ELISA as the method for the quantitative detection. It can detect Amoxicillin (AMO) in samples, such as cell-culture, serum, water etc. This kit is composed of ELISA Microtiter plate, HRP conjugate, antibody working solution, standard and other supplementary reagents. The microtiter plate provided in this kit has been pre-coated with coupled antigen. During the detection, AMO in the samples or standard competes with coupled antigen on the solid phase supporter for sites of anti-AMO antibody. Then Horseradish Peroxidase (HRP) conjugate is added to each well, and substrate reagent is added for color development. There is a negative correlation between the OD value of samples and the concentration of AMO. The concentration of AMO in the samples can be calculated by comparing the OD of the samples to the standard curve.

## Technical indicator

**Reaction mode** (Incubation time and temperature): 25°C ; 30 min, 30 min, 15 min.

**Detection limit:** Cell-culture, Serum, Water, Vaccine---10 ppb.

*Note: If you need to test other samples, Please contact your area sales manager.*

### Cross-reactivity:

Amoxicillin	100%
Ampicillin	>100%
Cloxacillin	>100%
Oxacillin	85%
Cephacetrile	5%
Cephazoline	5%

**Sample recovery rate:** 90% ±30%.

**Kits components**

Item	Specifications
ELISA Microtiter plate	96 wells
Standard Liquid	1 mL each (ppb=ng/mL=ng/g) (0 ppb, 0.2 ppb, 0.6 ppb, 1.8 ppb, 5.4 ppb, 16.2 ppb)
Antibody working Solution	7 mL
HRP Conjugate	12 mL
5×Concentrated Sample Solution	20 mL
Substrate Reagent A	6 mL
Substrate Reagent B	6 mL
Stop Solution	6 mL
20×Concentrated Wash Buffer	25 mL
Plate Sealer	3 pieces
Sealed Bag	1 piece
Manual	1 copy

Note: All reagent bottle caps must be tightened to prevent evaporation and microbial pollution

**Other materials required but not supplied**

**Instrument:** Microplate reader, Printer, Homogenizer, Oscillators, Vortex mixer, Centrifuge, Balance (sensitivity 0.01g).

**Micropipette:** Single channel (20-200 µL, 100-1000 µL), Multichannel (30-300 µL).

**Reagents:** Methanol

---

**Notes**

1. The overall OD value will be lower when reagents have not been brought to room temperature before use or room temperature is below 25°C.
2. If the wells turn dry during the washing procedure, it will lead to bad linear standard curve and poor repeatability. Operate the next step immediately after wash.
3. Mix thoroughly and wash the plate completely. The consistency of wash procedure can strongly affect the reproducibility of this ELISA kit.
4. FOR RESEARCH USE ONLY. ELISA Microtiter plate should be covered by plate sealer. Avoid the kit to strong light.
5. **Each reagent is optimized for use in the E-FS-SP001. Do not substitute reagents from any other manufacturer into the test kit. Do not combine reagents from other E-FS-SP001 with different lot numbers.**
6. Substrate Reagent should be abandoned if it turns blue color. When OD value of standard (concentration: 0) < 0.8 unit (A450nm < 0.8), it indicates the reagent may be deteriorated.
7. Stop solution is caustic, avoid contact with skin and eyes.
8. As the OD values of the standard curve may vary according to the conditions of the actual assay performance (e.g. operator, pipetting technique, washing technique or temperature effects), the operator should establish a standard curve for each test.
9. Even the same operator might get different results in two separate experiments. In order to get reproducible results, the operation of every step in the assay should be controlled.
10. **For mentioned sample fast and efficient extraction methods are included in the kit description. Please consult technical support for the applicability if other sample need to be tested.**
11. The kit is used for rapid screening of actual samples. If the test result is positive, the instrument method such as HPLC, LC/MS, etc. can be used for quantitative confirmation.

**Storage and expiry date**

Store the kit at 2-8°C. Do not freeze any test kit components.

Return any unused microwells to their original foil bag and reseal them together with the desiccant provided and further store at 2-8°C.

**Expiry date:** expiration date is on the packing box.

## Experimental preparation

Restore all reagents and samples to room temperature before use.

Open the microplate reader in advance, preheat the instrument, and set the testing parameters.

### 1. Sample pretreatment Notice:

Experimental apparatus should be clean, and the pipette should be disposable to avoid cross-contamination during the experiment.

### 2. Solution preparation

*Please prepare solution according to the number of samples. Don't use up all components in the kit at once!*

Solution 1: 70% Methanol (*for muscle, egg sample*)

**Methanol (V):** Deionized water (V) = 7:3.

Solution 2: Sample Solution (*for muscle, egg sample*)

Dilute **5×Concentrated Sample Solution** with deionized water. (5×Concentrated Sample Solution (V): Deionized water (V) = 1:4).

Solution 3: Wash Buffer

Dilute **20×Concentrated Wash Buffer** with deionized water. (20×Concentrated Wash Buffer (V): Deionized water (V) = 1:19).

### 3. Sample pretreatment procedure

#### 1.1 Pretreatment of serum (livestock), cell-culture, water sample:

- (1) Use the conventional method to prepare sample, the sample must be clear, no hemolysis and no pollution. Samples can be conserved at 2-8°C in 1 week, and it should be stored at -20°C for a long term storage.
- (2) Weigh 2±0.05 g of homogenate sample into a 50 mL centrifuge tube. Add 4 mL of **70% Methanol** (Solution 1), vortex for 1 min.
- (3) Centrifuge for 5 min at 4000 rpm.
- (4) Take 50 µL of supernatant to another centrifuge tube, then add 450 µL of **Sample Solution** (Solution 2), vortex for 10 s.
- (5) Take 50 µL of liquid for analysis.

**Note: Sample dilution factor: 20, detection limit: 10 ppb**

---

**Assay procedure**

Restore all reagents and samples to room temperature (25°C) before use. All the reagents should be mixed thoroughly by gently swirling before pipetting. Avoid foaming. The unused ELISA Microtiter plate should be sealed as soon as possible and stored at 2-8°C.

1. **Number:** number the sample and standard in order (multiple well), and keep a record of standard wells and sample wells. **Standard and Samples need test in duplicate.**
2. **Add Sample:** add 50 µL of **Standard** or **Sample** per well, add 50 µL of **Antibody Working Solution** to each well. Gently oscillate for 10 s to mix thoroughly and cover the plate with sealer. Incubate for 30 min at 25°C.
3. **Wash:** uncover the sealer carefully, remove the liquid in each well. Immediately add 260 µL of **Wash Buffer** (Solution 3) to each well and wash. Repeat the wash procedure for 4 times, 30 s intervals/time. Invert the plate and pat it against thick clean absorbent paper (If bubbles exist in the wells, clean tips can be used to prick them).
4. **HRP Conjugation:** add 100 µL of **HRP Conjugate** into each well, and incubate at 25°C for 30 min.
5. **Wash:** Repeat step 3 for washing.
6. **Color Development:** add 50 µL of **Substrate Reagent A** to each well, and then add 50 µL of **Substrate Reagent B**. Gently oscillate for 10 s to mix thoroughly. Incubate for 15 min at 25°C in the dark (The reaction time can be extended according to the actual color change).
7. **Stop Reaction:** add 50 µL of **Stop Solution** to each well, oscillate gently for 10 s to mix thoroughly.
8. **OD Measurement:** determine the optical density (OD value) of each well at 450 nm (reference wavelength 630 nm) with a microplate reader. This step should be finished in 5 min after stop reaction.

## Result analysis

### 1. Absorbance(%)= $A/A_0 \times 100\%$

A: Average absorbance of standard or sample

$A_0$ : Average absorbance of 0 ppb Standard

### 2. Drawing and calculation of standard curve

Create a standard curve by plotting the absorbance percentage of each standard on the y-axis against the log concentration on the x-axis to draw a semi-logarithmic plot. Add average absorbance value of sample to standard curve to get corresponding concentration. **If samples have been diluted, the concentration calculated from the standard curve must be multiplied by the dilution factor.**

For this kit, it is more convenient to use professional analysis form for accurate and fast analysis of batch samples.

