



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

(FOR RESEARCH USE ONLY. DO NOT USE IT IN CLINICAL DIAGNOSIS !)

2-step Plus Poly-HRP Anti Rabbit IgG Detection System (with DAB solution)

Catalog No: E-IR-R215

3 mL/ 6 mL/ 18 mL

Cat	product	3 mL	6 mL	18 mL	Storage
E-IR-R215A	Normal Goat Serum (Ready-to-Use)	3 mL	6 mL	18 mL	2~8 ℃
E-IR-R215B	Polyperoxidase-anti-Rabbit IgG (Ready-to-Use)	3 mL	6 mL	18 mL	2~8 ℃
E-IR-R215C	3% H ₂ O ₂	3 mL	6 mL	18 mL	2~8 ℃
E-IR-R215D	DAB Concentrate (20×)	150 μL	300 μL	900 μL	2~8 ℃
E-IR-R215E	DAB Substrate	3 mL	6 mL	18 mL	2~8 ℃
Manual		One Copy			

This manual must be read attentively and completely before using this product.

If you have any problems, please contact our Technical Service Center for help.

Phone: 240-252-7368(USA) 240-252-7376(USA)

Email: techsupport@elabscience.com

Website: www.elabscience.com

Please kindly provide us the lot number (on the outside of the box) of the kit for more efficient service.

Introduction

2-step plus is a two-step immunohistochemical broad spectrum detection reagent. It polymerizes monovalent Fab fragments of secondary antibody and enzyme, which replaces the secondary antibody and tertiary antibody in traditional method, can directly amplify the binding signal of antibody-antigen. This method not only retains the specific binding ability of antibody with antigen, but also can effectively avoid space steric hindrance caused by excessive polymer molecules. Compared with the traditional SP three-step method, this kit has the characteristics of simple, rapid, high-sensitivity. This system abandons the using of biotin, so it can avoid background staining by endogenous biotin. It can be used in IHC, in which the primary antibody is monoclonal/polyclonal antibody derived from rabbit.

Diluent for DAB concentrated solution has been provided in this kit to avoid the influence of the different water acidity and alkalinity on the DAB Chromogenic Agent.

Sample dyeing

1. Dewax and hydrate the paraffin section.
2. Make thermal repair or digestion treatment to antigen of the tissue section if necessary according to antigen/antibody situation.
3. Incubate with E-IR-R215C (3% H₂O₂) for 10 min to eliminate endogenous peroxidase activity. Wash with PBS or TBS, 2 min×3 times.
4. Add E-IR-R215A -Normal Goat Serum (Ready-to-Use), incubate at 37 °C for 30 min. Shake off any excess liquid.
5. Add primary antibody (Rabbit-IgG) with proper dilution ratio, incubate at 20~37 °C for 1~2h or at 4 °C overnight (then rewarm at 37 °C for 30 min). Wash with PBS or TBS, 2 min×3 times. Dry the section with absorbent paper.
6. Add E-IR-R215B (Polyperoxidase-anti-Rabbit IgG), incubate at room temperature or 37 °C for 20 min. Wash with PBS or TBS, 2 min×3 times.
7. Add 1 drop (approximately 50 μL) of E-IR-R215D (DAB Concentrate) into each 1 mL of E-IR-R215E (DAB Substrate), mix fully and the mixed reagent is the DAB Working Solution. Prepare fresh solution before use and the prepared solution should be stored in the dark. Fresh prepared DAB Working Solution is valid within 4 hours and the unused solution must be abandoned
8. Take control of the DAB coloration period, the color of tan or brownish yellow is the positive signal. Avoid of excessive reaction.
9. Wash the section with deionized water terminate the chromogenic reaction, then operate the procedures of counterstaining, dehydrating, transparentizing and sealing

Storage

Store at 2~8 °C, shading light. Avoid of freezing. Valid for 12 months. The reagents are valid within 6 months after opening.