

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten! See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere Liefer- und Versandbedingungen

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

linkedin.com/company/szaboscandic in





AFM1 (Aflatoxin M1) ELISA Kit

Catalog No: E-TO-E007

96T/96T*3

Version Number: V1.2
Replace version: V1.1

Revision Date: 2024.03.14

This manual must be read attentively and completely before using this product.

If you have any problems, please contact our Technical Service Center for help.

Toll-free: 1-888-852-8623 Tel: 1-832-243-6086 Fax: 1-832-243-6017

Email: techsupport@elabscience.com

Website: www.elabscience.com

Please kindly provide us the lot number (on the outside of the box) of the kit for more efficient service.



Test principle

This kit uses Competitive-ELISA as the method for the quantitative detection. It can detect Aflatoxin M1 (AFM1) in samples, such as milk, milk powder. This kit is composed of ELISA Microtiter plate, HRP conjugate, antibody working solution, standard and other supplementary reagents. The microtiter plate in this kit has been pre-coated with coupled antigen. During the reaction, AFM1 in the samples or standard competes with coupled antigen on the solid phase supporter for sites of anti-AFM1 antibody. Then Horseradish Peroxidase (HRP) conjugate is added to each microtiter plate well, and substrate reagent is added for color development. There is a negative correlation between the OD value of samples and the concentration of AFM1. The concentration of AFM1 in the samples can be calculated by comparing the OD of the samples to the standard curve.

Technical indicator

Reaction mode (Incubation time and temperature): 25°C; 30 min, 15 min. **Detection limit:** Milk---0.1 ppb; Milk powder---0.15 ppb; Urine---0.5 ppb

Cross-reactivity: Aflatoxin M1 ---100%

Sample recovery rate: Milk ---85% \pm 15%; Milk powder, Urine---80% \pm 15%.

Kits components

Item	Specifications
ELISA Microtiter plate	96 wells
Standard Liquid	1 mL each (ppb=ng/mL=ng/g) (0 ppb, 0.05 ppb, 0.15 ppb, 0.45 ppb, 1.35 ppb, 4.05 ppb)
HRP Conjugate	5.5 mL
Antibody Working Solution	5.5 mL
Substrate Reagent A	6 mL
Substrate Reagent B	6 mL
Stop Solution	6 mL
20×Concentrated Wash Buffer	40 mL
2×Concentrated Reconstitution Buffer	50 mL
Plate Sealer	3 pieces
Sealed Bag	1 piece
Manual	1 copy

Note: All reagent bottle caps must be tightened to prevent evaporation and microbial pollution.



Other materials required but not supplied

Instrument: Microplate reader, Printer, Homogenizer, Nitrogen Evaporators, Water bath, Vortex mixer, Centrifuge, Graduated pipette, Balance (sensibility 0.01 g).

High-precision transferpettor: Single channel (20-200 μL, 100-1000 μL), Multichannel (300 μL).

Reagents: Acetonitrile.

Notes

- 1. The overall OD value will be lower when reagents have not been brought to room temperature before use or room temperature is below 25°C.
- 2. If the wells turn dry during the washing procedure, it will lead to bad linear standard curve and poor repeatability. Operate the next step immediately after wash.
- 3. Mix thoroughly and wash the plate completely. The consistency of wash procedure can strongly affect the reproducibility of this ELISA kit.
- 4. FOR RESEARCH USE ONLY. ELISA Microtiter plate should be covered by plate sealer. Avoid the kit to strong light.
- 5. Each reagent is optimized for use in the E-TO-E007. Do not substitute reagents from any other manufacturer into the test kit. Do not combine reagents from other E-TO-E007 with different lot numbers.
- 6. Substrate Reagent should be abandoned if it turns blue color. When OD value of standard (concentration: 0 < 0.5 unit ($A_{450 \text{ nm}} < 0.5$), it indicates reagent is deteriorated.
- 7. Stop solution is caustic, avoid contact with skin and eyes.
- 8. As the OD values of the standard curve may vary according to the conditions of the actual assay performance (e.g. operator, pipetting technique, washing technique or temperature effects), the operator should establish a standard curve for each test.
- 9. Even the same operator might get different results in two separate experiments. In order to get reproducible results, the operation of every step in the assay should be controlled.
- 10. For mentioned sample fast and efficient extraction methods are included in the kit description. Please consult technical support for the applicability if other sample need to be tested.
- 11. The kit is used for rapid screening of actual samples. If the test result is positive, the instrument method such as HPLC, LC/MS, etc. can be used for quantitative confirmation.

Storage and expiry date

Store the kit at $2\text{-}8^{\circ}\text{C}$. Do not freeze any test kit components.

Return any unused microwells to their original foil bag and reseal them together with the desiccant provided and further store at $2-8\,^{\circ}\mathrm{C}$.

Expiry date: expiration date is on the packing box.



Experimental preparation

Restore all reagents and samples to room temperature before use.

Open the Microplate reader in advance, preheat the instrument, and set the testing parameters.

1. Sample pretreatment Notice:

Experimental apparatus should be clean, and the pipette should be disposable to avoid cross-contamination during the experiment.

2. Solution preparation

Please prepare solution according to the number of samples. Don't use up all components in the kit at once!

Solution 1: Reconstitution Buffer

Dilute the **2**×Concentrated Reconstitution Buffer with deionized water.

 $(2 \times Concentrated Reconstitution Buffer (V): Deionized water (V) = 1:1).$

Solution 2: 84% Acetonitrile (for milk powder sample)

Acetonitrile (V): Deionized water (V) =84:16.

Solution 3: Wash Buffer

Dilute the 20 × Concentrated Wash Buffer with deionized water.

 $(20 \times Concentrated Wash Buffer (V): Deionized water (V) = 1:19).$

3. Sample pretreatment procedure

3.1. Pretreatment of milk sample:

- (1) Take 1 mL of liquid milk into 50 mL centrifuge tube, add 4 mL of **Acetonitrile**, oscillate for 5 min, centrifuge at 4000 rpm for 10 min at room temperature.
- (2) Take 2.5 mL of supernatant to another centrifuge tube, dry with nitrogen evaporators or water bath at 50-60 °C. (Please do it in a ventilated environment)
- (3) Dissolve the residual with 1 mL of **Reconstitution Buffer** (Solution 1), vortex and mix fully.
- (4) Take 50 μL for analysis.

Note: Sample dilution factor: 2, detection limit: 0.1 ppb

3.2. Pretreatment of milk powder sample:

- (1) Weight 5 ± 0.05 g of milk powder into 50 mL centrifuge tube, add 20 mL of **84% Acetonitrile** (Solution 2), vortex for 5 min, centrifuge at 4000 rpm for 10 min at room temperature.
- (2) Take 1 mL clear liquid to another centrifuge tube, dry with nitrogen evaporators or water bath at 50-60°C.
- (3) Dissolve the residual with 750 µL of **Reconstitution Buffer** (Solution 1), vortex and mix fully.
- (4) Take 50 μL for analysis.

Note: Sample dilution factor: 3, detection limit: 0.15 ppb



3.3. Pretreatment of urine sample:

- (1) Take 100 μL of clarified fresh sample to detect (the turbid urine sample should be filtered or centrifuge at 4000 rpm for 5 min to get clear urine sample).
- (2) Add 900 μL of **Reconstitution Buffer** (Solution 1), mix fully. Samples can be stored at 2-8°C for one week. If samples will be not tested in a week, store them below -20°C, and please try to avoid freeze-thaw cycles.
- (3) Take $50 \mu L$ of the supernatant for analysis.

Note: Sample dilution factor: 10, detection limit: 0.5 ppb

Assay procedure

Restore all reagents and samples to room temperature (25°C) before use. All the reagents should be mixed thoroughly by gently swirling before pipetting. Avoid foaming. The unused ELISA Microtiter plate should be sealed together to sealed bag with the desiccant provided and stored at 2-8°C.

- 1. **Number:** number the sample and standard in order (multiple well), and keep a record of standard wells and sample wells. **Standard and Samples need test in duplicate.**
- 2. Add sample: add 50 μL of Standard or Sample per well, then add 50 μL of HRP Conjugate to each well, then add 50 μL of Antibody Working Solution, cover the plate with sealer, oscillate for 5 sec gently to mix thoroughly, incubate at 25 °C for 30 min in shading light.
- 3. **Wash:** uncover the sealer carefully, remove the liquid of each well. Immediately add 300 µL of **Wash Buffer** (Solution 3) to each well. Repeat wash procedure for 5 times, 30 s intervals/time. Invert the plate and pat it against thick clean absorbent paper (If bubbles exist in the wells, clean tips can be used to prick them).
- 4. **Color Development:** add 50 μL of **Substrate Reagent A** to each well, and then add 50 μL of **Substrate Reagent B**. Gently oscillate for 5 s to mix thoroughly. Incubate at 25°C for 15 min in shading light. (The reaction time can be extended according to the actual color change).
- 5. **Stop reaction:** add 50 μL of **Stop Solution** to each well, oscillate gently to mix thoroughly.
- 6. **OD Measurement:** determine the optical density (OD value) of each well at 450 nm (reference wavelength 630 nm) with a microplate reader. This step should be finished in 10 min after stop reaction.



Result analysis

1. Absorbance (%)= $A/A_0 \times 100\%$

A: Average absorbance of standard or sample

A₀: Average absorbance of 0 ppb Standard

2. Drawing and calculation of standard curve

Create a standard curve by plotting the absorbance percentage of each standard on the y-axis against the log concentration on the x-axis to draw a semi-logarithmic plot. Add average absorbance value of sample to standard curve to get corresponding concentration. **If samples have been diluted, the concentration calculated from the standard curve must be multiplied by the dilution factor.**

For this kit, it is more convenient to use professional analysis form for accurate and fast analysis on a large number of samples.



